



SERVING SCIENTISTS
SINCE 1984

CERTIFIED CLASS 7 CLEAN ROOM
MANUFACTURING

PRODUCTION IN
THE NETHERLANDS AND
THE UNITED STATES

SAMPLE PRESERVATION & STORAGE SOLUTIONS

Advancing research by serving scientists with sample preservation and storage

ABOUT MICRONIC

Our goal is to advance research by serving scientists in finding solutions that contribute to a higher quality of life. We develop and manufacture a range of Dutch-designed products to enhance the process of sample preservation and storage.

Micronic is an independent organization with its headquarters located in Lelystad, the Netherlands. At Micronic, we develop and manufacture our premium labware and precision engineered equipment in-house. We are committed to offering top-of-the-line products, delivering exceptional quality, reliability and durability. Above all, we stand behind our solutions and are confident that you will too. Our products are applied worldwide in the research laboratories of university hospitals, forensics, agricultural, veterinary and governmental institutes, as well as companies in biotech, food, chemical and pharmaceutical industries.

PRODUCTION AND QUALITY CHARACTERISTICS

- **Environmental commitment:** Micronic is aware of its interconnect-edness with the environment. We continuously think about how we can operate our business in a way that is friendly to this natural environment and sustainable to the earth, and we act accordingly. An example of our environmental commitment can be found in the ACT label. This label is the first environmental impact factor label for laboratory products. Micronic participates in this great program so that purchasers can make informed decisions about a product's environmental impact and to promote sustainability in the lab.
- **ISO 9001:2015 and 14001:2015 certified:** All Micronic's businesses, located in the Netherlands and the United States are certified for ISO 9001:2015. Micronic Manufacturing, which produces the sample storage labware is certified for ISO 14001:2015.



- **Class 7 clean room:** Micronic labware is manufactured and assembled in a NEN-EN-ISO 14644-1 Class 7 certified clean room environment (comparable to US Federal Standard 209E Class 10,000).
- **FDA approved and virgin material:** Materials are FDA approved and comply with the (USA) USP XXII tests, including class VI (Medical approved), and the European Pharmacopoeia-monograph 3.2.2 unless communicated otherwise.
- **Extractables:** The materials used for Micronic tubes are the result of extensive research. One selection criterion is that no measurable amount of extractables should leach into samples.
- **RNase and DNase:** By maintaining the highest possible hygiene level during our production processes, Micronic manufactured labware is free of any detectable RNase or DNase contamination.
- **Endotoxin (pyrogen):** By maintaining the highest possible hygiene level during its production processes, Micronic is able to limit the endotoxin level of produced and packaged labware to an acceptable minimum (< 0.01 EU/ml).
- **CE label:** Micronic labware products are developed and manufactured in compliance with the European Directive 98/79/EC, Annex III.
- **Sterilization:** Micronic products can be supplied gamma irradiated or ethylene oxide treated.



CONTENTS

| | |
|---|-----------|
| SAMPLE STORAGE TUBES | 4 |
| Superior features | 5 |
| Tube overview and specifications | 6 |
| Tubes with internal thread (96-well, 48-well and 24-well format) | 11 |
| Tubes with external thread (96-well, 48-well and 24-well format) | 18 |
| Tubes with external thread hybrid (96-well, 48-well and 24-well format) | 28 |
| Special tubes (tubestrip-8 and light sensitive tubes) | 35 |
| SAMPLE STORAGE RACKS | 44 |
| Superior features | 45 |
| Rack overview and specifications | 46 |
| 96-well format racks | 50 |
| 48-well format racks | 54 |
| 24-well format racks | 56 |
| 138-well format racks | 58 |
| SAMPLE STORAGE CAPS | 60 |
| Superior features | 61 |
| Cap overview and specifications | 62 |
| Screw Caps | 64 |
| TPE Push Caps | 66 |
| EVA Push Caps | 67 |
| SAMPLE HANDLING EQUIPMENT | 70 |
| Equipment Service & Warranty | 71 |
| Code Reader overview and specifications | 72 |
| Tube Readers - Single Tube | 74 |
| Rack Readers - Single Rack | 76 |
| Push Cap Cappers & Decapper | 78 |
| Screw Cap Recappers | 81 |
| Tube Handler | 85 |
| SAMPLE STORAGE STARTER PACKS | 90 |
| Overview Starter Packs | 92 |
| INNOVATION AT MICRONIC | 94 |



SERVING SCIENTISTS
SINCE 1984

CERTIFIED CLASS 7 CLEAN ROOM
MANUFACTURING

SAMPLE STORAGE TUBES

MICRONIC OFFERS A COMPLETE LINE OF SAMPLE STORAGE TUBES THAT FIT IN THE 138-WELL, 96-WELL, 48-WELL OR 24-WELL ANSI / SLAS STANDARD (FORMERLY RECOGNIZED AS THE ANSI / SBS STANDARD) RACK FORMAT. THE COMPREHENSIVE RANGE INCLUDES TUBES WITH VOLUMES RANGING FROM 0.30ML TO 6.00ML WHICH ARE SUITABLE FOR THE MOST COMMON LABORATORY STORAGE APPLICATIONS. THE TUBES ARE AVAILABLE NON-CODED, ALPHANUMERIC CODED OR WITH A UNIQUE 2D DATA-MATRIX CODE (ADHERE TO THE ECC200 STANDARD) ON THE TUBE BOTTOM, WHICH CAN BE COMBINED WITH A 1D BARCODE ON THE TUBE SIDE. IN ADDITION, MICRONIC OFFERS A VARIETY OF SPECIAL TUBES INCLUDING THE TUBESTRIP-8 AND AMBER TUBES FOR LIGHT SENSITIVE SAMPLES. THROUGH OUR LINE OF COMPATIBLE STORAGE RACKS, SEALING PRODUCTS AND EQUIPMENT MICRONIC IS ABLE TO OFFER YOU A COMPLETE SOLUTION.

SUPERIOR FEATURES OF MICRONIC SAMPLE STORAGE TUBES



UNIQUE MANUFACTURING PROCESS











All Micronic sample storage tubes are manufactured to industry leading strict tolerances. Due to the unique 2K injection molding technique, the highly transparent tube and the code surface are molded to one component leading to the following advantages:

- Code surface cannot be separated from the transparent tube enabling absolute traceability
- Solid and uniform walls allow maximum working volume
- No contamination during the production process


MICRONIC'S HIGH QUALITY AND STRONG SAMPLE STORAGE TUBES ARE ABLE TO WITHSTAND:

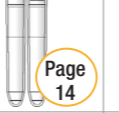- *Multiple freeze-thaw cycles.*
Extensive freeze-thaw tests are conducted with Micronic sample storage tubes and results show no leakages for both Micronic Push Cap tubes and Micronic Screw Cap tubes closed with TPE caps. Additionally, research also shows that no Micronic TPE caps pop off during the freeze-thaw process (reference: Test Report TR021901).
- *Grinding of samples by high speed bead beating for DNA/RNA extraction.*
When homogenizing samples, beads are often used to break open cell walls and membranes to release the desired DNA and RNA. The solid walls of Micronic tubes have been tested and are able to withstand the process of bead beating easily (reference: Test Report TR021801 and TR022701).

TUBE OVERVIEW AND SPECIFICATIONS - INTERNAL THREAD

| | 0.50 ML | 0.75 ML | 0.75 ML | 1.10 ML | 1.40 ML | 1.40 ML | 1.40 ML | 2.00 ML | 4.00 ML | 6.00 ML |
|--|---|---|---|--|---|---|---|---|---|---|
| |  Page 11 |  Page 12 |  Page 12 |  Page 13 |  Page 14 |  Page 14 |  Page 14 |  Page 15 |  Page 16 |  Page 17 |
| Storage rack | Loborack-96 | Loborack-96 | Loborack-96 | Roborack-96 | Comorack-96 | Roborack-96 | Roborack-96 | Roborack-96 | Micronic 48-4 | Micronic 24-4 |
| Ultra low temperature rack | Micronic 96-1/Q1 | Micronic 96-2 | Micronic 96-2 | Micronic 96-3 | Micronic 96-4 | Micronic 96-4 | Micronic 96-4 | Micronic 96-5 | | |
| Thread type | Internal | Internal | n/a | Internal | n/a | n/a | Internal | Internal | Internal | Internal |
| Tube type | Screw Cap tube | Screw Cap tube | Push Cap tube | Screw Cap tube | Push Cap tube | Push Cap tube | Screw Cap tube | Screw Cap tube | Screw Cap tube | Screw Cap tube |
| Inner tube shape | V | V | V or U | V | U | V or U | V or U | U | Flat | Flat |
| Working volume with Push Cap (+21°C)* | 0.30ml | 0.58ml | 0.56 or 0.59ml | 0.90ml | 1.14ml | 1.13 or 1.19ml | 1.15 or 1.20ml | 1.58ml | n/a | n/a |
| Working volume with Push Cap (-20°C)* | 0.30ml | 0.53ml | 0.53 or 0.53ml | 0.79ml | 1.01ml | 1.04 or 1.05ml | 1.01 or 1.05ml | 1.39ml | n/a | n/a |
| Working volume with Push Cap (-80°C)* | 0.30ml | 0.55ml | 0.51 or 0.55ml | 0.80ml | 0.95ml | 1.01 or 0.99ml | 0.95 or 1.06ml | 1.40ml**** | n/a | n/a |
| Working volume with Screw Cap (+21°C)* | 0.21ml | 0.48ml | n/a | 0.80ml | n/a | n/a | 1.04 or 1.10ml | 1.49ml | 2.70ml | 4.80ml |
| Working volume with Screw Cap (-20°C)* | 0.21ml | 0.48ml | n/a | 0.78ml | n/a | n/a | 1.00 or 1.06ml | 1.38ml | 2.65ml | 4.10ml |
| Working volume with Screw Cap (-80°C)* | 0.21ml | 0.48ml | n/a | 0.73ml | n/a | n/a | 1.01 or 0.98ml | 1.25ml**** | 2.45ml | 3.60ml |
| 2D Data-Matrix coded (bottom) | Yes | Yes | Yes | Yes | No | Yes | Yes | Yes | Yes | Yes |
| Alphanumeric coded (bottom) | Yes | Yes | Yes | No | Yes | Yes | Yes | Yes | Yes | No |
| Non-coded (bottom) | Yes | Yes | Yes | Yes | Yes | Yes | Yes | No | Yes | Yes |
| Human readable coded (bottom) | Yes | Yes | Yes | Yes | No | Yes | Yes | Yes | Yes | Yes |
| Capping options (SC = Screw Cap) | Low Profile SC Screw Cap Push Cap | Low Profile SC Screw Cap Push Cap | Push Cap | Low Profile SC Screw Cap Push Cap | Push Cap | Push Cap | Low Profile SC Screw Cap Push Cap | Low Profile SC Screw Cap Push Cap | Screw Cap | Screw Cap |
| Outer diameter | 9.0mm | 9.0mm | 9.0mm | 9.0mm | 9.0mm | 9.0mm | 9.0mm | 9.0mm | 13.0mm | 16.3mm |
| Tube height | 18.5mm | 26.0mm | 26.0mm | 36.5mm | 44.0mm | 44.0mm | 44.0mm | 59.0mm | 44.0mm | 43.4mm |
| Tube height with Push Cap** | 19.5mm | 27.0mm | 27.0mm | 37.5mm | 45.0mm | 45.0mm | 45.0mm | 60.0mm | n/a | n/a |
| Tube height with Screw Cap | 26.0mm | 33.5mm | n/a | 44.0mm | n/a | n/a | 51.5mm | 66.5mm | 51.0mm | 51.4mm |
| Available precapped | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Twist-lock | Yes | Yes | No | Yes | No | No | Yes | Yes | Yes | Yes |
| Snap tube option | Yes | Yes | Yes | Yes | No | Yes (V-bottom) | Yes | Yes | No | No |
| Sterilization option | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| DNase / RNase free | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Class 7 clean room production | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Minimum storage temperature*** | Vapor phase LN2 | Vapor phase LN2 | -80°C | Vapor phase LN2 | -80°C | -80°C | Vapor phase LN2 | -80°C**** | Vapor phase LN2 | Vapor phase LN2 |
| Availability | Bulk, Rack | Bulk, Refill, Rack | Bulk, Refill, Rack | Bulk, Rack | Bulk, Refill, Rack | Bulk, Refill, Rack | Bulk, Refill, Rack | Bulk, Rack | Bulk, Rack | Bulk, Rack |

TUBE RANGE INTERNAL THREAD

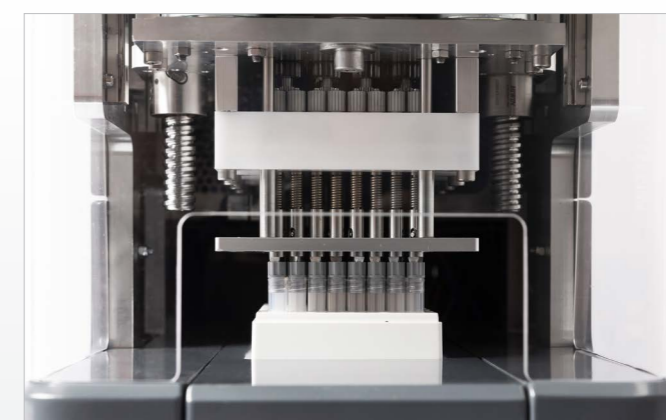











* If more sample volume is used in a tube than the specified working volume, it could be possible that deformation of the tube occurs. Request Test Report TR023201 for more information.

** Also applicable for the Low Profile Screw Cap.

*** Temperature range of the material of the tubes, caps and racks is approved from -80°C to +100°C. In practice tubes are used in lower temperatures until vapor phase LN2 (-150°C to -196°C). Micronic does not recommend to store tubes in liquid phase nitrogen (LN2). In practice tubes are only snap frozen in LN2 (without submerging the cap).

**** For -80°C storage of the 2.00ml tube its recommended to always use a rack with cover.

TUBE OVERVIEW AND SPECIFICATIONS - EXTERNAL THREAD

| | 0.30 ML | 0.40 ML | 0.75 ML | 0.80 ML | 1.10 ML | 1.40 ML | 2.00 ML | 1.00 ML | 4.00 ML | 3.50 ML |
|--|--|--|--|---|--|--|--|--|--|--|
| Ultra low temperature rack |  Page 18 Micronic 96-1 |  Page 19 Micronic 96-1 |  Page 20 Micronic 96-2 |  Page 21 Micronic 96-3 |  Page 22 Micronic 96-3 |  Page 23 Micronic 96-4 |  Page 24 Micronic 96-5 |  Page 25 Micronic 48-1 |  Page 26 Micronic 48-4 |  Page 27 Micronic 24-2 |
| Thread type | External | External | External | External | External | External | External | External | External | External |
| Tube type | Screw Cap tube | Screw Cap tube | Screw Cap tube | Screw Cap tube | Screw Cap tube | Screw Cap tube | Screw Cap tube | Screw Cap tube | Screw Cap tube | Screw Cap tube |
| Inner tube shape | V / Vr | V | V | V | V | V | V | Flat | Flat | Flat |
| Working volume with Push Cap (+21°C)* | 0.25ml / 0.20ml | 0.38ml | 0.57ml | 0.68ml | 0.85ml | 1.03ml | 1.48ml | n/a | n/a | n/a |
| Working volume with Screw Cap (+21°C)* | 0.32ml / 0.23ml | 0.40ml | 0.60ml | 0.75ml | 0.93ml | 1.11ml | 1.55ml | 0.85ml | 3.40ml | 3.35ml |
| Working volume with Screw Cap (-20°C)* | 0.25ml / 0.23ml | 0.38ml | 0.55ml | 0.65ml | 0.87ml | 1.05ml | 1.25ml | 0.85ml | 3.00ml | 3.05ml |
| Working volume with Screw Cap (-80°C)* | 0.25ml / 0.23ml | 0.36ml | 0.55ml | 0.65ml | 0.77ml | 0.94ml | 1.25ml | 0.85ml | 3.00ml | 3.05ml |
| 2D Data-Matrix coded (bottom) | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Alphanumeric coded (bottom) | No | No | No | No | No | No | No | No | No | No |
| Non-coded (bottom) | No | No | Yes | No | No | Yes | No | No | No | No |
| Human readable coded (bottom) | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Capping options | Screw Cap Screw Cap IG Push Cap | Screw Cap Screw Cap IG Push Cap | Screw Cap Screw Cap IG Push Cap | Screw Cap Screw Cap IG Push Cap | Screw Cap Screw Cap IG Push Cap | Screw Cap Screw Cap IG Push Cap | Screw Cap Screw Cap IG Push Cap | Screw Cap | Screw Cap | Screw Cap |
| Outer diameter | 8.8mm | 8.8mm | 9.0mm | 8.8mm | 9.0mm | 9.0mm | 9.0mm | 13.0mm | 13.0mm | 16.3mm |
| Tube height | 16.4mm | 21.8mm | 30.0mm | 34.8mm | 41.0mm | 48.0mm | 63.0mm | 16.0mm | 48.0mm | 30.0mm |
| Tube height with Push Cap | 17.8mm | 23.2mm | 31.4mm | 36.2mm | 42.7mm | 49.4mm | 64.4mm | n/a | n/a | n/a |
| Tube height with Screw Cap IG | 18.7mm | 24.1mm | 32.3mm | 37.0mm | 43.6mm | 50.3mm | 65.3mm | n/a | n/a | n/a |
| Tube height with Screw Cap | 19.9mm | 25.3mm | 33.5mm | 38.3mm | 44.6mm | 51.5mm | 66.5mm | 19.7mm | 51.5mm | 33.5mm |
| Available precapped | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Twist-lock | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Snap tube option | No | No | No | No | No | No | No | No | No | No |
| Sterilization option | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| DNase / RNase free | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Class 7 clean room production | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Minimum storage temperature** | Vapor phase LN2 | Vapor phase LN2 | Vapor phase LN2 | Vapor phase LN2 | Vapor phase LN2 | Vapor phase LN2 | -80°C*** | Vapor phase LN2 | Vapor phase LN2 | Vapor phase LN2 |
| Availability | Bulk, Rack | Bulk, Rack | Bulk, Rack | Bulk, Rack | Bulk, Rack | Bulk, Rack | Bulk, Rack | Bulk, Rack | Bulk, Rack | Bulk, Rack |

TUBE RANGE EXTERNAL THREAD










* If more sample volume is used in a tube than the specified working volume, it could be possible that deformation of the tube occurs. Request Test Report TR023201 for more information.

** Temperature range of the material of the tubes, caps and racks is approved from -80°C to +100°C. In practice tubes are used in lower temperatures until *vapor phase LN2* (-150°C to -196°C). Micronic does not recommend to store tubes in liquid phase nitrogen (LN2). In practice tubes are only snap frozen in LN2 (without submerging the cap).

*** For -80°C storage of the 2.00ml tube its recommended to always use a rack with cover.

TUBE OVERVIEW AND SPECIFICATIONS - EXTERNAL THREAD HYBRID

| | 0.75 ML | 1.40 ML | 2.00 ML | 3.00 ML | 1.50 ML | 3.00 ML | 6.00 ML |
|--|---|---|---|--|---|---|---|
| |  Page 28 |  Page 29 |  Page 30 |  Page 31 |  Page 32 |  Page 33 |  Page 34 |
| Ultra low temperature rack | Micronic 96-2 | Micronic 96-4 | Micronic 48-2 | Micronic 48-3 | Micronic 24-1 | Micronic 24-1 | Micronic 24-4 |
| Thread type | External | External | External | External | External | External | External |
| Tube type | Hybrid tube | Hybrid tube | Hybrid tube | Hybrid tube | Hybrid tube | Hybrid tube | Hybrid tube |
| Inner tube shape | U | U | Flat | Flat | Flat | Flat | Flat |
| Working volume with Push Cap (+21°C)* | 0.57ml | 1.03ml | n/a | n/a | n/a | n/a | n/a |
| Working volume with Screw Cap (+21°C)* | 0.62ml | 1.20ml | 1.95ml | 2.90ml | 1.45ml | 2.45ml | 6.00ml |
| Working volume with Screw Cap (-20°C)* | 0.62ml | 1.06ml | 1.80ml | 2.60ml | 1.40ml | 2.25ml | 5.55ml |
| Working volume with Screw Cap (-80°C)* | 0.60ml | 1.00ml | 1.65ml | 2.60ml | 1.30ml | 2.25ml | 5.30ml |
| 2D Data-Matrix coded (bottom) | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Alphanumeric coded (bottom) | No | No | No | No | No | No | No |
| Non-coded (bottom) | No | No | No | No | No | No | No |
| Human readable coded (bottom) | No | No | No | No | No | No | No |
| 1D barcoded (side) | Yes | Yes | Yes | Yes | No | No | Yes |
| Human readable coded (side) | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Capping options | Screw Cap Screw Cap IG Push Cap | Screw Cap Screw Cap IG Push Cap | Screw Cap | Screw Cap | Screw Cap | Screw Cap | Screw Cap |
| Outer diameter | 8.8mm | 8.8mm | 13.0mm | 13.0mm | 16.3mm | 16.3mm | 16.3mm |
| Tube height | 30.0mm | 48.0mm | 30.0mm | 41.0mm | 16.0mm | 23.0mm | 48.0mm |
| Tube height with Push Cap | 31.4mm | 49.4mm | n/a | n/a | n/a | n/a | n/a |
| Tube height with Screw Cap IG | 32.3mm | 50.3mm | n/a | n/a | n/a | n/a | n/a |
| Tube height with Screw Cap | 33.5mm | 51.5mm | 33.5mm | 44.5mm | 19.5mm | 26.5mm | 51.5mm |
| Available precapped | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Twist-lock | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Snap tube option | No | No | No | No | No | No | No |
| Sterilization option | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| DNase / RNase free | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Class 7 clean room production | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Minimum storage temperature** | Vapor phase LN2 | Vapor phase LN2 | Vapor phase LN2 | Vapor phase LN2 | Vapor phase LN2 | Vapor phase LN2 | Vapor phase LN2 |
| Availability | Bulk, Rack | Bulk, Rack | Bulk, Rack | Bulk, Rack | Bulk, Rack | Bulk, Rack | Bulk, Rack |

The 0.50ml Tubes with Internal Thread allow you to save between 35% and 50% of your freezer space and are ideal for low volume sample storage. The tube can be sealed with Push Caps and Screw Caps. Sealing 0.50ml Tubes with Low Profile Screw Caps provides an exceptionally efficient storage solution.









SUPERIOR FEATURES

- Internal thread provides the most optimal sealing for storage at ultra-low temperatures
- Available Non-coded, Alphanumeric coded or with a unique 2D Data-Matrix code (adhere to the ECC200 standard)
- Codes are permanently laser etched
- 2D codes also have a human-readable code
- Black code surface cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk, Lobarack-96, 96-1 and 96-Q1 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 2D codes
- Highly compact solution to save freezer space
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Snap tubes; tubes are locked into the rack wells to prevent sample loss from overturned racks
- Precapped with (Low Profile) Screw Cap or Push Cap
- Customized 2D Data-Matrix tube code and 1D rack barcode

SPECIFICATIONS

| | |
|---|--|
|  | Working volume at +21°C is 0.21 ml (Screw Cap) / 0.30ml (Push Cap) |
|  | V-bottom |
|  | Highest purity Polypropylene |
|  | Tube height is 18.5mm / Tube height in rack is 19.5mm (see page 51 for dim. drawing) |
|  | Resistant to Chemical solvents |
|  | Vapor phase LN2 (with Screw Cap) / -80°C (with Push Cap) to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 36.

0.75ML STORAGE TUBES INTERNAL THREAD

The popular 0.75ml Tubes with Internal Thread enable you to store your valuable samples in a safe and traceable way. The tubes are available in a Push Cap and Screw Cap version, and with a variety of codings. The alphanumeric and 2D codes are laser etched into the tube bottom to guarantee sample traceability.



96-WELL FORMAT

SUPERIOR FEATURES

- Internal thread provides the most optimal sealing for storage at ultra-low temperatures
- Available Non-coded, Alphanumeric coded or with a unique 2D Data-Matrix code (adhere to the ECC200 standard)
- Codes are permanently laser etched
- 2D codes also have a human-readable code
- Black code surface cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk, refill, Lobarack-96, 96-2 ULT Rack and 138-2 HD Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Amber tubes for light sensitive samples
- Sterilization using gamma irradiation or Eto treatment
- Snap tubes; tubes are locked into the rack wells to prevent sample loss from overturned racks
- Precapped with (Low Profile) Screw Cap or Push Cap
- Customized 2D Data-Matrix tube code and 1D rack barcode



SPECIFICATIONS

| | | |
|--|--|--|
| | | For the working volume, please see 'Tube Overview and Specs' on page 6-7 |
| | | V-bottom (Screw Cap tube) / V- or U-bottom (Push Cap tube) |
| | | Highest purity Polypropylene |
| | | Tube height is 26.0mm / Tube height in rack is 27.0mm (see page 51 for dim. drawing) |
| | | Resistant to Chemical solvents |
| | | Vapor phase LN2 (with Screw Cap) / -80°C (with Push Cap) to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 36.

1.10ML STORAGE TUBES INTERNAL THREAD

The traceable 1.10ml Tubes work equally well with either Screw Caps or Push Caps. The tubes are standard available in the Roborack-96 or Micronic 96-3 Rack. The 1D barcoded racks are based on the ANSI/SLAS standards for storage racks and are, just like the tubes, produced in a class 7 clean room.



96-WELL FORMAT

SUPERIOR FEATURES

- Internal thread provides the most optimal sealing for storage at ultra-low temperatures
- Available Non-coded, Alphanumeric coded or with a unique 2D Data-Matrix code (adhere to the ECC200 standard)
- Codes are permanently laser etched
- 2D codes also have a human-readable code
- Black code surface cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk, Roborack-96 and 96-3 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Snap tubes; tubes are locked into the rack wells to prevent sample loss from overturned racks
- Precapped with (Low Profile) Screw Cap or Push Cap
- Customized 2D Data-Matrix tube code and 1D rack barcode



Snap tubes

SPECIFICATIONS

| | | |
|--|--|--|
| | | Working volume at +21°C is 0.80ml (Screw Cap) / 0.90ml (Push Cap) |
| | | V-bottom |
| | | Highest purity Polypropylene |
| | | Tube height is 36.5mm / Tube height in rack is 37.5mm (see page 51 for dim. drawing) |
| | | Resistant to Chemical solvents |
| | | Vapor phase LN2 (with Screw Cap) / -80°C (with Push Cap) to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 37.

1.40ML STORAGE TUBES

INTERNAL THREAD

1.40ml Non-coded Tubes are proven to be an industry standard solution for many sample storage applications. The tubes provide an affordable way to improve your sample storage, handling and logistic processes. The 1.40ml tubes are also available with code surfaces that cannot be separated from the transparent tube due a unique 2K injection molding technique.



96-WELL FORMAT



SUPERIOR FEATURES

- Internal thread provides the most optimal sealing for storage at ultra-low temperatures
- Available Non-coded, Alphanumeric coded or with a unique 2D Data-Matrix code (adhere to the ECC200 standard)
- Codes are permanently laser etched
- 2D codes also have a human-readable code
- Black code surface cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk, refill, Comorack-96, Roborack-96 and 96-4 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Amber tubes for light sensitive samples
- Sterilization using gamma irradiation or Eto treatment
- Snap tubes; tubes are locked into the rack wells to prevent sample loss from overturned racks
- Precapped with (Low Profile) Screw Cap or Push Cap
- Customized 2D Data-Matrix tube code and 1D rack barcode



Suited for bead beating

SPECIFICATIONS

| | |
|--|--|
| | For the working volumes, please see 'Tube Overview and Specs' on page 6-7 |
| | V- or U-bottom (Screw Cap tube) / V- or U-bottom (Push Cap tube) |
| | Highest purity Polypropylene |
| | Tube height is 44.0mm / Tube height in rack is 45.0mm (see page 51 for dim. drawing) |
| | Resistant to Chemical solvents |
| | Vapor phase LN2 (with Screw Cap) / -80°C (with Push Cap) to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 37.

2.00ML STORAGE TUBES

INTERNAL THREAD

The 2.00ml Tubes combine high sample storage capacity and absolute sample traceability in an industry standard 96-well configuration. The storage tubes are provided with an alphanumeric or unique 2D Data-Matrix code laser encrypted on the tube bottom to enable easy identification and traceable storage of your samples.



96-WELL FORMAT

SUPERIOR FEATURES

- Internal thread provides the most optimal sealing for storage at low temperatures
- Available Alphanumeric coded or with a unique 2D Data-Matrix code (adhere to the ECC200 standard) on the tube bottom
- Codes are permanently laser etched
- 2D codes also have a human-readable code
- Black code surface cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk and 96-5 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Snap tubes; tubes are locked into the rack wells to prevent sample loss from overturned racks
- Precapped with (Low Profile) Screw Cap or Push Cap
- Customized 2D Data-Matrix tube code and 1D rack barcode



SPECIFICATIONS

| | |
|--|--|
| | Working volume at +21°C is 1.49ml (Screw Cap) / 1.58ml (Push Cap) |
| | U-bottom |
| | Highest purity Polypropylene |
| | Tube height is 59.0mm / Tube height in rack is 60.0mm (see page 51 for dim. drawing) |
| | Resistant to Chemical solvents |
| | -80°C* to +100°C |

* For -80°C storage of the 2.00ml tube it is recommended to always use a rack with cover

ORDERING INFORMATION

More information about product numbers and ordering, see page 38.

4.00ML STORAGE TUBES

INTERNAL THREAD

Our 4.00ml Tubes in 48-well format are ideal for storing or transporting larger volumes of liquid or solid biological samples. The tubes provide optimal sealing enabling high integrity storage at ultra-low temperatures. Additionally, the tubes are compatible with Micronic Screw Cap Recappers and automated systems.









SUPERIOR FEATURES

- Internal thread provides the most optimal sealing for storage at ultra-low temperatures
- Available Non-coded, Alphanumeric coded or with a unique 2D Data-Matrix code (adhere to the ECC200 standard)
- Codes are permanently laser etched
- 2D codes also have a large human-readable code
- Black code surface cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk and 48-4 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 2D codes
- Class 7 clean room production
- 48-format Screw Caps are available in 12 colors
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Precapped with Screw Cap
- Customized 2D Data-Matrix tube code and 1D rack barcode

SPECIFICATIONS

| | |
|---|--|
|  | Working volume at +21°C is 2.70ml (Screw Cap) |
|  | Flat bottom |
|  | Highest purity Polypropylene |
|  | Tube height is 44.0mm / Tube height in rack is 45.0mm (see page 55 for dim. drawing) |
|  | Resistant to Chemical solvents |
|  | Vapor phase LN2 to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 38.

6.00ML STORAGE TUBES

INTERNAL THREAD

6.00ml Sample Storage Tubes in 24-well format have a large working volume of 4.80ml at +21°C. These tubes are available Non-coded or with a 2D Data-Matrix code. The 2D Data-Matrix code permanently laser-etched on the tube bottom identifies the sample, its coordinates within a 24-tube storage rack, the particular rack and the location of the rack in your freezer.









SUPERIOR FEATURES

- Internal thread provides the most optimal sealing for storage at ultra-low temperatures
- Available Non-coded or with a unique 2D Data-Matrix code (adhere to the ECC200 standard) on the tube bottom
- 2D codes and 1D rack barcodes are permanently laser etched
- A large human-readable code next to the 2D code
- Black code surface cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk and 24-4 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 24-format Screw Caps are available in 12 colors
- 100% quality control on all 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Precapped with Screw Cap
- Customized 2D Data-Matrix tube code and 1D rack barcode

SPECIFICATIONS

| | |
|---|--|
|  | Working volume at +21°C is 4.80ml (Screw Cap) |
|  | Flat bottom |
|  | Highest purity Polypropylene |
|  | Tube height is 43.4mm / Tube height in rack is 44.0mm (see page 57 for dim. drawing) |
|  | Resistant to Chemical solvents |
|  | Vapor phase LN2 to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 38.

0.30ML STORAGE TUBES EXTERNAL THREAD

To address the need for low volume genomics, the 0.30ml Tube enables miniaturization of reaction volume required in genomic applications. The automation friendly tube cuts out intermediate (screening) plates and is more accessible for low volume liquid handlers. Ideal for DNA/RNA libraries, unique aliquoting and storage capacity with multi access.



96-WELL FORMAT

SUPERIOR FEATURES

- External thread for minimum cross contamination and maximum working volume
- Triple thread for optimal sealing quality
- Thick tube walls, sturdy design and excellent properties for ultra-low temperature storage
- Available with a unique 2D Data-Matrix code (adhere to the ECC standards) permanently laser etched on the tube bottom
- 2D codes also have a human-readable code
- Black code surface cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk and 96-1 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Precapped with Screw Cap (IG) or Push Cap
- Customized 2D Data-Matrix tube code and 1D rack barcode

Micronic 0.30ml Tubes External Thread

Christine Marchand
Status: Reviewer
Member since: 2019
Organization: Hoffmann-La Roche

Rating: 4.3 ★★★★★

Application Area: Storage of small aliquots of reference compounds
"A very convenient, space saving way of storage for small aliquots, storage in stackable 96er racks, identification made easy made by screw cap color coding"
Review date: 20 Jun 2019 (0.30ml Tubes External Thread)

Ease of use: ★★★★★
After sales service: ★★★★★
Value for money: ★★★★★

SPECIFICATIONS

| | |
|--|--|
| | For the working volumes, please see 'Tube Overview and Specs' on page 8-9 |
| | Vr- or V-bottom |
| | Highest purity Polypropylene |
| | Tube height is 16.4mm / Tube height in rack is 17.4mm (see page 51 for dim. drawing) |
| | Resistant to Chemical solvents |
| | Vapor phase LN2 to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 39.

0.40ML STORAGE TUBES EXTERNAL THREAD

The 0.40ml Tube enables efficient storage of low volumes of liquid samples. Due to thick tube walls, the tube has a sturdy design and excellent properties for ultra-low temperature storage. The 2D Data-Matrix codes are permanently laser-etched into the tube bottom so that they can never wear or fall off, ensuring absolute traceability.



96-WELL FORMAT

SUPERIOR FEATURES

- External thread for minimum cross contamination and maximum working volume
- Triple thread for optimal sealing quality
- Thick tube walls, sturdy design and excellent properties for ultra-low temperature storage
- Available with a unique 2D Data-Matrix code (adhere to the ECC standards) permanently laser etched on the tube bottom
- 2D codes also have a human-readable code
- Black code surface cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk and 96-1 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Precapped with Screw Cap (IG) or Push Cap
- Customized 2D Data-Matrix tube code and 1D rack barcode



SPECIFICATIONS

| | |
|--|--|
| | Working volume at +21°C is 0.40ml (Screw Cap) / 0.38ml (Push Cap) |
| | V-bottom |
| | Highest purity Polypropylene |
| | Tube height is 21.8mm / Tube height in rack is 22.8mm (see page 51 for dim. drawing) |
| | Resistant to Chemical solvents |
| | Vapor phase LN2 to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 39.

0.75ML STORAGE TUBES

EXTERNAL THREAD

Our externally threaded 0.75ml Tubes are provided with a triple start thread which gives an excellent closure to ensure long term sample preservation and storage. Using externally threaded storage tubes eliminates the possibility of the sample coming into contact with the screw thread, minimizing chance of cross-contamination while improving sample integrity.




96-WELL FORMAT

SUPERIOR FEATURES

- External thread for minimum cross contamination and maximum working volume
- Triple thread for optimal sealing quality
- Thick tube walls, sturdy design and excellent properties for ultra-low temperature storage
- Available with a unique 2D Data-Matrix code (adhere to the ECC standards) permanently laser etched on the tube bottom
- 2D codes also have a human-readable code
- Black code surface cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk, 96-2 ULT Rack and 138-2 HD Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS


- Sterilization using gamma irradiation or Eto treatment
- Precapped with Screw Cap (IG) or Push Cap
- Customized 2D Data-Matrix tube code and 1D rack barcode

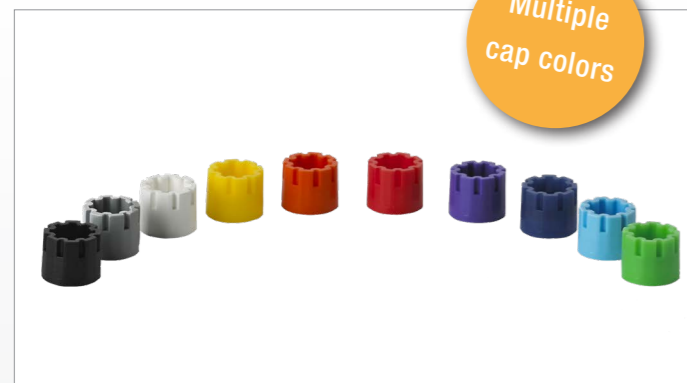

Multiple cap colors

SPECIFICATIONS

| | |
|--|--|
| | Working volume at +21°C is 0.60ml (Screw Cap) / 0.57ml (Push Cap) |
| | V-bottom |
| | Highest purity Polypropylene |
| | Tube height is 30.0mm / Tube height in rack is 31.0mm (see page 51 for dim. drawing) |
| | Resistant to Chemical solvents |
| | Vapor phase LN2 to +100°C |



ORDERING INFORMATION

More information about product numbers and ordering, see page 39.

0.80ML STORAGE TUBES

EXTERNAL THREAD

The space efficient 0.80ml Sample Tubes can hold a high (net) volume of biomaterial while occupying a relatively small amount of space, maximizing storage volume. Additionally, the code surface features a raised edge, which protects the 2D code against accidental scratches, ensuring absolute traceability.



96-WELL FORMAT

SUPERIOR FEATURES

- External thread for minimum cross contamination and maximum working volume
- Triple thread for optimal sealing quality
- Thick tube walls, sturdy design and excellent properties for ultra-low temperature storage
- Available with a unique 2D Data-Matrix code (adhere to the ECC standards) permanently laser etched on the tube bottom
- 2D codes also have a human-readable code
- Black code surface cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk, 96-3 ULT Rack and 138-2 HD Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free


ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Precapped with Screw Cap (IG) or Push Cap
- Customized 2D Data-Matrix tube code and 1D rack barcode



SPECIFICATIONS

| | |
|--|--|
| | Working volume at +21°C is 0.75ml (Screw Cap) / 0.68ml (Push Cap) |
| | V-bottom |
| | Highest purity Polypropylene |
| | Tube height is 34.8mm / Tube height in rack is 35.8mm (see page 51 for dim. drawing) |
| | Resistant to Chemical solvents |
| | Vapor phase LN2 to +100°C |

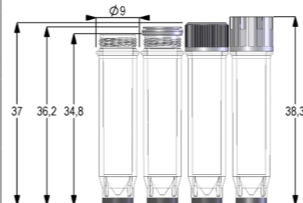

ORDERING INFORMATION

More information about product numbers and ordering, see page 39.

1.10ML STORAGE TUBES EXTERNAL THREAD

The triple thread of the 1.10ml Tubes External Thread gives an excellent closure and horizontal placement of the cap to ensure long-term sample preservation and storage. The code surface of the tubes with external thread features a raised edge, which protects the 2D code against accidental scratches, ensuring absolute traceability.



SUPERIOR FEATURES

- External thread for minimum cross contamination and maximum working volume
- Triple thread for optimal sealing quality
- Thick tube walls, sturdy design and excellent properties for ultra-low temperature storage
- Available with a unique 2D Data-Matrix code (adhere to the ECC standards) permanently laser etched on the tube bottom
- 2D codes also have a human-readable code
- Black code surface cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk and 96-3 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Precapped with Screw Cap (IG) or Push Cap
- Customized 2D Data-Matrix tube code and 1D rack barcode



SPECIFICATIONS

| | |
|--|--|
| | Working volume at +21°C is 0.93ml (Screw Cap) / 0.85ml (Push Cap) |
| | V-bottom |
| | Highest purity Polypropylene |
| | Tube height is 41.0mm / Tube height in rack is 42.3mm (see page 51 for dim. drawing) |
| | Resistant to Chemical solvents |
| | Vapor phase LN2 to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 40.

1.40ML STORAGE TUBES EXTERNAL THREAD

To anticipate on the trend to store samples in cryogenic temperatures, the 1.40ml Tubes with External Thread have a sturdy design and thick tube walls - excellent properties for ultra-low temperature storage. Using externally threaded storage tubes eliminates the possibility of the sample coming into contact with the screw thread, improving sample integrity.



SUPERIOR FEATURES

- External thread for minimum cross contamination and maximum working volume
- Triple thread for optimal sealing quality
- Thick tube walls, sturdy design and excellent properties for ultra-low temperature storage
- Available with a unique 2D Data-Matrix code (adhere to the ECC standards) permanently laser etched on the tube bottom
- 2D codes also have a human-readable code
- Black code surface cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk and 96-4 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Precapped with Screw Cap (IG) or Push Cap
- Customized 2D Data-Matrix tube code and 1D rack barcode



SPECIFICATIONS

| | |
|--|--|
| | Working volume at +21°C is 1.11ml (Screw Cap) / 1.03ml (Push Cap) |
| | V-bottom |
| | Highest purity Polypropylene |
| | Tube height is 48.0mm / Tube height in rack is 49.0mm (see page 51 for dim. drawing) |
| | Resistant to Chemical solvents |
| | Vapor phase LN2 to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 40.

2.00ML STORAGE TUBES

EXTERNAL THREAD

Our unmatched 2.00ml Tubes are ideal to store and handle high volume samples in an industry standard 96-well configuration. The triple thread of the tube with external thread gives an excellent closure and horizontal placement of the cap to ensure long-term sample preservation and storage.



96-WELL FORMAT

SUPERIOR FEATURES

- External thread for minimum cross contamination and maximum working volume
- Triple thread for optimal sealing quality
- Thick tube walls, sturdy design and excellent properties for low temperature storage
- Available with a unique 2D Data-Matrix code (adhere to the ECC standards) permanently laser etched on the tube bottom
- 2D codes also have a human-readable code
- Black code surface cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk and 96-5 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Precapped with Screw Cap (IG) or Push Cap
- Customized 2D Data-Matrix tube code and 1D rack barcode



SPECIFICATIONS

| | |
|--|--|
| | Working volume at +21°C is 1.55ml (Screw Cap) / 1.48ml (Push Cap) |
| | V-bottom |
| | Highest purity Polypropylene |
| | Tube height is 63.0mm / Tube height in rack is 63.8mm (see page 51 for dim. drawing) |
| | Resistant to Chemical solvents |
| | -80°C* to +100°C |

* For -80°C storage of the 2.00ml tube it is recommended to always use a rack with cover

ORDERING INFORMATION

More information about product numbers and ordering, see page 40.

1.00ML STORAGE TUBES

EXTERNAL THREAD

The Micronic 1.00ml Tissue Storage Tube with External Thread is specially designed for the storage of your valuable fresh frozen tissue samples. The 1.00ml tube can store solid samples up to a size of 0.40 x 0.40 x 0.70 cm. A wide shallow opening enables easy loading and removal of tissue samples using tweezers.



48-WELL FORMAT

SUPERIOR FEATURES

- External thread for minimum cross contamination and maximum working volume
- Triple thread for optimal sealing quality
- Thick tube walls, sturdy design and excellent properties for ultra-low temperature storage
- Available with a unique 2D Data-Matrix code (adhere to the ECC standards) permanently laser etched on the tube bottom
- 2D codes also have a human-readable code
- Transparent Screw Cap available for visual checks
- Black code surface cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk and 48-1 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Precapped with Screw Cap
- Customized 2D Data-Matrix tube code and 1D rack barcode



SPECIFICATIONS

| | |
|--|--|
| | Working volume at +21°C is 0.85ml (Screw Cap) |
| | Flat bottom |
| | Highest purity Polypropylene |
| | Tube height is 16.0mm / Tube height in rack is 17.0mm (see page 55 for dim. drawing) |
| | Resistant to Chemical solvents |
| | Vapor phase LN2 to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 40.

4.00ML STORAGE TUBES EXTERNAL THREAD

The 4.00ml Tube with External Thread has a massive working volume of 3.40ml (at +21°C). The sample storage tube is ideal for solid or large volumes of liquid samples. The 4.00ml tubes with a 2D Data-Matrix code permanently laser-etched on the bottom of the tube are the storage medium of choice for customers demanding long-term reliable code reading.



48-WELL FORMAT



SUPERIOR FEATURES

- External thread for minimum cross contamination and maximum working volume
- Triple thread for optimal sealing quality
- Thick tube walls, sturdy design and excellent properties for ultra-low temperature storage
- Available with a unique 2D Data-Matrix code (adhere to the ECC standards) permanently laser etched on the tube bottom
- 2D codes also have a human-readable code
- Transparent Screw Cap available for visual checks
- Black code surface cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk and 48-4 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free







ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Precapped with Screw Cap
- Customized 2D Data-Matrix tube code and 1D rack barcode



Ask about
code reading
options

SPECIFICATIONS

| | |
|---|--|
|  | Working volume at +21°C is 3.40ml (Screw Cap) |
|  | Flat bottom |
|  | Highest purity Polypropylene |
|  | Tube height is 48.0mm / Tube height in rack is 49.0mm (see page 55 for dim. drawing) |
|  | Resistant to Chemical solvents |
|  | Vapor phase LN2 to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 41.

3.50ML STORAGE TUBES EXTERNAL THREAD

The 3.50ml Tissue Tube with External Thread is specially designed for the storage of your fresh frozen tissue. The 3.50ml tube can store samples up to a size of 0.95 x 0.95 x 2.40 cm. With this tube, biobanks can take the next step to standardize and automate the storage of fresh frozen tissue samples.



24-WELL FORMAT

SUPERIOR FEATURES

- External thread for minimum cross contamination and maximum working volume
- Triple thread for optimal sealing quality
- Thick tube walls, sturdy design and excellent properties for ultra-low temperature storage
- Available with a unique 2D Data-Matrix code (adhere to the ECC standards) permanently laser etched on the tube bottom
- 2D codes also have a human-readable code
- Transparent Screw Cap available for visual checks
- Black code surface cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk and 24-2 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free







ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Precapped with Screw Cap
- Customized 2D Data-Matrix tube code and 1D rack barcode



Transparent
caps

SPECIFICATIONS

| | |
|---|--|
|  | Working volume at +21°C is 3.35ml (Screw Cap) |
|  | Flat bottom |
|  | Highest purity Polypropylene |
|  | Tube height is 30.0mm / Tube height in rack is 31.2mm (see page 57 for dim. drawing) |
|  | Resistant to Chemical solvents |
|  | Vapor phase LN2 to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 41.

0.75ML STORAGE TUBES

EXTERNAL THREAD HYBRID

The Tubes with External Thread Hybrid uniquely combine automated and visual sample identification. The unique combination of coding concepts on the 0.75ml Tube guarantees the complete traceability of samples. The high contrast 1D, 2D and human-readable codes are permanently laser-etched into the white surfaces, so that they can never wear or fall off.



SUPERIOR FEATURES

- 1D barcode/human readable code on the tube side wall for visual identification or verification
- Available with a 2D Data-Matrix code (adhere to the ECC standards) permanently laser etched on the tube bottom
- External thread for minimum cross contamination and maximum working volume
- Triple thread for optimal sealing quality
- Thick tube walls, sturdy design and excellent properties for ultra-low temperature storage
- White code surfaces cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk, 96-2 ULT Rack and 138-2 HD Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 1D and 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Precapped with Screw Cap or Push Cap
- Customized tube codes and rack barcode



SPECIFICATIONS

| | |
|--|--|
| | Working volume at +21°C is 0.62ml (Screw Cap) / 0.57ml (Push Cap) |
| | U-bottom |
| | Highest purity Polypropylene |
| | Tube height is 30.0mm / Tube height in rack is 31.0mm (see page 51 for dim. drawing) |
| | Resistant to Chemical solvents |
| | Vapor phase LN2 to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 41.

1.40ML STORAGE TUBES

EXTERNAL THREAD HYBRID

The transparent part of the tube walls enables easy visual checking of the sample, while the 4 coding concepts guarantee sample traceability. The 4 different coding concepts on the 1.40ml Tube enable researchers to verify information and keep identifying samples when automated or visual sample identification methods are challenging.




SUPERIOR FEATURES

- 1D barcode/human readable code on the tube side wall for visual identification or verification
- Available with a 2D Data-Matrix code (adhere to the ECC standards) permanently laser etched on the tube bottom
- External thread for minimum cross contamination and maximum working volume
- Triple thread for optimal sealing quality
- Thick tube walls, sturdy design and excellent properties for ultra-low temperature storage
- White code surfaces cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk and 96-4 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 1D and 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Precapped with Screw Cap or Push Cap
- Customized tube codes and rack barcode

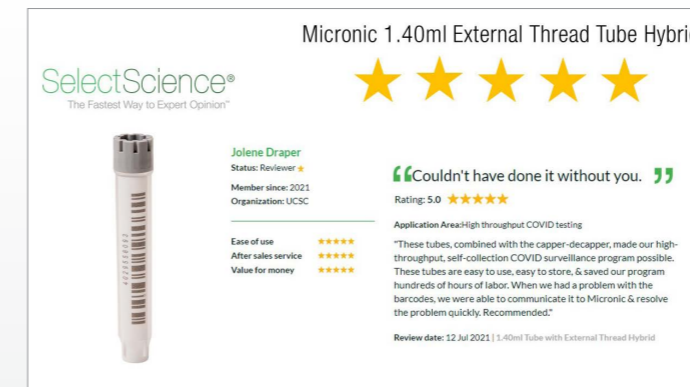

SPECIFICATIONS

| | |
|--|--|
| | Working volume at +21°C is 1.20ml (Screw Cap) / 1.03ml (Push Cap) |
| | U-bottom |
| | Highest purity Polypropylene |
| | Tube height is 48.0mm / Tube height in rack is 49.0mm (see page 51 for dim. drawing) |
| | Resistant to Chemical solvents |
| | Vapor phase LN2 to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 41 & 42.

2.00ML STORAGE TUBES

EXTERNAL THREAD HYBRID

The 2.00ml Tubes with External Thread Hybrid guarantee sample traceability and optimize working volume. The transparent part of the tube walls enable easy visual checking of the sample, while the 4 coding concepts guarantee sample traceability. Samples can be identified with the 2D Data-Matrix code, 1D barcode, numeric code and by written codes on the white tube surfaces.



SUPERIOR FEATURES

- 1D barcode/human readable code on the tube side wall for visual identification or verification
- Available with a 2D Data-Matrix code (adhere to the ECC standards) permanently laser etched on the tube bottom
- External thread for minimum cross contamination and maximum working volume
- Triple thread for optimal sealing quality
- Thick tube walls, sturdy design and excellent properties for ultra-low temperature storage
- White code surfaces cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk and 48-2 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 1D and 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Precapped with Screw Cap
- Customized tube codes and rack barcode



SPECIFICATIONS

| | | |
|--|--|--|
| | | Working volume at +21°C is 1.95ml (Screw Cap) |
| | | Flat bottom |
| | | Highest purity Polypropylene |
| | | Tube height is 30.0mm / Tube height in rack is 31.0mm (see page 55 for dim. drawing) |
| | | Resistant to Chemical solvents |
| | | Vapor phase LN2 to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 42.

3.00ML STORAGE TUBES

EXTERNAL THREAD HYBRID

The 3.00ml Tubes with External Thread Hybrid guarantee sample traceability, optimize working volume and increase overall sample storage efficiency. The unique combination of coding concepts and the dimensions of the tubes make the tubes extremely versatile. The tubes are compatible with ANSI / SLAS format racks, standard 9x9 or 9x10 cryo boxes, automated stores and liquid handlers.



SUPERIOR FEATURES

- 1D barcode/human readable code on the tube side wall for visual identification or verification
- Available with a 2D Data-Matrix code (adhere to the ECC standards) permanently laser etched on the tube bottom
- External thread for minimum cross contamination and maximum working volume
- Triple thread for optimal sealing quality
- Thick tube walls, sturdy design and excellent properties for ultra-low temperature storage
- White code surfaces cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk and 48-3 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 1D and 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free



ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Precapped with Screw Cap
- Customized tube codes and rack barcode

SPECIFICATIONS

| | | |
|--|--|--|
| | | Working volume at +21°C is 2.90ml (Screw Cap) |
| | | Flat bottom |
| | | Highest purity Polypropylene |
| | | Tube height is 41.0mm / Tube height in rack is 42.0mm (see page 55 for dim. drawing) |
| | | Resistant to Chemical solvents |
| | | Vapor phase LN2 to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 42.

1.50ML STORAGE TUBES

EXTERNAL THREAD HYBRID

Our 1.50ml Tube in 24-well format is a fully traceable solution to store your valuable biopsy, donor or transplant samples. The tube has an external screw thread, a wide opening and a flat inner bottom, enabling easy loading / removal of tissue samples using tweezers. The white code surface cannot be separated from the transparent tube due to the unique 2K injection molding technique.



24-WELL FORMAT

SUPERIOR FEATURES

- Numeric human readable code on the tube side wall for visual identification or verification
- Available with a 2D Data-Matrix code (adhere to the ECC standards) permanently laser etched on the tube bottom
- External thread for minimum cross contamination and maximum working volume
- Triple thread for optimal sealing quality
- Thick tube walls, sturdy design and excellent properties for ultra-low temperature storage
- White code surfaces cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk and 24-1 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 1D and 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Precapped with Screw Cap
- Customized tube codes and rack barcode



SPECIFICATIONS

| | |
|--|--|
| | Working volume at +21°C is 1.45ml (Screw Cap) |
| | Flat bottom |
| | Highest purity Polypropylene |
| | Tube height is 16.0mm / Tube height in rack is 17.0mm (see page 57 for dim. drawing) |
| | Resistant to Chemical solvents |
| | Vapor phase LN2 to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 42.

3.00ML STORAGE TUBES

EXTERNAL THREAD HYBRID

This tissue tube is ideal for the storage of solid tissue samples. With a unique 2D Data-Matrix code laser etched on the tube bottom, and a large human readable code on the tube sidewall, the 3.00ml Tissue Tube provides easy and unambiguous identification and storage of your samples.



24-WELL FORMAT

SUPERIOR FEATURES

- Numeric human readable code on the tube side wall for visual identification or verification
- Available with a 2D Data-Matrix code (adhere to the ECC standards) permanently laser etched on the tube bottom
- External thread for minimum cross contamination and maximum working volume
- Triple thread for optimal sealing quality
- Thick tube walls, sturdy design and excellent properties for ultra-low temperature storage
- White code surfaces cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk and 24-1 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all 1D and 2D codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Precapped with Screw Cap
- Customized tube codes and rack barcode



SPECIFICATIONS

| | |
|--|--|
| | Working volume at +21°C is 2.45ml (Screw Cap) |
| | Flat bottom |
| | Highest purity Polypropylene |
| | Tube height is 23.0mm / Tube height in rack is 24.0mm (see page 57 for dim. drawing) |
| | Resistant to Chemical solvents |
| | Vapor phase LN2 to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 43.

This 6.00ml Tube with External Thread Hybrid in 24-well format enables researchers to store tissue and high volumes of liquid samples. The combination of coding concepts and the dimensions of the tubes provide a secure solution to store samples at low temperatures while guaranteeing sample traceability.



SUPERIOR FEATURES

- 1D barcode/human readable code on the tube side wall for visual identification or verification
- Available with a 2D Data-Matrix code (adhere to the ECC standards) permanently laser etched on the tube bottom
- External thread for minimum cross contamination and maximum working volume
- Triple thread for optimal sealing quality
- Thick tube walls, sturdy design and excellent properties for ultra-low temperature storage
- White code surfaces cannot be separated from the transparent tube due to the 2K injection molding technique
- Available in bulk and 24-4 ULT Rack (based on the ANSI / SLAS standards for storage racks)
- 100% quality control on all codes
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Precapped with Screw Cap
- Customized tube codes and rack barcode


SPECIFICATIONS

| | | |
|--|--|--|
| | | Working volume at +21°C is 6.00ml |
| | | Flat bottom |
| | | Highest purity Polypropylene |
| | | Tube height is 48.0mm / Tube height in rack is 49.0mm (see page 57 for dim. drawing) |
| | | Resistant to Chemical solvents |
| | | Vapor phase LN2 to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 43.

We offer two special storage solutions. First, an economical, Non-coded solution for storing up to 8 samples in a linked row (1.40ml). The Tubestrip-8 have the same features and benefits as other Micronic tubes. Second, we offer amber tinted polypropylene storage tubes. Used in conjunction with a screw cap or push cap, the Amber Tubes ensure the integrity of light sensitive biological samples.

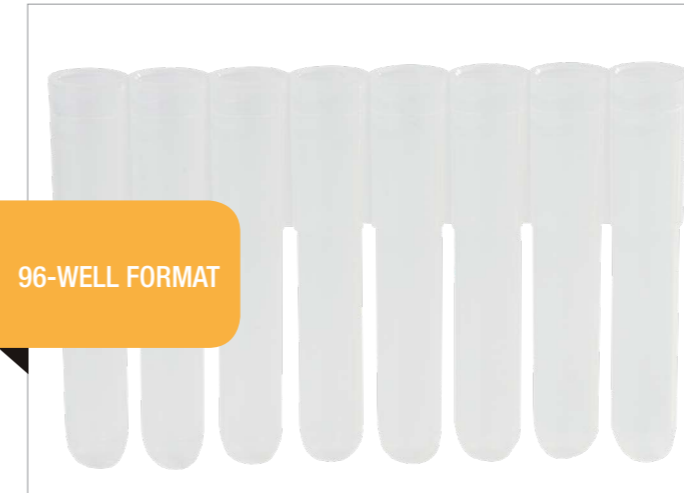

SUPERIOR FEATURES

Tubestrip-8

- Standard available Non-coded
- Class 7 clean room production
- RNase/DNase and Pyrogen free
- Additional options: sterilization using gamma irradiation or Eto treatment, snap tubestrips
- Low cost solution for storing up to 8 samples in a linked row

Amber tubes

- For light sensitive samples
- Available Non-coded or with unique laser etched 2D Data-Matrix code (adhere to the ECC200 standard) on the tube bottom
- 2D codes also have a human-readable code
- Black code surface cannot be separated from the transparent tube due to the unique 2K injection molding technique
- Available in Lobarack-96, Roborack-96 and ULT Rack
- Class 7 clean room production
- RNase/DNase and Pyrogen free
- Additional options: sterilization using gamma irradiation or Eto treatment, snap tubes, precapped, customized 2D Data-Matrix code and 1D rack barcode

SPECIFICATIONS

| | | |
|--|--|--|
| | | Working volume at +21°C is 1.20ml |
| | | U-bottom |
| | | Highest purity Polypropylene |
| | | Tube height is 45.4mm / Tube height with Push Cap rack is 46.4mm (EVA) |
| | | Resistant to Chemical solvents |
| | | -80°C to +100°C |

*For specifications of the Amber tubes see page 11 (0.75ml) or page 13 (1.40ml)

ORDERING INFORMATION

More information about product numbers and ordering, see page 43.

ORDERING INFORMATION - TUBES WITH INTERNAL THREAD

0.75ML STORAGE TUBES

| Catalog number | Description | Case Size |
|---|---|---------------|
| 0.50ml Non coded Screw Cap tubes with Internal Thread V-bottom without caps | | |
| MP32040 | 0.50ml Non coded Screw Cap tubes with Internal Thread in Bulk | 5 x 960 tubes |
| MPW32041BC3 | 0.50ml Non coded Screw Cap tubes with Internal Thread in Lobarack-96 (standard cover) | 4 x 10 racks |
| MPW32068BC3 | 0.50ml Non coded Screw Cap tubes with Internal Thread in Lobarack-96 (low cover) | 4 x 10 racks |
| MP32100 | 0.50ml Non coded Screw Cap tubes with Internal Thread in Micronic 96-1 Rack with low cover | 4 x 10 racks |
| MP32101 | 0.50ml Non coded Screw Cap tubes with Internal Thread in Micronic 96-Q1 Rack with high cover | 4 x 10 racks |
| 0.50ml Alphanumeric coded Screw Cap tubes with Internal Thread V-bottom without caps | | |
| MPW42049BC3 | 0.50ml Alphanumeric coded Screw Cap tubes with Internal Thread in Lobarack-96 (standard cover) | 4 x 10 racks |
| MPW42072BC3 | 0.50ml Alphanumeric coded Screw Cap tubes with Internal Thread in Lobarack-96 (low cover) | 4 x 10 racks |
| MP42100 | 0.50ml Alphanumeric coded Screw Cap tubes with Internal Thread in Micronic 96-1 Rack with low cover | 4 x 10 racks |
| MP42101 | 0.50ml Alphanumeric coded Screw Cap tubes with Internal Thread in Micronic 96-Q1 Rack with high cover | 4 x 10 racks |
| 0.50ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread V-bottom without caps | | |
| MP52326 | 0.50ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Bulk | 5 x 960 tubes |
| MPW52325BC3 | 0.50ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Lobarack-96 (standard cover) | 4 x 10 racks |
| MPW52337BC3 | 0.50ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Lobarack-96 (low cover) | 4 x 10 racks |
| MP52500 | 0.50ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Micronic 96-1 Rack with low cover | 4 x 10 racks |
| MP52501 | 0.50ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Micronic 96-Q1 Rack with high cover | 4 x 10 racks |

Also available in a Rack with Bottom ID or precapped with (Low Profile) Screw Cap or Push Cap.

0.75ML STORAGE TUBES

| Catalog number | Description | Case Size |
|---|--|----------------|
| 0.75ml Non coded Screw Cap tubes with Internal Thread V-bottom without caps | | |
| MP32069 | 0.75ml Non coded Screw Cap tubes with Internal Thread in Bulk | 5 x 960 tubes |
| MP32070 | 0.75ml Non coded Screw Cap tubes with Internal Thread in refill carrier-96 | 5 x 10 refills |
| MPW32071BC3 | 0.75ml Non coded Screw Cap tubes with Internal Thread in Lobarack-96 | 4 x 10 racks |
| MP32110 | 0.75ml Non coded Screw Cap tubes with Internal Thread in Micronic 96-2 Rack with low cover | 4 x 10 racks |
| MP32111 | 0.75ml Non coded Screw Cap tubes with Internal Thread in Micronic 96-2 Rack with high cover | 4 x 10 racks |
| 0.75ml Alphanumeric coded Screw Cap tubes with Internal Thread V-bottom without caps | | |
| MP42067 | 0.75ml Alphanumeric coded Screw Cap tubes with Internal Thread in refill carrier-96 | 5 x 10 refills |
| MPW42068BC3 | 0.75ml Alphanumeric coded Screw Cap tubes with Internal Thread in Lobarack-96 | 4 x 10 racks |
| MP42110 | 0.75ml Alphanumeric coded Screw Cap tubes with Internal Thread in Micronic 96-2 Rack with low cover | 4 x 10 racks |
| MP42111 | 0.75ml Alphanumeric coded Screw Cap tubes with Internal Thread in Micronic 96-2 Rack with high cover | 4 x 10 racks |
| 0.75ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread V-bottom without caps | | |
| MP52349 | 0.75ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Bulk | 5 x 960 tubes |
| MP52350 | 0.75ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in refill carrier-96 | 5 x 10 refills |
| MPW52351BC3 | 0.75ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Lobarack-96 | 4 x 10 racks |
| MP52519 | 0.75ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Micronic 96-2 Rack with low cover | 4 x 10 racks |
| MP52520 | 0.75ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Micronic 96-2 Rack with high cover | 4 x 10 racks |

Other item numbers are available for 0.75ml Push Cap tubes V- or U-bottom. Also available in a Rack with Bottom ID, 138-2 HD Rack or precapped with (Low Profile) Screw Cap or Push Cap.

1.10ML STORAGE TUBES

| Catalog number | Description | Case Size |
|---|--|---------------|
| 1.10ml Non Coded Screw Cap tubes with Internal Thread V-bottom without caps | | |
| MP32033 | 1.10ml Non coded Screw Cap tubes with Internal Thread in Bulk | 4 x 960 tubes |
| MP32033L | 1.10ml Non coded Screw Cap tubes with Internal Thread Snap in Bulk | 4 x 960 tubes |
| MPW32034BC3 | 1.10ml Non coded Screw Cap tubes with Internal Thread in Roborack-96 | 4 x 10 racks |
| MPW32034LBC3 | 1.10ml Non coded Screw Cap tubes with Internal Thread Snap in Roborack-96 | 4 x 10 racks |
| MP32120 | 1.10ml Non coded Screw Cap tubes with Internal Thread in Micronic 96-3 Rack with low cover | 4 x 10 racks |
| MP32121 | 1.10ml Non coded Screw Cap tubes with Internal Thread in Micronic 96-3 Rack with high cover | 4 x 10 racks |
| 1.10ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread V-bottom without caps | | |
| MP52318 | 1.10ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Bulk | 4 x 960 tubes |
| MP52318L | 1.10ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread Snap in Bulk | 4 x 960 tubes |
| MPW52319BC3 | 1.10ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Roborack-96 | 4 x 10 racks |
| MPW52319LBC3 | 1.10ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread Snap in Roborack-96 | 4 x 10 racks |
| MP52540 | 1.10ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Micronic 96-3 Rack with low cover | 4 x 10 racks |
| MP52541 | 1.10ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Micronic 96-3 Rack with high cover | 4 x 10 racks |

Also available in a Rack with Bottom ID or precapped with (Low Profile) Screw Cap or Push Cap.

1.40ML STORAGE TUBES

| Catalog number | Description | Case Size |
|--|--|----------------|
| 1.40ml Standard Push Cap tubes U-bottom without caps (round bottom) | | |
| MP32022 | 1.40ml Non coded Push Cap tubes in Bulk | 4 x 960 tubes |
| MPW42002BC3 | 1.40ml Non coded Push Cap tubes in Roborack-96 | 4 x 10 racks |
| MP22502 | 1.40ml Non coded Push Cap tubes in Comorack-96 (polycarbonate) | 4 x 10 racks |
| 1.40ml Screw Cap tubes with Internal Thread V-bottom without caps | | |
| MP32074 | 1.40ml Non coded Screw Cap tubes with Internal Thread in Bulk | 4 x 960 tubes |
| MPW32076BC3 | 1.40ml Non coded Screw Cap tubes with Internal Thread in Roborack-96 | 2 x 10 racks |
| MP32130 | 1.40ml Non coded Screw Cap tubes with Internal Thread in Micronic 96-4 Rack with low cover | 4 x 10 racks |
| MP32131 | 1.40ml Non coded Screw Cap tubes with Internal Thread in Micronic 96-4 Rack with high cover | 2 x 10 racks |
| MP42070 | 1.40ml Alphanumeric coded Screw Cap tubes with Internal Thread in refill carrier-96 | 5 x 10 refills |
| MPW42071BC3 | 1.40ml Alphanumeric coded Screw Cap tubes with Internal Thread in Roborack-96 | 2 x 10 racks |
| MP42130 | 1.40ml Alphanumeric coded Screw Cap tubes with Internal Thread in Micronic 96-4 Rack with low cover | 4 x 10 racks |
| MP42131 | 1.40ml Alphanumeric coded Screw Cap tubes with Internal Thread in Micronic 96-4 Rack with high cover | 2 x 10 racks |
| MP52353 | 1.40ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Bulk | 4 x 960 tubes |
| MPW52355BC3 | 1.40ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Roborack-96 | 2 x 10 racks |
| MP52550 | 1.40ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Micronic 96-4 Rack with low cover | 4 x 10 racks |
| MP52551 | 1.40ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Micronic 96-4 Rack with high cover | 2 x 10 racks |
| 1.40ml Screw Cap tubes with Internal Thread U-bottom without caps | | |
| MP32062 | 1.40ml Non coded Screw Cap tubes with Internal Thread in Bulk | 4 x 960 tubes |
| MPW32042BC3 | 1.40ml Non coded Screw Cap tubes with Internal Thread in Roborack-96 | 2 x 10 racks |
| MPW42063BC3 | 1.40ml Alphanumeric coded Screw Cap tubes with Internal Thread in Roborack-96 | 2 x 10 racks |
| MP52345 | 1.40ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Bulk | 4 x 960 tubes |
| MPW52330BC3 | 1.40ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Roborack-96 | 2 x 10 racks |

Other item numbers are available for 1.40ml Push Cap tubes V- or U-bottom. Also available in a Rack with Bottom ID precapped with (Low Profile) Screw Cap or Push Cap.

from page 13

from page 14

2.00ML STORAGE TUBES

| Catalog number | Description | Case Size |
|---|--|---------------|
| 2.00ml Screw Cap tubes with Internal Thread U-bottom without caps | | |
| MP42150 | 2.00ml Alphanumeric coded Screw Cap tubes with Internal Thread in Micronic 96-5 Rack with low cover | 4 x 6 racks |
| MP42151 | 2.00ml Alphanumeric coded Screw Cap tubes with Internal Thread in Micronic 96-5 Rack with high cover | 4 x 6 racks |
| MP52359 | 2.00ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Bulk | 4 x 576 tubes |
| MP52570 | 2.00ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Micronic 96-5 Rack with low cover | 4 x 6 racks |
| MP52571 | 2.00ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Micronic 96-5 Rack with high cover | 4 x 6 racks |
| 2.00ml Screw Cap tubes with Internal Thread U-bottom standard precapped with Grey (Low Profile) Screw Caps | | |
| MP42150-V20 | 2.00ml Alphanumeric coded Screw Cap tubes with Internal Thread precapped with Grey Low Profile Screw Caps in Micronic 96-5 Rack with low cover | 4 x 6 racks |
| MP42151-Z20 | 2.00ml Alphanumeric coded Screw Cap tubes with Internal Thread precapped with Grey Screw Caps in Micronic 96-5 Rack with high cover | 4 x 6 racks |
| MP52359-Z20 | 2.00ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread precapped with Grey Screw Caps in Bulk | 4 x 576 tubes |
| MP52570-V20 | 2.00ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread precapped with Grey Low Profile Screw Caps in Micronic 96-5 Rack with low cover | 4 x 6 racks |
| MP52571-Z20 | 2.00ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread precapped with Grey Screw Caps in Micronic 96-5 Rack with high cover | 4 x 6 racks |
| 2.00ml Screw Cap tubes with Internal Thread U-bottom standard precapped with TPE Caps Natural | | |
| MP52359-X01 | 2.00ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread precapped with TPE Caps Natural in Bulk | 4 x 576 tubes |
| MP52570-X01 | 2.00ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread precapped with TPE Caps Natural in Micronic 96-5 Rack with low cover | 4 x 6 racks |

from page 15

4.00ML STORAGE TUBES

| Catalog number | Description | Case Size |
|--|---|---------------|
| 4.00ml Screw Cap tubes with Internal Thread without caps | | |
| MP32200 | 4.00ml Non coded Screw Cap tubes with Internal Thread Bulk | 4 x 480 tubes |
| MP32201 | 4.00ml Non coded Screw Cap tubes with Internal Thread in Micronic 48-4 Rack with high cover | 2 x 10 racks |
| MP42200 | 4.00ml Alphanumeric coded Screw Cap tubes with Internal Thread in Micronic 48-4 Rack with high cover | 2 x 10 racks |
| MP52440 | 4.00ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Bulk | 4 x 480 tubes |
| MP52441 | 4.00ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Micronic 48-4 Rack with high cover | 2 x 10 racks |
| 4.00ml Screw Cap tubes with Internal Thread standard precapped with Grey Screw Caps | | |
| MP52440-Z20 | 4.00ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread precapped with Grey Screw Caps in Bulk | 4 x 480 tubes |
| MP52441-Z20 | 4.00ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread precapped with Grey Screw Caps in Micronic 48-4 Rack | 2 x 10 racks |

from page 16

6.00ML STORAGE TUBES

| Catalog number | Description | Case Size |
|--|---|---------------|
| 6.00ml Screw Cap tubes with Internal Thread without caps | | |
| MP32300 | 6.00ml Non coded Screw Cap tubes with Internal Thread Bulk | 4 x 240 tubes |
| MP32301 | 6.00ml Non coded Screw Cap tubes with Internal Thread in Micronic 24-4 Rack with high cover | 2 x 10 racks |
| MP52400 | 6.00ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Bulk | 4 x 240 tubes |
| MP52410 | 6.00ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread in Micronic 24-4 Rack with high cover | 2 x 10 racks |
| 6.00ml Screw Cap tubes with Internal Thread standard precapped with Grey Screw Caps | | |
| MP52400-Z20 | 6.00ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread precapped with Grey Screw Caps in Bulk | 4 x 240 tubes |
| MP52410-Z20 | 6.00ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread precapped with Grey Screw Caps in Micronic 24-4 Rack | 2 x 10 racks |

from page 17

ORDERING INFORMATION - TUBES WITH EXTERNAL THREAD

0.30ML STORAGE TUBES

| Catalog number | Description | Case Size |
|--|--|---------------|
| 0.30ml Non Coded Screw Cap tubes with External Thread without caps | | |
| MP32146 | 0.30ml Non coded Screw Cap tubes with External Thread V-bottom Bulk | 5 x 960 tubes |
| MP32147 | 0.30ml Non coded Screw Cap tubes with External Thread V-bottom in Micronic 96-1 Rack with low cover | 4 x 10 racks |
| 0.30ml 2D Data-Matrix coded Screw Cap tubes with External Thread without caps | | |
| MP52701 | 0.30ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom Bulk | 5 x 960 tubes |
| MP52703 | 0.30ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom in Micronic 96-1 Rack with low cover | 4 x 10 racks |

from page 18

Also available in a Rack with Bottom ID or precapped with Screw Cap (IG) or Push Cap

0.40ML STORAGE TUBES

| Catalog number | Description | Case Size |
|--|--|---------------|
| 0.40ml Screw Cap tubes with External Thread without caps | | |
| MP52716 | 0.40ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom Bulk | 5 x 960 tubes |
| MP52718 | 0.40ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom in Micronic 96-1 Rack with high cover | 4 x 10 racks |
| 0.40ml Screw Cap tubes with External Thread standard precapped with Grey Screw Cap IG | | |
| MP52716-G71 | 0.40ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom precapped with Grey Screw Cap IG Bulk | 5 x 960 tubes |
| MP52718-G71 | 0.40ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom precapped with Grey Screw Cap IG in Micronic 96-1 Rack with high cover | 4 x 10 racks |

from page 19

Also available in a Rack with Bottom ID or precapped with Screw Cap or Push Cap

0.75ML STORAGE TUBES

| Catalog number | Description | Case Size |
|--|---|---------------|
| 0.75ml Non Coded Screw Cap tubes with External Thread without caps | | |
| MP32150 | 0.75ml Non coded Screw Cap tubes with External Thread V-bottom Bulk | 5 x 960 tubes |
| MP32151 | 0.75ml Non coded Screw Cap tubes with External Thread V-bottom in Micronic 96-2 Rack with high cover | 4 x 10 racks |
| 0.75ml 2D Data-Matrix coded Screw Cap tubes with External Thread without caps | | |
| MP52704 | 0.75ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom Bulk | 4 x 960 tubes |
| MP52706 | 0.75ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom in Micronic 96-2 Rack with high cover | 4 x 10 racks |

from page 20

Also available in a Rack with Bottom ID or 138-2 HD Rack or precapped with Screw Cap (IG) or Push Cap

0.80ML STORAGE TUBES

| Catalog number | Description | Case Size |
|--|---|---------------|
| 0.80ml Screw Cap tubes with External Thread without caps | | |
| MP52720 | 0.80ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom Bulk | 4 x 960 tubes |
| MP52722 | 0.80ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom in Micronic 96-3 Rack with low cover | 4 x 10 racks |
| 0.80ml Screw Cap tubes with External Thread standard precapped with Grey Screw Cap IG | | |
| MP52720-G71 | 0.80ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom precapped with Grey Screw Cap IG Bulk | 4 x 960 tubes |
| MP52722-G71 | 0.80ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom precapped with Grey Screw Cap IG in Micronic 96-3 Rack with low cover | 4 x 10 racks |

from page 21

Also available in a Rack with Bottom ID or 138-2 HD Rack or precapped with Screw Cap or Push Cap

1.10ML STORAGE TUBES

| Catalog number | Description | Case Size |
|---|--|---------------|
| 1.10ml 2D Data-Matrix coded Screw Cap tubes with External Thread without caps | | |
| MP52730 | 1.10ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom Bulk | 4 x 960 tubes |
| MP52732 | 1.10ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom in Micronic 96-3 Rack with high cover | 4 x 10 racks |
| 1.10ml 2D Data-Matrix coded Screw Cap tubes with External Thread standard precapped with Grey Screw Cap IG | | |
| MP52730-G71 | 1.10ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom precapped with Grey Screw Cap IG Bulk | 4 x 960 tubes |
| MP52732-G71 | 1.10ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom precapped with Grey Screw Cap IG in Micronic 96-3 Rack with high cover | 4 x 10 racks |

Also available in a Rack with Bottom ID or precapped with Screw Cap or Push Cap

1.40ML STORAGE TUBES

| Catalog number | Description | Case Size |
|--|---|---------------|
| 1.40ml Non Coded Screw Cap tubes with External Thread without caps | | |
| MP32152 | 1.40ml Non coded Screw Cap tubes with External Thread V-bottom Bulk | 4 x 960 tubes |
| MP32153 | 1.40ml Non coded Screw Cap tubes with External Thread V-bottom in Micronic 96-4 Rack with high cover | 2 x 10 racks |
| 1.40ml 2D Data-Matrix coded Screw Cap tubes with External Thread without caps | | |
| MP52708 | 1.40ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom Bulk | 4 x 960 tubes |
| MP52710 | 1.40ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom in Micronic 96-4 Rack with high cover | 2 x 10 racks |

Also available in a Rack with Bottom ID or precapped with Screw Cap (IG) or Push Cap

2.00ML STORAGE TUBES

| Catalog number | Description | Case Size |
|--|--|---------------|
| 2.00ml Screw Cap tubes with External Thread without caps | | |
| MP52712 | 2.00ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom Bulk | 4 x 576 tubes |
| MP52714 | 2.00ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom in Micronic 96-5 Rack with high cover | 4 x 6 racks |
| 2.00ml Screw Cap tubes with External Thread standard precapped with Grey Screw Cap IG | | |
| MP52712-G71 | 2.00ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom precapped with Grey Screw Cap IG Bulk | 4 x 576 tubes |
| MP52714-G71 | 2.00ml 2D Data-Matrix coded Screw Cap tubes with External Thread V-bottom precapped with Grey Screw Cap IG in Micronic 96-5 Rack with high cover | 4 x 6 racks |

1.00ML STORAGE TUBES

| Catalog number | Description | Case Size |
|--|---|---------------|
| 1.00ml Screw Cap tubes with External Thread without caps | | |
| MP52800 | 1.00ml 2D Data-Matrix coded Screw Cap tubes with External Thread Flat bottom Bulk | 5 x 480 tubes |
| MP52802 | 1.00ml 2D Data-Matrix coded Screw Cap tubes with External Thread Flat bottom in Micronic 48-1 Rack with low cover | 4 x 10 racks |
| 1.00ml Screw Cap tubes with External Thread standard precapped with Grey Screw Caps | | |
| MP52800-Y20 | 1.00ml 2D Data-Matrix coded Screw Cap tubes with External Thread Flat bottom precapped with Grey Screw Cap Bulk | 5 x 480 tubes |
| MP52802-Y20 | 1.00ml 2D Data-Matrix coded Screw Cap tubes with External Thread Flat bottom precapped with Grey Screw Cap in Micronic 48-1 Rack with low cover | 4 x 10 racks |

Also available precapped with Screw Cap or Push Cap

4.00ML STORAGE TUBES

| Catalog number | Description | Case Size |
|---|--|---------------|
| 4.00ml Screw Cap tubes with External Thread without caps | | |
| MP52808 | 4.00ml 2D Data-Matrix coded Screw Cap tubes with External Thread Flat bottom Bulk | 4 x 480 tubes |
| MP52810 | 4.00ml 2D Data-Matrix coded Screw Cap tubes with External Thread Flat bottom in Micronic 48-4 Rack with high cover | 2 x 10 racks |
| 4.00ml Screw Cap tubes with External Thread precapped with Grey Screw Caps | | |
| MP52808-Y20 | 4.00ml 2D Data-Matrix coded Screw Cap tubes with External Thread Flat bottom precapped with Grey Screw Cap Bulk | 4 x 480 tubes |
| MP52810-Y20 | 4.00ml 2D Data-Matrix coded Screw Cap tubes with External Thread Flat bottom precapped with Grey Screw Cap in Micronic 48-4 Rack with high cover | 2 x 10 racks |

3.50ML STORAGE TUBES

| Catalog number | Description | Case Size |
|---|--|---------------|
| 3.50ml Screw Cap tubes with External Thread without caps | | |
| MP52904 | 3.50ml 2D Data-Matrix coded Screw Cap tubes with External Thread Flat bottom Bulk | 4 x 240 tubes |
| MP52906 | 3.50ml 2D Data-Matrix coded Screw Cap tubes with External Thread Flat bottom in Micronic 24-2 Rack with high cover | 4 x 10 racks |
| 3.50ml Screw Cap tubes with External Thread precapped with Grey Screw Caps | | |
| MP52904-Y20 | 3.50ml 2D Data-Matrix coded Screw Cap tubes with External Thread Flat bottom precapped with Grey Screw Cap Bulk | 4 x 240 tubes |
| MP52906-Y20 | 3.50ml 2D Data-Matrix coded Screw Cap tubes with External Thread Flat bottom precapped with Grey Screw Cap in Micronic 24-2 Rack with high cover | 4 x 10 racks |

ORDERING INFORMATION - TUBES WITH EXTERNAL THREAD HYBRID

0.75ML STORAGE TUBES HYBRID

| Catalog number | Description | Case Size |
|--|---|---------------|
| 0.75ml Screw Cap tubes with External Thread Hybrid without caps | | |
| MP52755 | 0.75ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D U-bottom Bulk | 4 x 960 tubes |
| MP52762 | 0.75ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W00 U-bottom Bulk | 4 x 960 tubes |
| MP52757 | 0.75ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D U-bottom in Micronic 96-2 Rack with high cover | 4 x 10 racks |
| MP52763 | 0.75ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W00 U-bottom in Micronic 96-2 Rack with high cover | 4 x 10 racks |
| 0.75ml Screw Cap tubes with External Thread Hybrid precapped with Grey Screw Cap IG | | |
| MP52755-G71 | 0.75ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D U-bottom precapped with Grey Screw Cap IG Bulk | 4 x 960 tubes |
| MP52757-G71 | 0.75ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D U-bottom precapped with Grey Screw Cap IG in Micronic 96-2 Rack with high cover | 4 x 10 racks |

Also available in a Rack with Bottom ID or a 138-2 HD Rack or precapped with Screw Cap or Push Cap

1.40ML STORAGE TUBES HYBRID

| Catalog number | Description | Case Size |
|--|--|---------------|
| 1.40ml Screw Cap tubes with External Thread Hybrid without caps | | |
| MP52759 | 1.40ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D U-bottom Bulk | 4 x 960 tubes |
| MP52764 | 1.40ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W00 U-bottom Bulk | 4 x 960 tubes |
| MP52761 | 1.40ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D U-bottom in Micronic 96-4 Rack with high cover | 2 x 10 racks |
| MP52765 | 1.40ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W00 U-bottom in Micronic 96-4 Rack with high cover | 2 x 10 racks |

1.40ML STORAGE TUBES HYBRID

from page 29

| Catalog number | Description | Case Size |
|--|---|---------------|
| 1.40ml Screw Cap tubes with External Thread Hybrid precapped with Grey Screw Cap IG | | |
| MP52759-G71 | 1.40ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D U-bottom precapped with Grey Screw Cap IG Bulk | 4 x 960 tubes |
| MP52761-G71 | 1.40ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D U-bottom precapped with Grey Screw Cap IG in Micronic 96-4 Rack with high cover | 2 x 10 racks |

Also available in a Rack with Bottom ID or precapped with Screw Cap or Push Cap

2.00ML STORAGE TUBES HYBRID

from page 30

| Catalog number | Description | Case Size |
|--|---|---------------|
| 2.00ml Screw Cap tubes with External Thread Hybrid without caps | | |
| MP52855 | 2.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D Flat bottom Bulk | 4 x 480 tubes |
| MP52856 | 2.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W00 Flat bottom Bulk | 4 x 480 tubes |
| MP52858 | 2.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D Flat bottom in Micronic 48-2 Rack with high cover | 4 x 10 racks |
| MP52859 | 2.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W00 Flat bottom in Micronic 48-2 Rack with high cover | 4 x 10 racks |
| 2.00ml Screw Cap tubes with External Thread Hybrid precapped with Grey Screw Caps | | |
| MP52855-Y20 | 2.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D Flat bottom precapped with Grey Screw Cap Bulk | 4 x 480 tubes |
| MP52858-Y20 | 2.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D Flat bottom precapped with Grey Screw Cap in Micronic 48-2 Rack with high cover | 4 x 10 racks |

3.00ML STORAGE TUBES HYBRID

from page 31

| Catalog number | Description | Case Size |
|--|---|---------------|
| 3.00ml Screw Cap tubes with External Thread Hybrid without caps | | |
| MP52861 | 3.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D Flat bottom Bulk | 4 x 480 tubes |
| MP52862 | 3.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W00 Flat bottom Bulk | 4 x 480 tubes |
| MP52864 | 3.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D Flat bottom in Micronic 48-3 Rack with high cover | 4 x 10 racks |
| MP52865 | 3.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W00 Flat bottom in Micronic 48-3 Rack with high cover | 4 x 10 racks |
| 3.00ml Screw Cap tubes with External Thread Hybrid precapped with Grey Screw Caps | | |
| MP52861-Y20 | 3.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D Flat bottom precapped with Grey Screw Cap Bulk | 4 x 480 tubes |
| MP52864-Y20 | 3.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D Flat bottom precapped with Grey Screw Cap in Micronic 48-3 Rack with high cover | 4 x 10 racks |

1.50ML STORAGE TUBES HYBRID

from page 32

| Catalog number | Description | Case Size |
|--|--|---------------|
| 1.50ml Screw Cap tubes with External Thread Hybrid without caps | | |
| MP52951 | 1.50ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid WHR Flat bottom Bulk | 5 x 240 tubes |
| MP52952 | 1.50ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W00 Flat bottom Bulk | 5 x 240 tubes |
| MP52954 | 1.50ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid WHR Flat bottom in Micronic 24-1 Rack with low cover | 4 x 10 racks |
| MP52955 | 1.50ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W00 Flat bottom in Micronic 24-1 Rack with low cover | 4 x 10 racks |
| 1.50ml Screw Cap tubes with External Thread Hybrid precapped with Grey Screw Caps | | |
| MP52951-Y20 | 1.50ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid WHR Flat bottom precapped with Grey Screw Cap Bulk | 5 x 240 tubes |
| MP52954-Y20 | 1.50ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid WHR Flat bottom precapped with Grey Screw Cap in Micronic 24-1 Rack with low cover | 4 x 10 racks |

3.00ML STORAGE TUBES HYBRID

from page 33

| Catalog number | Description | Case Size |
|--|---|---------------|
| 3.00ml Screw Cap tubes with External Thread Hybrid without caps | | |
| MP52957 | 3.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid WHR Flat bottom Bulk | 4 x 240 tubes |
| MP52958 | 3.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W00 Flat bottom Bulk | 4 x 240 tubes |
| MP52960 | 3.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid WHR Flat bottom in Micronic 24-1 Rack with high cover | 4 x 10 racks |
| MP52961 | 3.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W00 Flat bottom in Micronic 24-1 Rack with high cover | 4 x 10 racks |
| 3.00ml Screw Cap tubes with External Thread Hybrid precapped with Grey Screw Caps | | |
| MP52957-Y20 | 3.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid WHR Flat bottom precapped with Grey Screw Cap Bulk | 4 x 240 tubes |
| MP52960-Y20 | 3.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid WHR Flat bottom precapped with Grey Screw Cap in Micronic 24-1 Rack with high cover | 4 x 10 racks |

6.00ML STORAGE TUBES HYBRID

from page 34

| Catalog number | Description | Case Size |
|--|---|---------------|
| 6.00ml Screw Cap tubes with External Thread Hybrid without caps | | |
| MP52963 | 6.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D Flat bottom Bulk | 4 x 240 tubes |
| MP52964 | 6.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W00 Flat bottom Bulk | 4 x 240 tubes |
| MP52966 | 6.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D Flat bottom in Micronic 24-4 Rack with high cover | 2 x 10 racks |
| MP52967 | 6.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W00 Flat bottom in Micronic 24-4 Rack with high cover | 2 x 10 racks |
| 6.00ml Screw Cap tubes with External Thread Hybrid precapped with Grey Screw Caps | | |
| MP52963-Y20 | 6.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D Flat bottom precapped with Grey Screw Cap Bulk | 4 x 240 tubes |
| MP52966-Y20 | 6.00ml 2D Data-Matrix coded Screw Cap tubes with Ext. Thread Hybrid W1D Flat bottom precapped with Grey Screw Cap in Micronic 24-4 Rack with high cover | 2 x 10 racks |

ORDERING INFORMATION - SPECIAL TUBES

TUBESTRIP-8 & AMBER TUBES FOR LIGHT SENSITIVE SAMPLES

from page 35


| Catalog number | Description | Case Size |
|---|---|----------------|
| 1.40ml Non coded Push Cap Tubestrip U-bottom | | |
| MP32067 | 1.40ml Non coded Push Cap tubes in strip of 8-tubes | 4 x 120 strips |
| MP32067L | 1.40ml Non coded Push Cap tubes in strip of 8-tubes Snap | 4 x 120 strips |
| Products for Light Sensitive Samples - 0.75ml / V-bottom | | |
| MP32087 | 0.75ml Non coded Screw Cap tubes with Internal Thread Amber V-bottom Bulk | 5 x 960 tubes |
| MPW32078BC3 | 0.75ml Non coded Push Cap tubes Amber V-bottom in Lobarack-96 White | 4 x 10 racks |
| MPW32079BC3 | 0.75ml Non coded Screw Cap tubes with Internal Thread Amber V-bottom in Lobarack-96 White | 4 x 10 racks |
| MP32091 | 0.75ml Non coded Screw Cap tubes with Internal Thread Amber V-bottom in Micronic 96-2 Rack with high cover | 4 x 10 racks |
| MP32090 | 0.75ml Non-coded Push Cap tubes Amber V-bottom in Micronic 96-2 Rack with low cover | 4 x 10 racks |
| MPW52648BC3 | 0.75ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread Amber V-bottom in Lobarack-96 White | 4 x 10 racks |
| Products for Light Sensitive Samples - 1.40ml / U- or V-bottom | | |
| MP32088 | 1.40ml Non coded Screw Cap tubes with Internal Thread Amber V-bottom Bulk | 4 x 960 tubes |
| MPW32086BC3 | 1.40ml Non coded Screw Cap tubes with Internal Thread Amber V-bottom in Roborack-96 | 2 x 10 rack |
| MPW52649BC3 | 1.40ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread Amber V-bottom in Roborack-96 | 2 x 10 racks |
| MP52560 | 1.40ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread Amber V-bottom in Micronic 96-4 Rack with low cover | 4 x 10 racks |
| MP52561 | 1.40ml 2D Data-Matrix coded Screw Cap tubes with Internal Thread Amber V-bottom in Micronic 96-4 Rack with high cover | 2 x 10 racks |



SERVING SCIENTISTS
SINCE 1984


CERTIFIED CLASS 7 CLEAN ROOM
MANUFACTURING

SUPERIOR FEATURES OF MICRONIC SAMPLE STORAGE RACKS


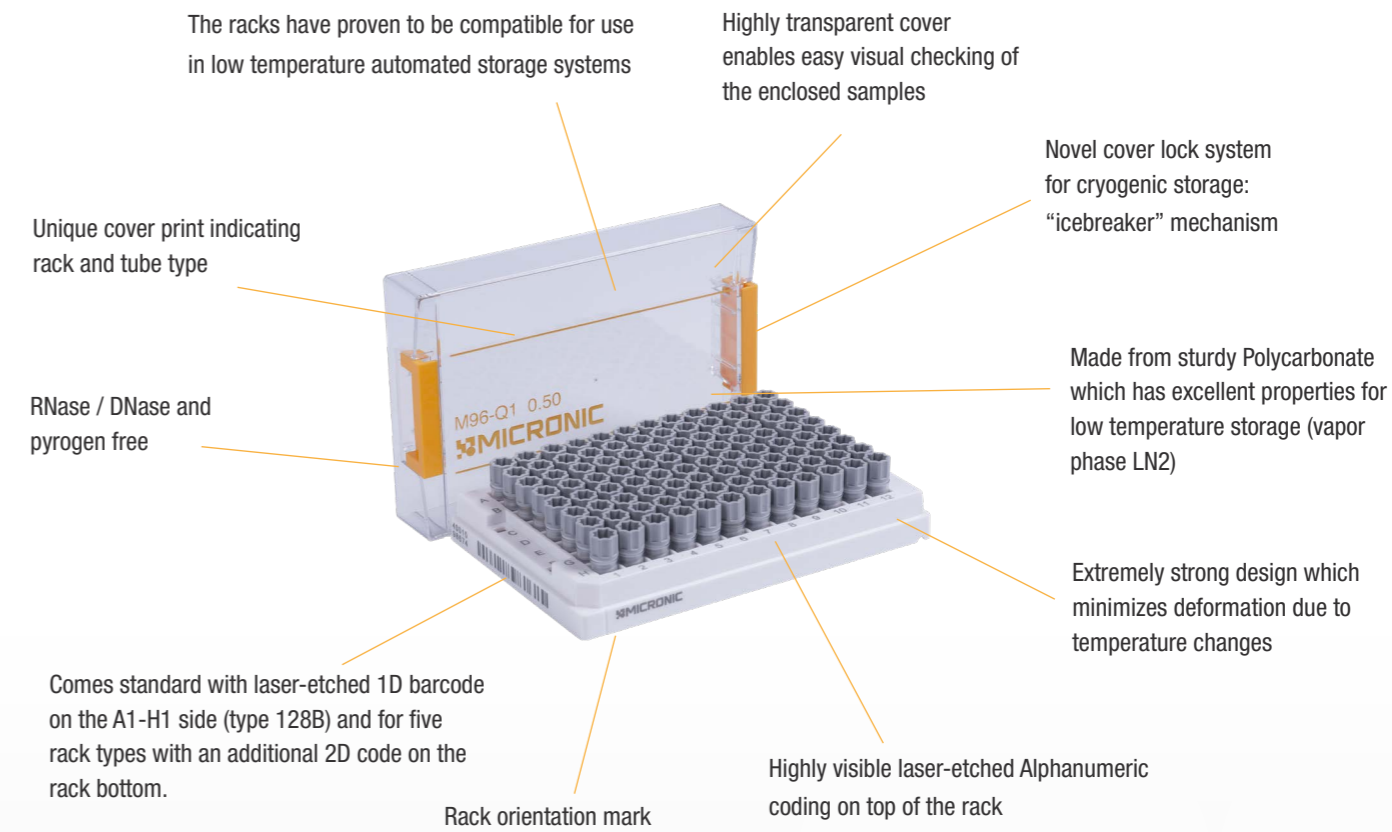






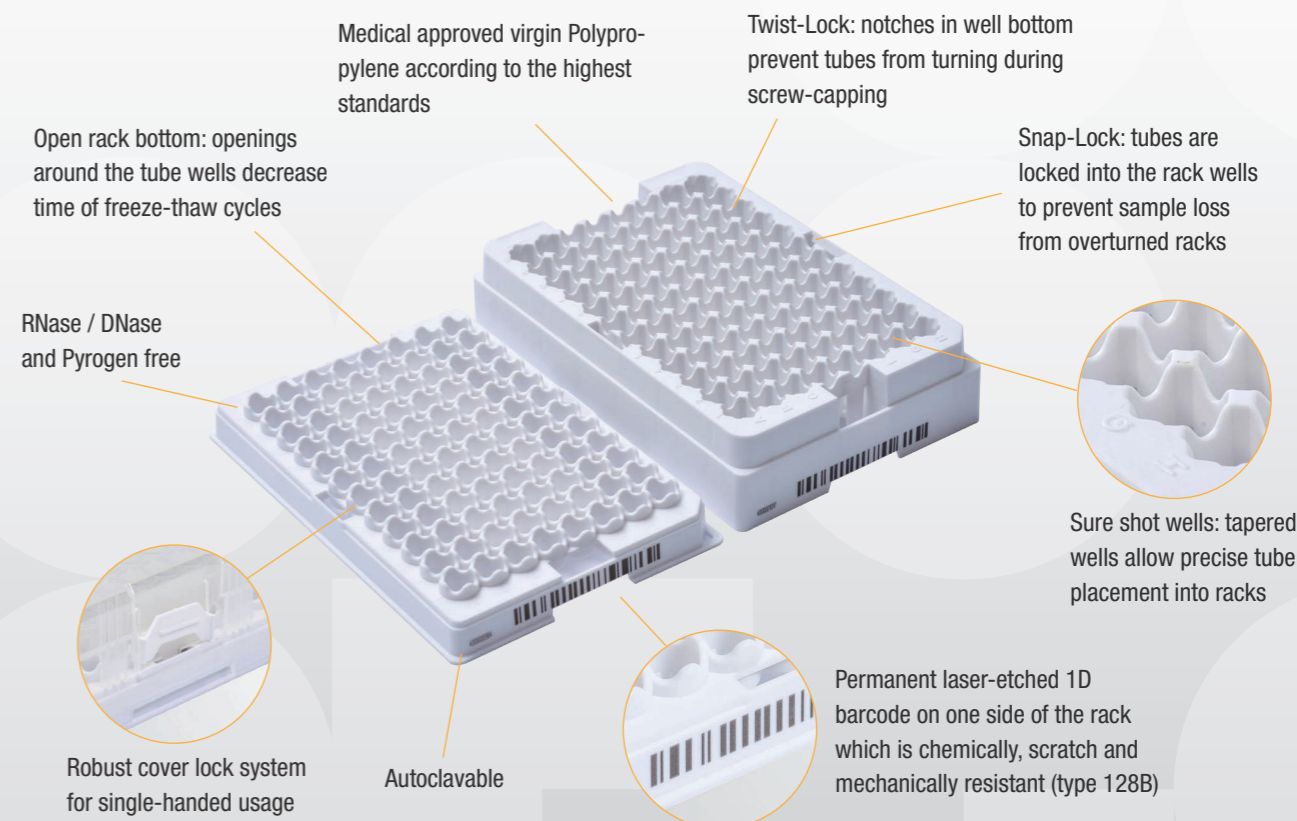

SAMPLE STORAGE RACKS

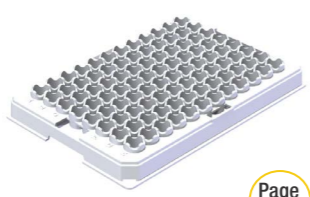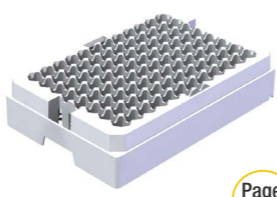ALL MICRONIC STORAGE RACKS ARE BASED ON THE GLOBAL RECOGNIZED ANSI / SLAS 1-2004 THROUGH ANSI / SLAS 4-2004 STANDARDS (FORMERLY RECOGNIZED AS THE ANSI / SBS STANDARDS) AND PROVIDE AN EASY AND AUTOMATION FRIENDLY METHOD TO STORE YOUR SAMPLE STORAGE TUBES. FOR EACH SAMPLE STORAGE TUBE FORMAT MICRONIC OFFERS A ROBUST AND USER FRIENDLY STORAGE RACK. THE RACKS FEATURE A LASER ETCHED 1D RACK BARCODE ON ONE SIDE OF THE RACK AND FOR FIVE RACKS AN ADDITIONAL 2D DATA-MATRIX CODE ON THE BOTTOM. FURTHERMORE, IT FEATURES A 'TWIST-LOCK' TO PREVENT TUBES FROM TURNING DURING SCREW (DE)CAPPING. BESIDES THE STANDARD BARCODED RACKS, MICRONIC ALSO OFFERS THE SERVICE TO SUPPLY LASER-ETCHED 1D BARCODES IN CUSTOMIZED FORMATS.












RACK OVERVIEW AND SPECIFICATIONS

| | LOBORACK-96 | ROBORACK-96 | COMORACK-96 | MICRONIC 96-1 | MICRONIC 96-Q1 | MICRONIC 96-2 | MICRONIC 96-3 | MICRONIC 96-4 |
|---|--|--|--|---|---|---|---|---|
| |  Page 52 |  Page 52 |  Page 52 |  Page 50 |  Page 50 |  Page 50 |  Page 50 |  Page 50 |
| Rack material | PP | PP | PC | PC | PC | PC | PC | PC |
| Quantity of tubes in rack | 96 | 96 | 96 | 96 | 96 | 96 | 96 | 96 |
| Available for tube volumes (nominal) | 0.50ml and 0.75ml (internal thread) | 1.10ml, 1.40ml and 2.50ml (internal thread) | 1.40ml (standard tube) (internal thread) | 0.30ml, 0.40ml and 0.50ml (internal/external thread) | 0.50ml (internal thread) | 0.75ml (internal/external thread) | 0.80ml and 1.10ml (internal/external thread) | 1.40ml (internal/external thread) |
| Rack height without cover | 15.0mm | 28.4mm | 31.1mm | 12.0mm | 16.0mm | 19.0mm | 22.7mm | 30.0mm |
| Tube height in rack without cover <i>(*Tubes with Push Caps and tubes with Low Profile Screw Caps are identical in height)</i> | <i>Internal Thread</i> 21.0mm (0.50ml without cap) 22.0mm (0.50ml with Push Cap*) 28.5mm (0.50ml with Screw Cap) 28.5mm (0.75ml without cap) 29.5mm (0.75ml Push Cap*) 36.0mm (0.75ml Screw Cap) | <i>Internal Thread</i> 39.0mm (1.10ml without cap) 40.0mm (1.10ml with Push Cap*) 46.5mm (1.10ml with Screw Cap) 46.5mm (1.40ml without cap) 47.5mm (1.40ml Push Cap*) 54.5mm (1.40ml Screw Cap) 85.0mm (2.50ml without cap) 86.0mm (2.50ml with Push Cap) | <i>Standard Tube</i> 46.5mm (1.40ml without cap) 47.5mm (1.40ml with Push Cap) | <i>Internal Thread</i> 19.5mm (0.50ml without cap) 20.5mm (0.50ml with Push Cap*) 26.5mm (0.50ml with Screw Cap) <i>External Thread</i> 17.4mm (0.30ml without cap) 18.8mm (0.30ml with Push Cap) 19.6mm (0.30ml with Screw Cap IG) 20.8mm (0.30ml with Screw Cap) 22.8mm (0.40ml without cap) 24.2mm (0.40ml with Push Cap) 25.0mm (0.40ml with Screw Cap IG) 26.3mm (0.40ml with Screw Cap) | <i>Internal Thread</i> 19.5mm (without cap) 20.5mm (with Push Cap*) 27.2mm (with Screw Cap) <i>External Thread</i> 31.0mm (without cap) 32.4mm (with Push Cap) 33.3mm (with Screw Cap IG) 34.5mm (with Screw Cap) | <i>Internal Thread</i> 27.0mm (without cap) 28.0mm (with Push Cap*) 34.7mm (with Screw Cap) <i>External Thread</i> 31.0mm (without cap) 32.4mm (with Push Cap) 33.3mm (with Screw Cap IG) 34.5mm (with Screw Cap) | <i>Internal Thread</i> 37.5mm (1.10ml without cap) 38.5mm (1.10ml with Push Cap*) 45.0mm (1.10ml with Screw Cap) <i>External Thread</i> 35.8mm (0.80ml without cap) 37.2mm (with Push Cap) 38.0mm (with Screw Cap IG) 39.3mm (0.80ml with Screw Cap) 42.3mm (1.10ml without cap) 43.7mm (1.10ml with Push Cap) 44.6mm (1.10ml with Screw Cap IG) 45.8mm (1.10 with Screw Cap) | <i>Internal Thread</i> 45.0mm (without cap) 46.0mm (with Push Cap*) 53.0mm (with Screw Cap) <i>External Thread</i> 49.0mm (without cap) 50.4mm (with Push Cap) 51.3mm (with Screw Cap IG) 52.5mm (with Screw Cap) |
| Rack height with low cover | 23.8mm | 41.2mm | n/a | 23.8mm | 30.8mm | 30.8mm | 41.5mm | 48.8mm |
| Rack height with standard cover | 31.7mm | 49.1mm | 52.0mm | n/a | n/a | n/a | n/a | n/a |
| Rack height with high cover | 38.7mm | 56.1mm | n/a | 30.8mm | n/a | 37.8mm | 48.5mm | 55.8mm |
| Rack height with extra high cover | n/a | 88.0mm | n/a | n/a | n/a | n/a | n/a | n/a |
| Available barcoded | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Available laser etched barcoded | Yes | Yes | No | Yes | Yes | Yes | Yes | Yes |
| Twist-Lock | Yes | Yes | No | Yes | Yes | Yes | Yes | Yes |
| Sterilization option | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| DNase / RNase free | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Sure shot design | Yes | Yes | No | No | No | No | No | No |
| Rack locking system (cover) | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Autoclavable** | Yes | Yes | No | No | No | No | No | No |
| Class 7 clean room production | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Temperature range*** | Vapor phase LN2 to +121°C | Vapor phase LN2 to +121°C | -80°C to +140°C | Vapor phase LN2 to +140°C | Vapor phase LN2 to +140°C | Vapor phase LN2 to +140°C | Vapor phase LN2 to +140°C | Vapor phase LN2 to +140°C |
| Chemical resistance | +++ | +++ | + | ++/+ | ++/+ | ++/+ | ++/+ | ++/+ |

** The racks can be autoclaved in +121°C. Please request more information about the number of cycles per rack type.

*** Temperature range of the material of the tubes, caps and racks is approved from -80°C to +100°C. In practice racks are used in higher temperatures for autoclaving (+121°C) and in lower temperatures until vapor phase LN2 (-150°C to -196°C) for storage. Micronic does not recommend to store racks in liquid phase nitrogen (LN2).

RACK OVERVIEW AND SPECIFICATIONS

| | MICRONIC 96-5 | MICRONIC 48-1 | MICRONIC 48-2 | MICRONIC 48-3 | MICRONIC 48-4 | MICRONIC 24-1 | MICRONIC 24-2 | MICRONIC 24-4 | MICRONIC 138-2 |
|---|---|--|---|--|--|--|--|--|---|
| |  Page 50 |  Page 54 |  Page 54 |  Page 54 |  Page 54 |  Page 56 |  Page 56 |  Page 56 |  Page 58 |
| Rack material | PC | PC | PC | PC | PC | PC | PC | PC | PC |
| Quantity of tubes in rack | 96 | 48 | 48 | 48 | 48 | 24 | 24 | 24 | 138 |
| Available for tube volumes (nominal) | 2.00ml (internal/external thread) | 1.00ml (external thread) | 2.00ml (external thread hybrid) | 3.00ml (external thread hybrid) | 4.00ml (internal/external thread) | 1.50ml and 3.00ml (external thread hybrid) | 3.50ml (external thread hybrid) | 6.00ml (internal/external thread hybrid) | 0.75ml and 0.80ml (internal/external thread hybrid) |
| Rack height without cover | 41.0mm | 12.0mm | 19.0mm | 22.7mm | 30.0mm | 12.0mm | 19.0mm | 30.0mm | 20.2mm |
| Tube height in rack without cover <i>(*Tubes with Push Caps and tubes with Low Profile Screw Caps are identical in height)</i> | <i>Internal Thread</i> 60.0mm (without cap) 61.0mm (with Push Cap*) 67.5mm (with Screw Cap) <i>External Thread</i> 64.0mm (without cap) 65.4mm (with Push Cap) 66.3mm (with Screw Cap IG) 67.5mm (with Screw Cap) | <i>External Thread</i> 17.0mm (without cap) 20.7mm (with Screw Cap) | <i>External Thread</i> 31.0mm (without cap) 34.5mm (with Screw Cap) | <i>External Thread</i> 42.0mm (without cap) 45.5mm (with Screw Cap) | <i>Internal Thread</i> 45.0mm (without cap) 52.4mm (with Screw Cap) <i>External Thread</i> 49.0mm (without cap) 52.5mm (with Screw Cap) | <i>External Thread</i> 17.0mm (1.50ml without cap) 20.5mm (1.50ml with Screw Cap) 24.0mm (3.00ml without cap) 27.5mm (3.00ml with Screw Cap) | <i>External Thread</i> 31.2mm (without cap) 34.7mm (with Screw Cap) | <i>Internal Thread</i> 44.0mm (without cap) 51.9mm (with Screw Cap) <i>External Thread</i> 49.0mm (without cap) 52.5mm (with Screw Cap) | <i>Internal Thread</i> 27.0mm (0.75ml without cap) 28.0mm (0.75ml with Push Cap*) 34.5mm (0.75ml with Screw Cap) <i>External Thread</i> 31.0mm (0.75ml without cap) 32.4mm (0.75ml with Push Cap) 33.3mm (0.75ml with Screw Cap IG) 34.5mm (0.75ml with Screw Cap) 35.8mm (0.80ml without cap) 37.2mm (0.80ml with Push Cap) 38.1mm (0.80ml with Screw Cap IG) 39.3mm (0.80ml with Screw Cap) |
| Rack height with low cover | 66.8mm | 23.8mm | n/a | n/a | n/a | 23.8mm | n/a | n/a | n/a |
| Rack height with high cover | 74.3mm | n/a | 30.8mm | 41.5mm | 55.8mm | 30.8mm | 37.8mm | 55.8mm | 55.8mm |
| Available barcoded | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Available laser etched barcoded | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Twist-Lock | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Sterilization option | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| DNase / RNase free | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Sure shot design | No | No | No | No | No | No | No | No | No |
| Rack locking system (cover) | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Autoclavable** | No | No | No | No | No | No | No | No | No |
| Class 7 clean room production | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Temperature range*** | Vapor phase LN2 to +140°C | Vapor phase LN2 to +140°C | Vapor phase LN2 to +140°C | Vapor phase LN2 to +140°C | Vapor phase LN2 to +140°C | Vapor phase LN2 to +140°C | Vapor phase LN2 to +140°C | Vapor phase LN2 to +140°C | Vapor phase LN2 to +140°C |
| Chemical resistance | ++/+ | ++/+ | ++/+ | ++/+ | ++/+ | ++/+ | ++/+ | ++/+ | ++/+ |

** The racks can be autoclaved in +121°C. Please request more information about the number of cycles per rack type.

*** Temperature range of the material of the tubes, caps and racks is approved from -80°C to +100°C. In practice racks are used in higher temperatures for autoclaving (+121°C) and in lower temperatures until vapor phase LN2 (-150°C to -196°C) for storage. Micronic does not recommend to store racks in liquid phase nitrogen (LN2).

The Micronic ULT Rack Range is developed to securely store samples down to ultra-low temperatures. Absolute traceability and reproducibility on the racks is ensured through novel laser-etched alphanumeric visual location aids on top of the racks and a unique laser etched 1D barcode on the side of the racks. The open bottom design of the storage racks facilitate quick defrosting of samples.



ULT RACK RANGE



Resistance to extreme conditions

SUPERIOR FEATURES

- Dimensions based on the Society for Laboratory Automation and Screening (ANSI / SLAS) standards for storage racks
- Extremely strong design which minimizes deformation due to temperature changes
- Manufactured from Polycarbonate which has excellent properties for low temperatures (vapor phase LN2)
- Highly transparent cover
- Standard available with laser etched 1D rack barcode
- Unique laser etched Alphanumeric coding on top
- Rack orientation marks on bottom and side
- Secure and safe locking system
- Twist-Lock
- Open rack bottom
- Compatible with the Micronic Rack Readers
- Class 7 clean room production
- RNase/DNase and Pyrogen free
- 2D Data-Matrix code on the rack bottom


ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Customized 1D rack barcode

SPECIFICATIONS

| | | |
|---------------|--|--|
| | | Compatible with tubes from 0.30ml to 2.00ml |
| | | 12 rows of 8 tubes (96 tubes) |
| | | Polycarbonate |
| | | Length: 127.8mm (5.03 inches), width: 85.5mm (3.37 inches) |
| | | Resistant to Chemical solvents |
| See next page | | Vapor phase LN2 to +140°C |

DIMENSIONS

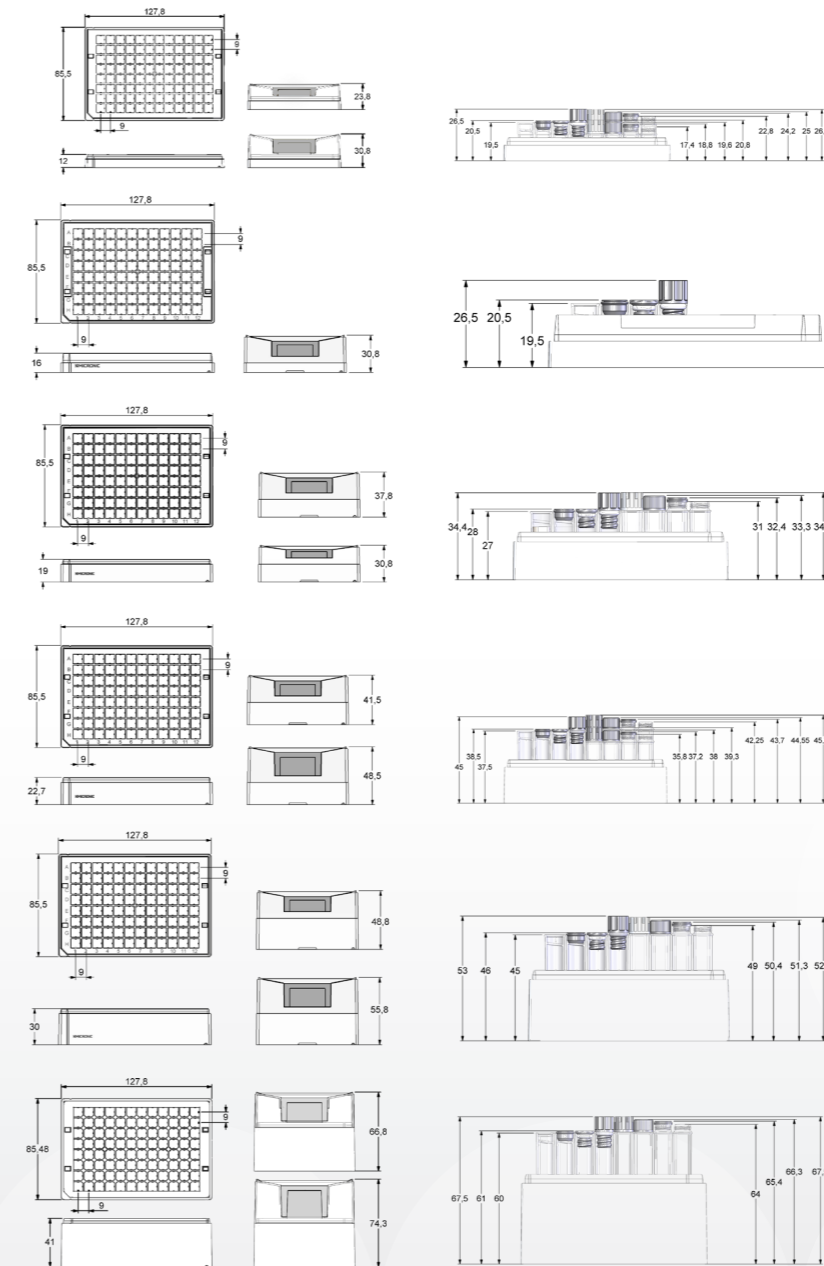

Micronic 96-1 rack dimensions and tube in rack dimensions
Left: 0.50ml internally threaded tubes with TPE Push Caps / Low Profile Screw Caps or Screw Caps. Right: 0.30ml / 0.40ml externally threaded tubes with Screw Caps / Screw Caps IG or Push Caps.

Micronic 96-Q1 rack dimensions and tube in rack dimensions
0.50ml internally threaded tubes with TPE Push Caps / Low Profile Screw Caps or Screw Caps.

Micronic 96-2 rack dimensions and tube in rack dimensions
Left: 0.75ml internally threaded tubes with TPE Push Caps / Low Profile Screw Caps or Screw Caps. Right: 0.75ml externally threaded tubes (Hybrid) with Screw Caps / Screw Caps IG or Push Caps.

Micronic 96-3 rack dimensions and tube in rack dimensions
Left: 1.10ml internally threaded tubes with TPE Push Caps / Low Profile Screw Caps or Screw Caps. Right: 0.80ml / 1.10ml externally threaded tubes with Screw Caps / Screw Caps IG or Push Caps.

Micronic 96-4 rack dimensions and tube in rack dimensions
Left: 1.40ml internally threaded tubes with TPE Push Caps / Low Profile Screw Caps or Screw Caps. Right: 1.40ml externally threaded tubes (Hybrid) with Screw Caps / Screw Caps IG or Push Caps.

Micronic 96-5 rack dimensions and tube in rack dimensions
Left: 2.00ml internally threaded tubes with TPE Push Caps / Low Profile Screw Caps or Screw Caps. Right: 2.00ml externally threaded tubes with Screw Caps / Screw Caps IG or Push Caps.

ORDERING INFORMATION

| Catalog number | Description | Case Size |
|--|--|--------------|
| Tube storage racks in 96-well ANSI/SLAS standard format (empty) | | |
| MP51200 | Micronic 96-1 Rack with low cover for 0.30ml tubes capped with Screw Caps | 4 x 10 racks |
| MP51201 | Micronic 96-1 Rack with low cover for 0.50ml tubes capped with Push Caps or (Low Profile) Screw Caps | 4 x 10 racks |
| MP51204 | Micronic 96-2 Rack with low cover for 0.75 ml tubes capped with Push Caps or Low Profile Screw Caps | 4 x 10 racks |
| MP51217 | Micronic 96-3 Rack with low cover for 0.80ml tubes capped with Screw Caps | 4 x 10 racks |
| MP51207 | Micronic 96-3 Rack with low cover for 1.10ml tubes capped with Push Caps or Low Profile Screw Caps | 4 x 10 racks |
| MP51210 | Micronic 96-4 Rack with low cover for 1.40ml tubes capped with Push Caps or Low Profile Screw Caps | 4 x 10 racks |
| MP51215 | Micronic 96-5 Rack with low cover for 2.00ml tubes capped with Push Caps or Low Profile Screw Caps | 4 x 6 racks |

Other item numbers are available with high cover and bottom rack ID

Precision manufactured in a certified class 7 clean room production facility, these racks are based on the Society for Laboratory Automation and Screening (ANSI/SLAS) standards for storage racks. The racks are designed to hold 96 individual storage tubes, also known as a 96-well configuration.



SUPERIOR FEATURES

- Dimensions based on the Society for Laboratory Automation and Screening (ANSI / SLAS) standards for storage racks
- Standard available with laser etched 1D rack barcode (Loborack-96 and Roborack-96)
- Secure and safe locking system
- Twist-Lock (Loborack-96 and Roborack-96)
- Sure shot design: tapered wells allow precise tube placement into racks (Loborack-96 and Roborack-96)
- Manufactured from Polypropylene (Loborack-96 and Roborack-96) or Polycarbonate (Comorack-96)
- Open rack bottom
- Compatible with the Micronic Rack Readers
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Customized 1D rack barcode
- Pre-labelled rack barcode (Comorack-96)




SPECIFICATIONS


| | | |
|----------------------|--|---|
| <p>See next page</p> | | Compatible with tubes from 0.50ml to 2.00ml |
| | | 12 rows of 8 tubes (96 tubes) |
| | | Polypropylene (Loborack-96 and Roborack-96) / Polycarbonate (Comorack-96) |
| | | Length: 127.8mm (5.03 inches), width: 85.5mm (3.37 inches) |
| | | Resistant to Chemical solvents |
| | | Vapor phase LN2 to +121°C |

DIMENSIONS


Loborack-96 dimensions and tube in rack dimensions: **0.50ml (right) and 0.75ml (left) internally threaded tubes.**

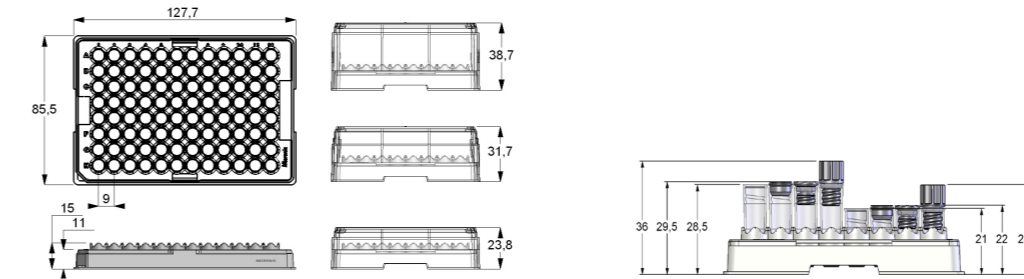

Roborack-96 dimensions and tube in rack dimensions: **1.10ml (lower right), 1.40ml (left) and 2.50ml (upper right) internally threaded tubes.**

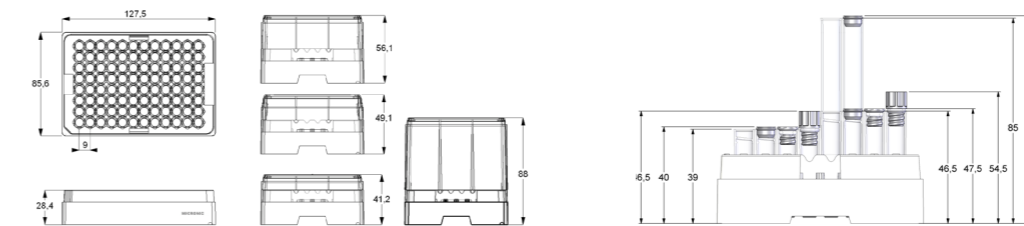

Comorack-96 dimensions and tube in rack dimensions: **standard 1.40ml internally threaded tubes.**

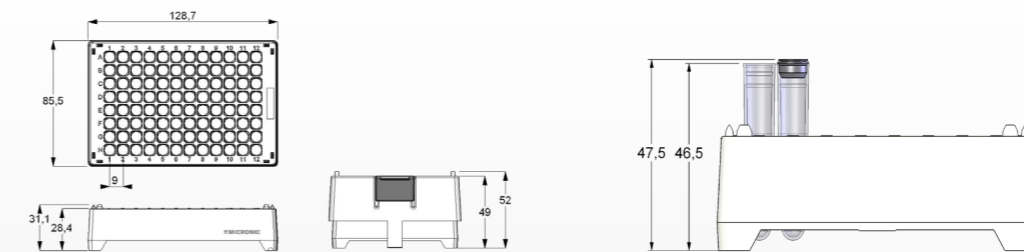

ORDERING INFORMATION

| Catalog number | Description | Case Size |
|--|---|--------------|
| Tube storage racks in 96-well ANSI/SLAS standard format (empty) | | |
| MPW51013BC3 | Loborack-96 White, Laser etched (including low cover) (barcoded A1-H1 side) | 4 x 10 racks |
| MPW51004BC3 | Loborack-96 White, Laser etched (including standard cover) (barcoded A1-H1 side) | 4 x 10 racks |
| MPW51015BC3 | Loborack-96 White, Laser etched (including high cover) (barcoded A1-H1 side) | 4 x 10 racks |
| MPW51001BC3 | Roborack-96 White, Laser etched (including standard cover) (barcoded A1-H1 side) | 4 x 10 racks |
| MPW51012BC3 | Roborack-96 White, Laser etched (including high cover) for 1.40ml tubes precapped with Screw Caps (barcoded A1-H1 side) | 2 x 10 racks |
| MPW51016BC3 | Roborack-96 White, Laser etched (including extra high cover) for 2.50ml tubes (barcoded A1-H1 side) | 4 x 6 racks |
| MPW51017BC3 | Roborack-96 White, Laser etched (including low cover) (barcoded A1-H1 side) | 4 x 10 racks |
| MP22504 | Comorack-96 (including cover) (polycarbonate) | 4 x 10 racks |

The Micronic 48-well format ULT Racks can hold 6x8 (48) individual tubes. Next to the standard ANSI/SLAS footprint, the ULT Racks are provided with more features ideal for automation. For example: an outstanding Twist-Lock design which prevents tubes from turning during screw (de)capping, easy lead-in tube and cover placement and closed side walls with minimal notches.



SUPERIOR FEATURES

- Dimensions based on the Society for Laboratory Automation and Screening (ANSI / SLAS) standards for storage racks
- Extremely strong design which minimizes deformation due to temperature changes
- Manufactured from Polycarbonate which has excellent properties for low temperatures (vapor phase LN2)
- Highly transparent cover
- Standard available with laser etched 1D rack barcode
- Unique laser etched Alphanumeric coding on top
- Rack orientation marks on bottom and side
- Secure and safe locking system
- Twist-Lock
- Open rack bottom
- Compatible with the Micronic Rack Readers
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Customized 1D rack barcode




SPECIFICATIONS

| | | |
|--|--|--|
| | | Compatible with tubes from 1.00ml to 4.00ml |
| | | 8 rows of 6 tubes (48 tubes) |
| | | Polycarbonate |
| | | Length: 127.8mm (5.03 inches), width: 85.5mm (3.37 inches) |
| | | Resistant to Chemical solvents |
| | | Vapor phase LN2 to +140°C |

See next page

DIMENSIONS

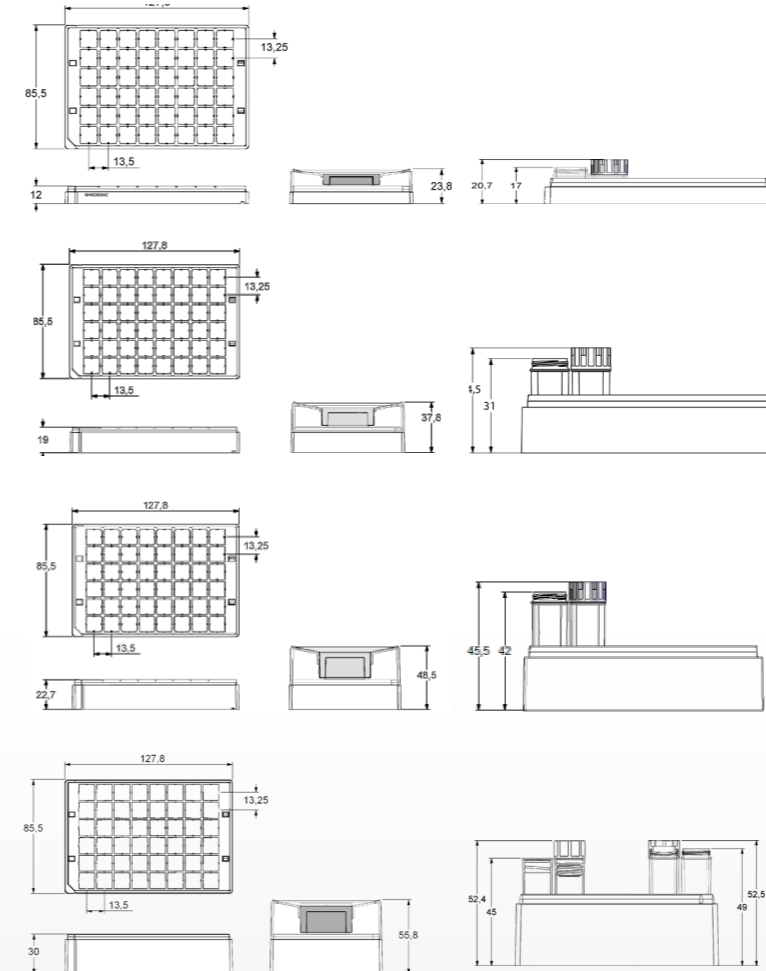

Micronic 48-1 rack dimensions and tube in rack dimensions **1.00ml externally threaded tubes with Screw Caps.**

Micronic 48-2 rack dimensions and tube in rack dimensions **2.00ml externally threaded Hybrid tubes with Screw Caps.**

Micronic 48-3 rack dimensions and tube in rack dimensions **3.00ml externally threaded Hybrid tubes with Screw Caps.**

Micronic 48-4 rack dimensions and tube in rack dimensions **Left: 4.00ml internally threaded tubes with Screw Caps. Right: 4.00ml externally threaded tubes with Screw Caps.**

ORDERING INFORMATION

| Catalog number | Description | Case Size |
|--|--|--------------|
| Tube storage racks in 48-well ANSI/SLAS standard format (empty) | | |
| MP51304 | Micronic 48-1 Rack with low cover for 1.00ml tubes capped with Screw Caps | 4 x 10 racks |
| MP51303 | Micronic 48-2 Rack with high cover for 2.00ml tubes capped with Screw Caps | 4 x 10 racks |
| MP51302 | Micronic 48-3 Rack with high cover for 3.00ml tubes capped with Screw Caps | 4 x 10 racks |
| MP51301 | Micronic 48-4 Rack with high cover for 4.00ml tubes capped with Screw Caps | 2 x 10 racks |

Micronic ULT Rack

The Fastest Way to Expert Opinion™

★★★★☆

Sandra Story
Status: Reviewer
Member since: 2020
Organization: NUBAD, LLC

Ease of use ★★★★★
After sales service ★★★★★
Value for money ★★★★★

“Everyone loves this product that helps efficiency and organization.”

Rating: 4.3 ★★★★★

Application Area: Drug development stability storage

“The multitube rack is a perfect fit for our liquid handler. Keeping track of inventory is no longer a complicated chore and we have confidence in the long term storage stability and that our product can be shipped safely.”

Review date: 27 Mar 2020 | Micronic ULT 96-4 Rack

The Micronic 24-well format racks are made from Polycarbonate and have excellent properties for low temperature conditions. The material used to manufacture the racks is selected on quality and the highest purity levels. The racks have an extremely strong design which minimizes deformation due to temperature changes.



SUPERIOR FEATURES

- Dimensions based on the Society for Laboratory Automation and Screening (ANSI / SLAS) standards for storage racks
- Extremely strong design which minimizes deformation due to temperature changes
- Manufactured from Polycarbonate which has excellent properties for low temperatures (vapor phase LN2)
- Highly transparent cover
- Standard available with laser etched 1D rack barcode
- Unique laser etched Alphanumeric coding on top
- Rack orientation marks on bottom and side
- Secure and safe locking system
- Twist-Lock
- Open rack bottom
- Compatible with the Micronic Rack Readers
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS


- Sterilization using gamma irradiation or Eto treatment
- Customized 1D rack barcode



SPECIFICATIONS

| | | |
|---------------|--|--|
| | | Compatible with tubes from 1.50ml to 6.00ml |
| | | 6 rows of 4 tubes (24 tubes) |
| | | Polycarbonate |
| | | Length: 127.8mm (5.03 inches), width: 85.5mm (3.37 inches) |
| | | Resistant to Chemical solvents |
| See next page | | Vapor phase LN2 to +140°C |

DIMENSIONS

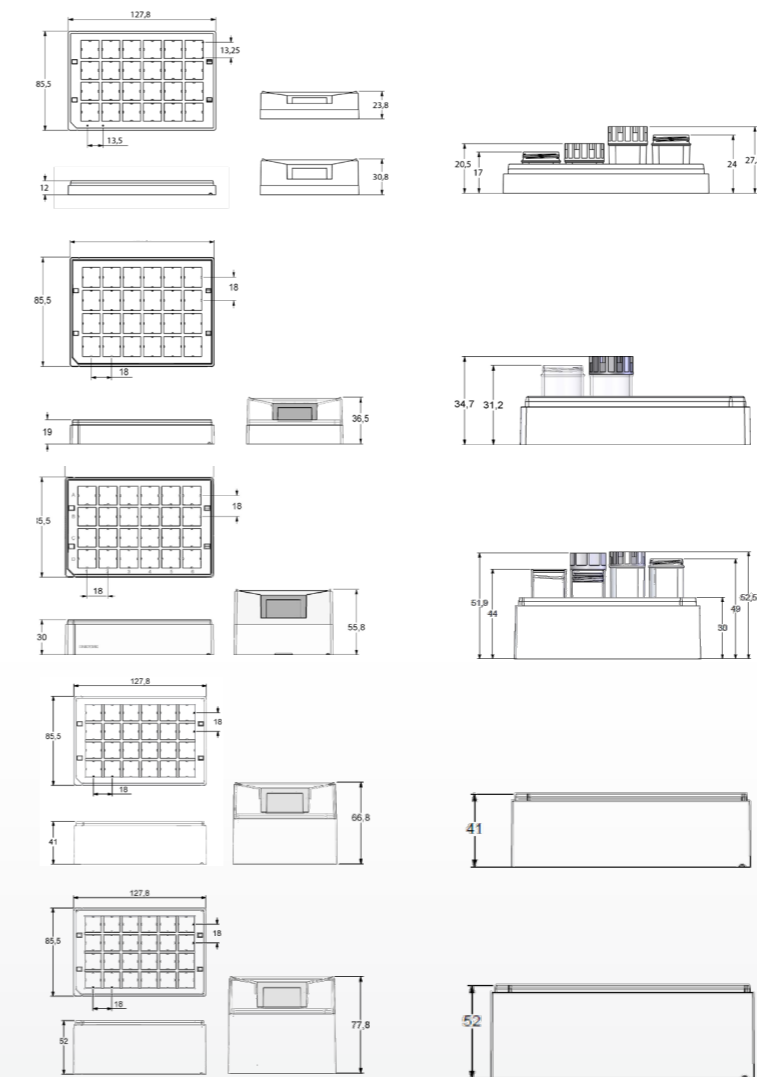

Micronic 24-1 rack dimensions and tube in rack dimensions
Left: 1.50ml externally threaded (Hybrid) tubes with Screw Caps. Right: 3.00ml externally threaded (Hybrid) tubes with Screw Caps.

Micronic 24-2 rack dimensions and tube in rack dimensions
3.50ml externally threaded tubes with Screw Caps.

Micronic 24-4 rack dimensions and tube in rack dimensions
Left: 6.00ml internally threaded tubes with Screw Caps. Right: 6.00ml externally threaded (Hybrid) tubes with Screw Caps.

Micronic 24-5 rack dimensions
High volume tubes.

Micronic 24-6 rack dimensions
High volume tubes.

ORDERING INFORMATION

| Catalog number | Description | Case Size |
|--|--|--------------|
| Tube storage racks in 24-well ANSI/SLAS standard format (empty) | | |
| MP51404 | Micronic 24-1 Rack with low cover for 1.50ml tubes capped with Screw Caps | 4 x 10 racks |
| MP51406 | Micronic 24-1 Rack with high cover for 3.00ml tubes capped with Screw Caps | 4 x 10 racks |
| MP51409 | Micronic 24-2 Rack with high cover for 3.50ml tubes capped with Screw Caps | 4 x 10 racks |
| MP51403 | Micronic 24-4 Rack with high cover for 6.00ml tubes capped with Screw Caps | 2 x 10 racks |
| MP51407 | Micronic 24-5 Rack with low cover for tubes capped with Screw Caps | 4 x 6 racks |
| MP51408 | Micronic 24-6 Rack with low cover for tubes capped with Screw Caps | 4 x 6 racks |

The Micronic 138-well format HD rack is made to store more samples in an ANSI/SLAS format rack, and is especially designed to be compatible with automated processes. The rack walls are closed for easy gripping with most robotic arms. In addition, the rack uses a unique open bottom design, which permits tubes to be manipulated by robotic systems.



SUPERIOR FEATURES








- Dimensions based on the Society for Laboratory Automation and Screening (ANSI / SLAS) standards for storage racks
- Increases the number of samples that can be stored up to 43% compared to a standard rack in the ANSI/SLAS footprint
- Extremely strong design which minimizes deformation due to temperature changes
- Manufactured from Polycarbonate which has excellent properties for low temperatures (vapor phase LN2)
- Rack side walls are closed for easy and various gripping of most robotic arms
- Standard available with laser etched 1D rack barcode
- Twist-Lock
- Open rack bottom
- Compatible with automated processes
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Customized 1D rack barcode

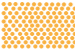


SPECIFICATIONS

| | | |
|--|---|--|
|  |  | Compatible with tubes from 0.75ml to 0.80ml |
| |  | 6 rows of 13 tubes and 5 rows of 12 tubes (138 tubes) |
| |  | Polycarbonate |
| |  | Length: 127.8mm (5.03 inches), width: 85.5mm (3.37 inches) |
| |  | Resistant to Chemical solvents |
| |  | Vapor phase LN2 to +100°C |

See next page

DIMENSIONS

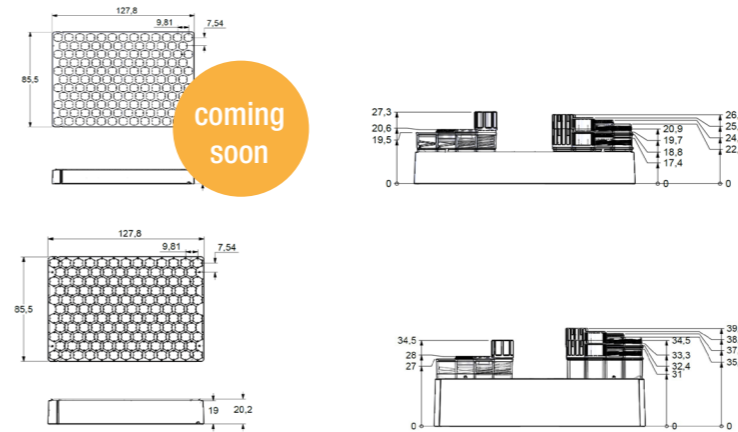

Micronic 138-1 rack dimensions and tube in rack dimensions
Left: 0.50ml internally threaded tubes with TPE Push Caps / Low Profile Screw Caps or Screw Caps. Right: 0.30ml / 0.40ml externally threaded tubes with TPE Push Caps / Screw Caps IG or Screw Caps.

Micronic 138-2 rack dimensions and tube in rack dimensions
Left: 0.75ml internally threaded tubes with TPE Push Caps / Low Profile Screw Caps or Screw Caps. Right: 0.75ml / 0.80ml externally threaded (hybrid) tubes with TPE Push Caps / Screw Caps IG or Screw Caps.

ORDERING INFORMATION

| Catalog number | Description | Case Size |
|----------------|---|--------------|
| MP51420 | Micronic 138-2 HD Rack without cover for 0.75ml tubes (empty) (Barcoded A1-H1 side) | 4 x 10 racks |



SERVING SCIENTISTS
SINCE 1984

CERTIFIED CLASS 7 CLEAN ROOM
MANUFACTURING

SAMPLE STORAGE CAPS

MICRONIC OFFERS THE MOST EXTENSIVE RANGE OF TUBE CAPPING PRODUCTS, INCLUDING SCREW CAPS AND PUSH CAPS, IN THE MARKET. FOR EFFECTIVE CRYOGENIC STORAGE, A SCREW CAP IS RECOMMENDED SINCE IT ALLOWS FOR THE TIGHTEST SEAL POSSIBLE. FOR STORAGE FROM -80°C AND ABOVE, EITHER A PUSH CAP (TPE CAPCLUSTER) OR A SCREW CAP CAN BE USED. ALL TUBES HAVE AN UPPER SCREW THREAD OR DOUBLE O-RING DESIGN PROVIDING A PERFECT SEAL AND CLOSURE ENSURING MAXIMUM SAMPLE INTEGRITY. THE MICRONIC PUSH CAPS ARE AVAILABLE IN 96-FORMAT AND THE SCREW CAPS IN 96-, 48- AND 24-FORMAT. ALL SEALING PRODUCTS ARE AVAILABLE IN A RANGE OF COLORS!

Industry
leading cap
range

SUPERIOR FEATURES OF MICRONIC SAMPLE STORAGE CAPS

All Screw Caps are manufactured from virgin plastic raw material, RNase/DNase and pyrogen free, automation friendly and have a temperature range of: Vapor phase LN2 to +100°C

High quality Silicon O-ring allows for optimal sealing for tubes with internal thread

Available in 12 colors



Available in 12 colors



Triple start thread allows for the tightest seal possible for tubes with external thread

TPE Push Caps for tubes with internal thread & external thread

Multi-pierceable without losing sealing integrity

Available with split septum to pierce with a pipette tip

RNase / DNase and Pyrogen free



Temperature range: -80°C to +100°C

TPE Caps in Natural: manufactured from materials for medical or non-medical applications

Available in Natural + 12 colors



STANDARD MICRONIC CAP COLOR CODING GUIDE FOR SAMPLE STORAGE TUBES





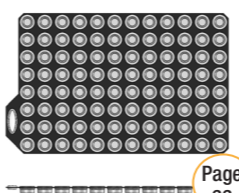
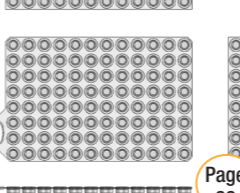
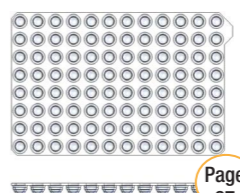
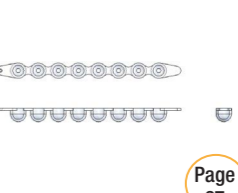
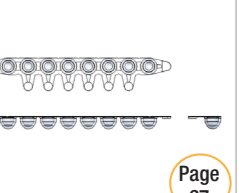
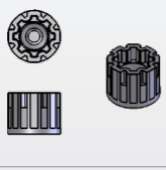
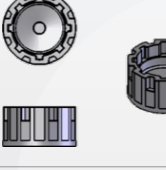
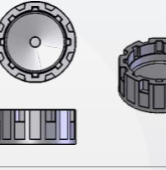
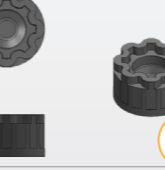
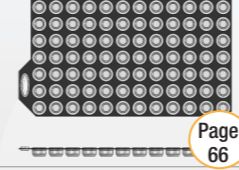
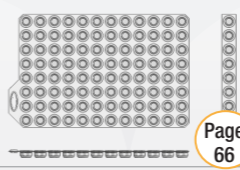
During the collection, processing and storage of samples, sample traceability is crucial. While most laboratories use 1D and 2D barcodes on sample tubes to ensure traceability of samples, cap color inspection will further promote the tube identification process. Cap color coding in laboratories is particularly useful when used to distinguish between different types of research purposes. Color provides a visual clue that works faster than text alone. By adding cap color coding next to 1D and 2D barcodes, users can speed up sample storage and retrieval time, eliminate human errors, and reduce the risk of lost samples.

To ensure that aliquots, which are transferred from sample collection tubes (vacuum tubes) into smaller sample storage tubes can still be identified by cap color inspection, Micronic offers a standard cap color coding solution for 96-, 48- and 24-format sample storage tubes. This solution is related to the cap colors of vacuum tubes. To each type of blood sample (i.e. additive which is present to pre-treat or preserve the blood), a cap color is assigned. Next to blood samples, the cap color solution can also be used for other biological samples (e.g. DNA, tissue, urine, feces, plants, or seeds).

For more information about the standard cap color coding guide, please see Micronic's Tech Note 'Cap Color Coding' on our website www.micronic.com!



CAP OVERVIEW AND SPECIFICATIONS
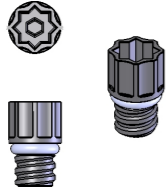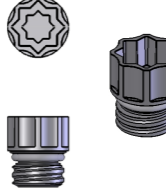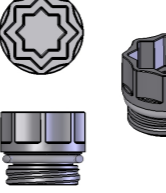
| | SCREW CAP-96 | SCREW CAP-48 | SCREW CAP-24 | LOW PROFILE SC | TPE CAPCLUSTER | TPE CAPMAT | EVA CAPMAT | EVA CAPBAND | EVA BREAK-OFF |
|--|--|--|---|--|--|--|---|--|--|
| Caps for tubes with internal thread |  Page 64 |  Page 64 |  Page 64 |  Page 64 |  Page 66 |  Page 66 |  Page 67 |  Page 67 |  Page 67 |
| Cap material | PP + Silicon O-ring | PP + Silicon O-ring | PP + Silicon O-ring | PP + Silicon O-ring | TPE | TPE (LYO) | EVA | EVA | EVA |
| Capping for # tubes | 1 | 1 | 1 | 1 | 96 or 1 | 96, 12 or 8 | 96 | 8 | 8 or 1 |
| Format | Single cap for 96-format tubes | Single cap for 48-format tubes | Single cap for 24-format tubes | Single cap for 96-format tubes | Mat of 96 individual caps | Mat of 96 caps | Mat of 96 caps | Strip of 8 caps | Strip of 8 caps |
| Compatible tubes | 0.50ml, 0.75ml, 1.10ml, 1.40ml, 2.00ml | 4.00ml | 6.00ml | 0.50ml, 0.75ml, 1.10ml, 1.40ml, 2.00ml | 0.50ml, 0.75ml, 1.10ml, 1.40ml, 2.00ml, 2.50ml | 0.50ml, 0.75ml, 1.10ml, 1.40ml, 2.00ml, 2.50ml | 0.50ml, 0.75ml, 1.10ml, 1.40ml, 2.00ml, 2.50ml | 0.50ml, 0.75ml, 1.10ml, 1.40ml, 2.00ml, 2.50ml | 0.50ml, 0.75ml, 1.10ml, 1.40ml, 2.00ml, 2.50ml |
| Colors* | 12 different colors | 12 different colors | 12 different colors | 12 different colors | Natural + 12 different colors | Natural + 7 different colors | Transparent | Transparent | Transparent |
| Pierceable | No | No | No | No | Yes | Yes | No | No | No |
| Available as pre-split | No | No | No | No | Yes | Yes | No | No | No |
| Re-usable | Yes | Yes | Yes | Yes | No | No | No | No | No |
| Sterilization | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Available DNase / RNase free | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Class 7 clean room production | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes | Yes |
| Minimum storage temperature** | Vapor phase LN2 | Vapor phase LN2 | Vapor phase LN2 | Vapor phase LN2 | -80°C | -80°C | -40°C | -40°C | -40°C |
| Chemical resistance | ++ | ++ | ++ | ++ | +++ | +++ | ++ | ++ | ++ |
| | SCREW CAP-96 | SCREW CAP-48 | SCREW CAP-24 | SCREW CAP IG | TPE CAPCLUSTER | TPE CAPMAT | <p>* Please be aware that great care was taken to develop and thoroughly test the caps to secure the stability of the colors. However, as it is not possible to cover all applications, we recommend users to test the colored caps within their own applications.</p> <p>** Temperature range of the material of the tubes, caps and racks is approved from -80°C to +100°C. In practice Screw Caps are used in lower temperatures until <i>vapor phase LN2</i> (-150°C to -196°C). Micronic does not recommend to store tubes in liquid phase nitrogen. In practice tubes are only snap frozen in LN2 (without submerging the cap).</p> | | |
| Caps for tubes with external thread |  Page 65 |  Page 65 |  Page 65 |  Page 65 |  Page 66 |  Page 66 | | | |
| Cap material | PE | PE | PE | PE | TPE | TPE | | | |
| Capping for # tubes | 1 | 1 | 1 | 1 | 96 or 1 | 96 or 1 | | | |
| Format | Single cap for 96-format tubes | Single cap for 48-format tubes | Single cap for 24-format tubes | Single cap for 96-format tubes | Mat of 96 individual caps | Mat of 96 caps | | | |
| Compatible tubes | 0.30ml, 0.40ml, 0.75ml, 0.80ml, 1.40ml, 2.00ml | 1.00ml, 2.00ml, 3.00ml 4.00ml | 1.50ml, 3.00ml, 3.50ml, 6.00ml | 0.30ml, 0.40ml, 0.75ml, 0.80ml, 1.40ml, 2.00ml | 0.30ml, 0.40ml, 0.75ml, 0.80ml, 1.40ml, 2.00ml | 0.30ml, 0.40ml, 0.75ml, 0.80ml, 1.40ml, 2.00ml | | | |
| Colors* | 12 different colors and transparent | 12 different colors and transparent | 12 different colors and transparent | 12 different colors and transparent | Natural + 12 different colors | Natural + 7 different colors | | | |
| Re-usable | Yes | Yes | Yes | Yes | No | No | | | |
| Sterilization option | Yes | Yes | Yes | Yes | Yes | Yes | | | |
| DNase / RNase free | Yes | Yes | Yes | Yes | Yes | Yes | | | |
| Class 7 clean room production | Yes | Yes | Yes | Yes | Yes | Yes | | | |
| Minimum storage temperature** | Vapor phase LN2 | Vapor phase LN2 | Vapor phase LN2 | Vapor phase LN2 | -80°C | -80°C | | | |
| Chemical resistance | ++ | ++ | ++ | ++ | +++ | +++ | | | |



Micronic Screw Caps-96 provide a secure closure for long-term sample storage at ultra-low temperatures or a safe sealing during the transportation of your samples. To meet the diverse needs of different researchers, Micronic offers a range of sealing solutions. Tubes with Internal thread can be sealed with the standard Screw Cap, the Screw Cap Ultra and the Low Profile Screw Cap.



INTERNAL THREAD

SUPERIOR FEATURES

- Low Profile Screw Caps save freezer space up to 33%!
- Screw Caps feature a high quality Silicon O-ring
- Available in 12 different colors: Grey, White, Yellow, Orange, Red, Pink, Purple, Blue, Light Blue, Light Green, Green and Black
- Automation friendly
- Ideal for sample storage at ultra-low temperatures or for sample transportation
- Compatible with Micronic internally threaded Screw Cap tubes in 96-, 48- or 24-format
- Small packaging of 960, 480, 240 caps
- Resists chemical solvents
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Pre-capped on Micronic Screw Cap tubes



SPECIFICATIONS

| | | | |
|--|--|--|---|
| | | | Compatible with Screw Cap tubes from 0.50ml to 6.00ml |
| | | | Compatible with tubes in 96-, 48- and 24-well format |
| | | | Polypropylene with Silicon O-ring |
| | | | See drawings on the left |
| | | | Resistant to Chemical solvents |
| | | | Vapor phase LN2 to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 68.

Tubes with External Thread can be sealed by standard Screw Caps or the Screw Caps IG for Tubes with External Thread. Using externally threaded storage tubes eliminates the possibility of the sample coming into contact with the screw thread, minimizing chance of cross-contamination while improving sample integrity.



EXTERNAL THREAD

SUPERIOR FEATURES

- Triple start thread for optimal sealing quality
- Available in transparent and 12 different colors: Grey, White, Yellow, Orange, Red, Pink, Purple, Blue, Light Blue, Light Green, Green and Black
- Automation friendly
- Screw Cap IG has an improved gripping for both manually and automated handling
- The Screw Cap IG has a newly designed lead-in on top of the Screw Cap which ensures a better tube placement
- Ideal for sample storage at ultra-low temperatures or for sample transportation
- Compatible with Micronic externally threaded Screw Cap tubes in 96-, 48- or 24-format
- Small packaging of 960, 480 or 240 caps
- Resists chemical solvents
- Class 7 clean room production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Pre-capped on Micronic Screw Cap tubes



SPECIFICATIONS

| | | | |
|--|--|--|---|
| | | | Compatible with Screw Cap tubes from 0.30ml to 6.00ml |
| | | | Compatible with tubes in 96-, 48- and 24-well format |
| | | | Polyethylene |
| | | | See drawings on the left |
| | | | Resistant to Chemical solvents |
| | | | Vapor phase LN2 to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 68.

TPE PUSH CAPS

INTERNAL AND EXTERNAL THREAD

The TPE Push Caps are compatible with tubes in 96-well format (internal and external thread). Several TPE cap formats are available: the Capcluster-96 which leaves each tube individually capped, the Capmat-96 holds 96 caps in a fixed foil in which the tubes are sealed together, and the Capband-8 or -12 comes in a mat format with tear off lines to obtain cap strips.




INTERNAL THREAD & EXTERNAL THREAD

SUPERIOR FEATURES

- Caps can be pierced repeatedly without losing seal integrity
- Available in Natural and 12 different colors: Grey, White, Yellow, Orange, Red, Pink, Purple, Blue, Light Blue, Light Green, Green and Black (Capcluster-96)
- TPE Caps in Natural: manufactured from materials for medical or non-medical applications
- Capmats and Capclusters (patented) are compatible with tubes with internal and external thread in 96-format
- Economical solution storage at temperatures down to -80°C
- Resists chemical solvents
- Class 7 clean room production
- RNase/DNase and Pyrogen free
- Special TPE push cap for tubes with internal thread: Lyo Caps-96, which are specially developed for the lyophilization process as vents in the caps allow vapor to escape from the tubes

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment
- Pre-capped on Micronic tubes
- Split septum to pierce with a pipette tip

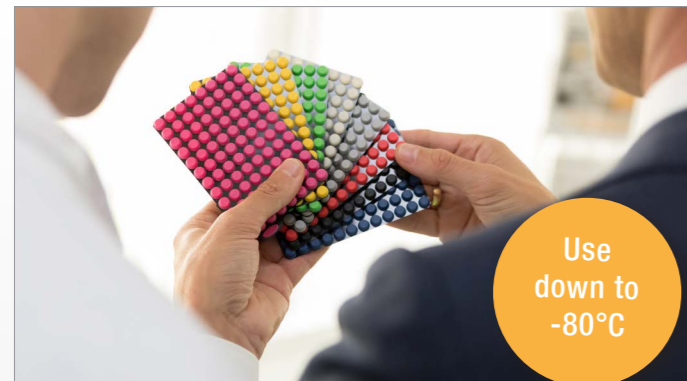

Use down to -80°C

SPECIFICATIONS

| | | |
|--|--|--|
| | | Compatible with Push Cap and Screw Cap tubes from 0.30ml to 2.00ml |
| | | Compatible with tubes in 96-well format |
| | | Thermo Plastic Elastomer (Virgin) |
| | | See drawing on the left |
| | | Resistant to Chemical solvents |
| | | -80°C to +100°C (without retaining foil) |

ORDERING INFORMATION

More information about product numbers and ordering, see page 69.

EVA PUSH CAPS

INTERNAL THREAD

The EVA Push Caps are a cost effective way to securely close sample storage tubes in a 96-well format. The EVA Push Caps for tubes in 96-well format are compatible with the 0.50ml, 0.75ml, 1.10ml, 1.40ml and 2.00ml. The caps are compatible with Micronic Capping Equipment.



INTERNAL THREAD

SUPERIOR FEATURES

- Easy solution
- Available in Transparent
- Compliant with the US and European Food contact regulations
- Compatible with Micronic tubes in 96-well format
- Resists chemical solvents
- Clean room Class 7 production
- RNase/DNase and Pyrogen free

ADDITIONAL OPTIONS

- Sterilization using gamma irradiation or Eto treatment



MANUAL DECAPPERS

We developed a variety of decappers for the manual removal of Push Caps and Screw Caps. Our Manual Decappers secure a quick and easy removal the caps, and are compatible with a variety of brands

SPECIFICATIONS

| | | |
|--|--|---|
| | | Compatible with tubes from 0.50ml to 2.00ml |
| | | Compatible with tubes in 96-well format |
| | | Ethyl Vinyl Acetate |
| | | See drawing on the left |
| | | Resistant to Chemical solvents |
| | | -40°C to +100°C |

ORDERING INFORMATION

More information about product numbers and ordering, see page 69.

ORDERING INFORMATION - CAPS

SCREW CAPS INTERNAL THREAD

| Catalog number | Description | Case Size |
|--|--|--------------|
| Screw Caps Classic for internally threaded tubes in 96-well format* | | |
| MP53120 | Screw Caps for internally threaded tubes in 96-well format - Grey | 1 x 960 caps |
| MP53101 | Screw Caps for internally threaded tubes in 96-well format - White | 1 x 960 caps |
| MP53104 | Screw Caps for internally threaded tubes in 96-well format - Yellow | 1 x 960 caps |
| MP53130 | Screw Caps for internally threaded tubes in 96-well format - Orange | 1 x 960 caps |
| MP53102 | Screw Caps for internally threaded tubes in 96-well format - Red | 1 x 960 caps |
| MP53121 | Screw Caps for internally threaded tubes in 96-well format - Pink | 1 x 960 caps |
| MP53109 | Screw Caps for internally threaded tubes in 96-well format - Purple | 1 x 960 caps |
| MP53103 | Screw Caps for internally threaded tubes in 96-well format - Blue | 1 x 960 caps |
| MP53124 | Screw Caps for internally threaded tubes in 96-well format - Light Blue | 1 x 960 caps |
| MP53123 | Screw Caps for internally threaded tubes in 96-well format - Light Green | 1 x 960 caps |
| MP53105 | Screw Caps for internally threaded tubes in 96-well format - Green | 1 x 960 caps |
| MP53122 | Screw Caps for internally threaded tubes in 96-well format - Black | 1 x 960 caps |
| Screw Caps Ultra for internally threaded tubes in 96-well format | | |
| MP53354 | Screw Caps Ultra for internally threaded tubes in 96-well format - Yellow | 1 x 960 caps |
| MP53352 | Screw Caps Ultra for internally threaded tubes in 96-well format - Red | 1 x 960 caps |
| MP53353 | Screw Caps Ultra for internally threaded tubes in 96-well format - Blue | 1 x 960 caps |
| MP53373 | Screw Caps Ultra for internally threaded tubes in 96-well format - Light Green | 1 x 960 caps |

* Other item numbers are available for Screw Caps Low Profile in 96-format, Screw Caps in 48-format and Screw Caps in 24-format.

SCREW CAPS EXTERNAL THREAD

| Catalog number | Description | Case Size |
|--|--|--------------|
| Screw Caps for externally threaded tubes in 96-well format* | | |
| MP53825 | Screw Caps for externally threaded tubes in 96-well format - Transparent | 1 x 960 caps |
| MP53820 | Screw Caps for externally threaded tubes in 96-well format - Grey | 1 x 960 caps |
| MP53801 | Screw Caps for externally threaded tubes in 96-well format - White | 1 x 960 caps |
| MP53804 | Screw Caps for externally threaded tubes in 96-well format - Yellow | 1 x 960 caps |
| MP53800 | Screw Caps for externally threaded tubes in 96-well format - Orange | 1 x 960 caps |
| MP53802 | Screw Caps for externally threaded tubes in 96-well format - Red | 1 x 960 caps |
| MP53821 | Screw Caps for externally threaded tubes in 96-well format - Pink | 1 x 960 caps |
| MP53809 | Screw Caps for externally threaded tubes in 96-well format - Purple | 1 x 960 caps |
| MP53803 | Screw Caps for externally threaded tubes in 96-well format - Blue | 1 x 960 caps |
| MP53806 | Screw Caps for externally threaded tubes in 96-well format - Light Blue | 1 x 960 caps |
| MP53823 | Screw Caps for externally threaded tubes in 96-well format - Light Green | 1 x 960 caps |
| MP53805 | Screw Caps for externally threaded tubes in 96-well format - Green | 1 x 960 caps |
| MP53822 | Screw Caps for externally threaded tubes in 96-well format - Black | 1 x 960 caps |
| Screw Caps IG for externally threaded tubes in 96-well format** | | |
| MP53870 | Screw Caps for externally threaded tubes in 96-well format - Transparent | 1 x 960 caps |
| MP53871 | Screw Caps for externally threaded tubes in 96-well format - Grey | 1 x 960 caps |
| MP53872 | Screw Caps for externally threaded tubes in 96-well format - White | 1 x 960 caps |

* Other item numbers are available for Screw Caps for externally threaded tubes in 48-format and in 24-format.

** Other colors and item numbers are available for Screw Caps IG for externally threaded tubes.

TPE PUSH CAPS INTERNAL AND EXTERNAL THREAD

from page 66

| Catalog number | Description | Case Size |
|---|--|------------------|
| Pierceable TPE Capclusters for capping 96 individual tubes with External thread* | | |
| MP53970 | Pierceable TPE Capcluster Natural for capping 96 individual tubes with External Thread | 4 x 10 clusters |
| Pierceable TPE Capclusters for capping 96 individual tubes with Internal thread | | |
| MP53065 | Pierceable TPE Capcluster Natural for capping 96 individual tubes with Internal Thread | 4 x 10 clusters |
| MP53068 | Pierceable TPE Capcluster Grey for capping 96 individual tubes with Internal Thread | 4 x 10 clusters |
| MP53073 | Pierceable TPE Capcluster White for capping 96 individual tubes with Internal Thread | 4 x 10 clusters |
| MP53072 | Pierceable TPE Capcluster Yellow for capping 96 individual tubes with Internal Thread | 4 x 10 clusters |
| MP53043 | Pierceable TPE Capcluster Orange for capping 96 individual tubes with Internal Thread | 4 x 10 clusters |
| MP53044 | Pierceable TPE Capcluster Red for capping 96 individual tubes with Internal Thread | 4 x 10 clusters |
| MP53069 | Pierceable TPE Capcluster Pink for capping 96 individual tubes with Internal Thread | 4 x 10 clusters |
| MP53077 | Pierceable TPE Capcluster Purple for capping 96 individual tubes with Internal Thread | 4 x 10 clusters |
| MP53066 | Pierceable TPE Capcluster Blue for capping 96 individual tubes with Internal Thread | 4 x 10 clusters |
| MP53045 | Pierceable TPE Capcluster Light Blue for capping 96 individual tubes with Internal Thread | 4 x 10 clusters |
| MP53067 | Pierceable TPE Capcluster Light Green for capping 96 individual tubes with Internal Thread | 4 x 10 clusters |
| MP53046 | Pierceable TPE Capcluster Green for capping 96 individual tubes with Internal Thread | 4 x 10 clusters |
| MP53071 | Pierceable TPE Capcluster Black for capping 96 individual tubes with Internal Thread | 4 x 10 clusters |
| Pierceable TPE Capmats for capping 96 tubes with External thread* | | |
| MP53990 | Pierceable TPE Capmat Natural for capping 96 tubes with External Thread | 4 x 10 clusters |
| Pierceable TPE Capmats for capping 96 tubes with Internal thread* | | |
| MP53081 | Pierceable TPE Capmat Natural for capping 96 tubes with Internal Thread | 4 x 10 mats |
| MP53084 | Pierceable TPE Capmat Grey for capping 96 tubes with Internal Thread | 4 x 10 mats |
| Pierceable TPE Capbands for capping tubes with Internal thread* | | |
| MP53089 | Pierceable TPE Capband Natural in Capmat format for capping 8 tubes with Internal Thread | 4 x 10 mats |
| MP53098 | Pierceable TPE Capband Natural in Capmat format for capping 12 tubes with Internal Thread | 4 x 10 mats |
| TPE Push Caps RB for capping tubes with Internal thread | | |
| MP53925 | Pierceable TPE Capcluster Natural RB for capping 96 individual tubes with Internal Thread | 24 x 10 clusters |
| MP53930 | Split TPE Capcluster Natural RB for capping 96 individual tubes with Internal Thread | 24 x 10 clusters |
| MP53940 | Split TPE Capband Natural RB for capping 8 tubes with Internal Thread | 24 x 10 clusters |
| Lyo Capclusters for capping 96 individual tubes with Internal Thread (large pack)* | | |
| MP53042 | Lyo Capclusters for capping 96 individual tubes with Internal Thread | 24 x 10 clusters |

* Other colors and item numbers are available for TPE Push Caps

EVA PUSH CAPS INTERNAL THREAD

from page 67

| Catalog number | Description | Case Size |
|----------------------------------|--|----------------|
| EVA Capping Range | | |
| MP82605 | EVA Capband break-off for capping 1 tube | 4 x 1000 bands |
| MP2271B1 | EVA Capband for capping 8 tubes | 24 x 120 bands |
| MP33000 | EVA Capmat for capping 96 tubes | 24 x 10 mats |
| MP33001 | EVA Capmat for capping 96 tubes for Comorack-96 only (polycarbonate) | 24 x 10 mats |
| Manual Decapping Products | | |
| MP54000 | Manual Decapper-1 | 10 pieces |
| MP54001 | Manual Decapper-8 | 10 pieces |
| MP65010 | Manual Decapper-8 (Aluminum) | 1 unit |
| MP54004 | Screw Cap Remover | 10 pieces |
| MP54006 | Low Profile Screw Cap remover | 10 pieces |

Designed
in-house

EQUIPMENT SERVICE & WARRANTY

MICRONIC EQUIPMENT ASSEMBLY CENTER

All Micronic equipment is designed and assembled in-house. We have well-trained personnel on hand to provide service on equipment. These services range from annual maintenance, installation support and customization, to repairs.

At Micronic, high quality and a high level of service are of paramount importance. By assembling equipment in-house, Micronic's equipment assembly personnel and engineers have the first-hand knowledge and skills necessary to perform services quickly and smoothly.



ASSEMBLED IN
THE NETHERLANDS
AND THE UNITED STATES

CONTINUOUS INNOVATION IN
SAMPLE HANDLING EQUIPMENT



SAMPLE HANDLING EQUIPMENT

TO ENABLE FAST, ACCURATE AND EFFICIENT SAMPLE IDENTIFICATION, MICRONIC OFFERS AN EXTENSIVE CODE READER RANGE WHICH CONSISTS OF SINGLE TUBE SCANNERS AND WHOLE RACK SCANNERS. THE SCANNERS CAN READ CODED 24-, 48- AND 96-WELL FORMAT TUBES. MICRONIC OFFERS ALSO A WIDE RANGE OF MANUAL AND AUTOMATED CAPPING AND DECAPPING EQUIPMENT. THESE DEVICES ENABLE SIGNIFICANT PRODUCTIVITY GAINS TO BE MADE IN THE OPENING AND CLOSING OF MULTIPLE SAMPLE STORAGE TUBES. IN ADDITION, SEVERAL SAMPLE MANAGEMENT SOLUTIONS ARE AVAILABLE THAT SUIT ALL THROUGHPUT AND BUDGET NEEDS.

MICRONIC WARRANTY ON EQUIPMENT

The Micronic products are designed and manufactured to the highest quality standards to deliver outstanding performance, ease of use and ease of installation. In case a user encounters any difficulties while installing or using the product, please let us know, we're glad to help.

Micronic warrants that its products are of satisfactory quality and conform to applicable EU legislation. If the product turns out to be defective, Micronic will arrange to repair or replace the product during the warranty period. The warranty period starts on the day of Micronic's invoice of the original product and ends after five (5) years with respect to Micronic Code Reading equipment and after one (1) year with respect to any other equipment.

The Micronic Warranty does not affect consumer statutory rights under applicable local laws, nor consumer rights against the seller arising from the purchase transaction. For more information about Micronic Warranty on equipment or about its applicability, please request the 'Micronic Warranty Statement'.

Service needed?

To obtain service please contact Micronic America / Asia / Europe or a local Micronic supplier.

CODE READER OVERVIEW AND SPECIFICATIONS

| | TUBE READER DT300 | TUBE READER DT510 | RACK READER DR505 | RACK READER DR700 | RACK READER DR710 |
|----------------------------------|--|---|--|--|--|
| |  Page 74 |  Page 75 |  Page 76 |  Page 77 |  Page 77 |
| Basic Features | | | | | |
| Imaging technology | CIS | CMOS | CIS | CCD | CCD |
| Decode time | < 1 second | < 1 second | 1 second | 1 second | 1 second |
| Scan time | < 1 second | < 1 second | 5 seconds | 7 seconds | 7 seconds |
| Dimensions (mm) | 150 x 40 x 40 | 90 x 80 x 95.4 | 265 x 186 x 70 | 309 x 230 x 87 | 309 x 230 x 87 |
| Decoding software | n/a | n/a | Micronic Code Reader Software | Micronic Code Reader Software | Micronic Code Reader Software |
| Software compatibility | | | | | |
| Windows 10 and 7* | Yes | Yes | Yes | Yes | Yes |
| Connection | USB x1 | USB x1 | USB x2 | USB x2 | USB x2 |
| Output options | | | | | |
| Output text file | Yes | Yes | Yes | Yes | Yes |
| Configure output file | Yes | Yes | Yes | Yes | Yes |
| Export to Excel | Yes | Yes | Yes | Yes | Yes |
| ODBC Database push | n/a | n/a | Yes | Yes | Yes |
| User account independent | Yes | n/a | Yes | Yes | Yes |
| Network ready | n/a | n/a | Yes | Yes | Yes |
| TCP/IP Remote control | n/a | n/a | Yes | Yes | Yes |
| Tube / rack compatibility | | | | | |
| 24-well format tube / rack | Yes | Yes | Yes | Yes | Yes |
| 48-well format tube / rack | Yes | Yes | Yes | Yes | Yes |
| 96-well format tube / rack | Yes | Yes | Yes | Yes | Yes |
| Read all 2D Data-Matrix codes | Yes | Yes | Yes | Yes | Yes |
| Other features | | | | | |
| No tube, no code feature | n/a | n/a | Yes | Yes | Yes |
| Single tube reading | Yes | Yes | Yes | Yes | Yes |
| Rack templates | n/a | n/a | Yes | Yes | Yes |
| Anti-frost system | n/a | No | Yes | Yes | Yes |
| Robot gripper friendly | n/a | No | Yes | Yes | Yes |
| Linear barcode reading | Yes, 1D tube and rack barcodes | Yes, 1D tube and rack barcodes | Optional, 1D rack barcodes | Optional, 1D rack barcodes | Optional, 1D rack barcodes |
| Rack bottom ID (2D code) reading | n/a | Yes | Yes | Yes | Yes |

*Please contact us for compatibility with other Windows versions.




MICRONIC
INNOVATION IN SAMPLE STORAGE

BROCHURE

- SERVING SCIENTISTS SINCE 1964
- INDUSTRY LEADING 5-YEAR WARRANTY ON CODE READERS
- ASSEMBLY IN THE NETHERLANDS AND THE UNITED STATES

MICRONIC CODE READER RANGE
Single Tube and Whole Rack Scanners

The Micronic Tube Reader DT300 enables users to reliably read 1D and 2D tube and rack codes almost instantly. This small lightweight device is practical to carry around, easy-to-use and can read any 1D barcode or 2D Data-Matrix code on 24-, 48- and 96-well format tubes and racks. The DT300 offers convenient, high performance scanning anywhere you might need it.

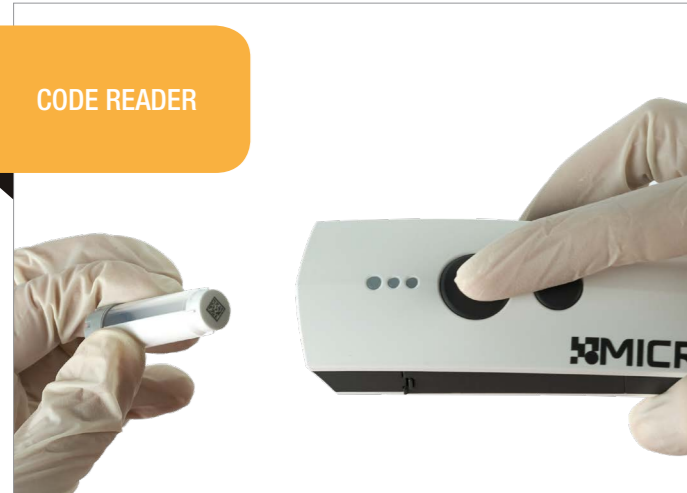

SUPERIOR FEATURES

- Instantly reads 2D and 1D codes
- Compatible with iOS, Android and Windows and Bluetooth dongles attached to any laptop
- Storage of approximately 15,000 2D scans and 17,000 1D scans
- Personalized configuration with time, date, user, project name, and other useful data
- Reads directly into a LIMS system, Excel spreadsheet or notepad
- Long-lasting battery through 10-12 hours of continuous usage



REQUEST DEMO

Experience what the DT300 has to offer by requesting a demo. E-mail us at sales@micronic.com or call +1 484 480 3372 (Micronic America) / +65 85 028 878 (Micronic Asia) / +31 (0)320 277 099 (Micronic Europe).

SPECIFICATIONS

| | | |
|--|--|--|
| | | Compatible with tubes from 0.30ml to 6.00ml |
| | | Compatible with Micronic tubes and tubes from other brands |
| | | Scan time is < 1 second |
| | | 150 x 40 x 40 mm / 5.91 x 1.57 x 1.57 inch |
| | | 80 grams / 0.17 lbs |
| | | Via USB cable |

ORDERING INFORMATION

More information about product numbers and ordering, see page 87.

The Micronic Tube Reader DT510 is high-speed code reading at its best. The DT510 is the most versatile single tube reader in the market - it reads 1D rack barcodes and 1D barcoded and 2D Data-Matrix coded tubes in less than a second. It is a plug-and-play reader that requires no software installation and is very easy to use.



SUPERIOR FEATURES

- High-end sophisticated decoding engine
- Fully automated scan and decode feature
- No software installation required
- Scans and decodes 1D & 2D rack codes and 1D & 2D Data-Matrix coded tubes in less than a second
- Compatible with tubes in 96-, 48- and 24-well format
- Small footprint
- Keyboard wedge functionality
- Correct code reads are indicated with a bleep
- USB connection (no power supply needed)
- Compatible with other suppliers of 1D and 2D barcoded tubes



SPECIFICATIONS

| | | |
|--|--|--|
| | | Compatible with tubes from 0.30ml to 6.00ml |
| | | Compatible with Micronic tubes and tubes from other brands |
| | | Scan time is < 1 second |
| | | 90 x 80 x 95.4 mm / 3.54 x 3.15 x 3.76 inch |
| | | 321 grams / 0.7 lbs |
| | | 5VDC 0.3A |

ORDERING INFORMATION

More information about product numbers and ordering, see page 87.

The Micronic Rack Reader DR505 reads whole racks filled with 2D Data-Matrix coded tubes quickly, cost effective and with limited lab space. The compact reader is built for high throughput and automated systems. It enables a reliable interface with your sample organization.




SUPERIOR FEATURES






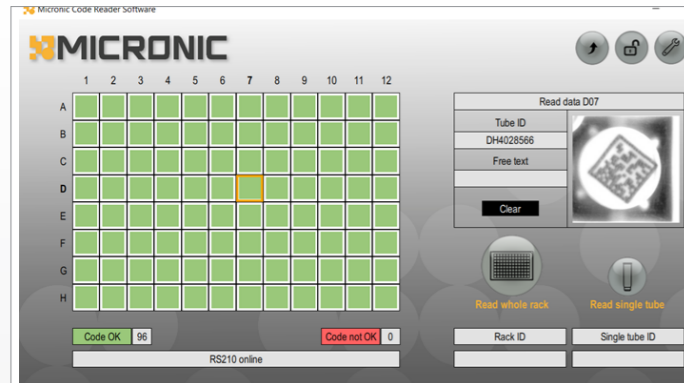

- Reads a whole rack (in 96-, 48- or 24-well format) in less than 5 seconds and a single 96-well format tube in less than 2 seconds
- Anti-frost system which minimizes condensation of frozen samples on the scanner plate
- Scans and decodes 2D Data-Matrix coded tubes
- Small bench top solution
- Easy-to-use Code Reader Software (latest version)
- Outputs CSV-files (text)
- Compatible with most brands of 2D Data-Matrix coded tubes
- 5-year warranty
- Integration with LIMS systems

ADDITIONAL OPTIONS

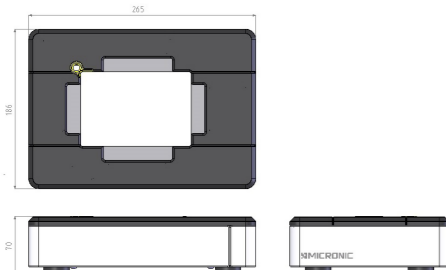- Integrated Micronic Side Barcode Reader for DR500 line



SPECIFICATIONS

| | | |
|---|---|--|
|  |  | Compatible with tubes from 0.30ml to 6.00ml |
| |  | Compatible with Micronic tubes and tubes from other brands |
| |  | 5 seconds (whole rack) / 2 seconds (single tube) |
| |  | 265 x 186 x 70 mm / 10.4 x 7.3 x 2.8 inch |
| |  | 3.4 kg / 7.5 lbs |
| |  | 115 - 230 V |

ORDERING INFORMATION

More information about product numbers and ordering, see page 87.

The high-end Micronic Rack Readers DR700 and DR710 read whole racks filled with 2D Data-Matrix coded tubes. The readers feature a CCD image sensor ensuring the highest accuracy 2D code reading. The DR710 has a uniquely designed active anti-frost system, which prevents the condensation of the scanning plate so that codes from frozen samples can be read.



SUPERIOR FEATURES








- CCD image sensor for high resolution image quality
- Active anti-frost system which prevents condensation of frozen samples on the scanner plate (DR710 model)
- Reads racks filled with tubes (in 96-, 48- or 24-well format) in 7 seconds
- Reads a single tube (in 96-, 48- or 24-well format) in less than 2 seconds
- Scans and decodes 2D Data-Matrix codes
- Small bench top solution
- Easy-to-use Code Reader Software (latest version)
- Outputs CSV-files (text)
- Compatible with most brands of 2D Data-Matrix coded tubes
- 5-year warranty

ADDITIONAL OPTIONS

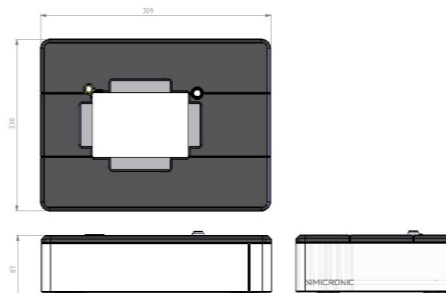- Integrated Micronic Side Barcode Reader for DR700 line



SPECIFICATIONS

| | | |
|---|---|--|
|  |  | Compatible with tubes from 0.30ml to 6.00ml |
| |  | Compatible with Micronic tubes and tubes from other brands |
| |  | 7 seconds (whole rack) / 2 seconds (single tube) |
| |  | 309 x 230 x 87 mm / 12.2 x 9.1 x 3.4 inch |
| |  | 5.1 kg / 11.2 lbs (estimated) |
| |  | 115 - 230 V |

ORDERING INFORMATION

More information about product numbers and ordering, see page 87.

The Manual Micronic Push Cap Capper CP400 offers a cost effective way of applying uniform pressure to Push Caps in an entire 96 tubes rack in only a single action. Built to meet the varying needs of different users, the Push Cap Capper CP400 can seal full or partially full racks with the complete range of Micronic TPE or EVA style capmats, capbands and capclusters.



SUPERIOR FEATURES

- Caps tubes (up to 1.40 ml) in all ANSI/SLAS format racks up to maximum height of 48.0mm
- Caps an entire 96-format rack in one single action with minimum force required
- Compact design and easy to use - no installation required
- Safe operation and even capping guaranteed
- Integrity of sample ensured
- Minimizes cross-contamination
- Works on full and partially-full racks
- Height adapters match customer's specifications
- Cost effective
- Robust and durable material
- Compatible with all Micronic push caps (TPE or EVA style capmats, capbands and capclusters)
- Compatible with other brands of tubes, caps, and racks

ADDITIONAL OPTIONS

- Adapters available for tubes from other brands



SPECIFICATIONS

| | | |
|--|--|--|
| | | Compatible with tubes from 0.30ml to 1.40ml |
| | | Compatible with Micronic tubes and tubes from other brands |
| | | +/- 5 seconds (whole rack) |
| | | 262 x 185 x 164 mm / 10.4 x 7.3 x 2.8 inch |
| | | 7 Kg / 15 lbs |
| | | n/a |

ORDERING INFORMATION

More information about product numbers and ordering, see page 87.

The Push Cap Cappers CP600 and CP610 can cap Micronic Push Caps on 96-format tubes in rack. For capping, simply place the rack filled with tubes in the sample holder and place a capmat, capband or capcluster on the tubes. A controlled pressure system on the Electric Cappers presses all the caps down evenly to ensure uniform, high integrity storage of your samples in capped tubes.



SUPERIOR FEATURES

- No adapters required
- Caps an entire rack of 96 tubes in an average of 7 seconds (depends on the tube height)
- Operates without any buttons
- Compact and robust design
- Controlled pressure system for an evenly capping result
- Works on full and partially full racks (minimum amount of 10 tubes)
- Compatible with the Micronic range of 96-format tubes and racks
- Works with tube sizes up to 1.40ml (CP600) or with tube sizes up to 2.00ml (CP610)
- Also compatible with other brands of tubes and caps



"The capping equipment and TPE push caps have been super efficient and easy to use, I use them as often as I can. The reliability of the equipment has been great with continuous usage."

- Bristol Myers Squibb

SPECIFICATIONS

| | | |
|--|--|---|
| | | Compatible with tubes up to 2.00ml (CP610) / up to 1.40ml (CP600) |
| | | Compatible with Micronic tubes and tubes from other brands |
| | | 7 seconds (whole rack) |
| | | 312 x 209 x 286 mm / 12.3 x 8.2 x 11.3 inch (CP600) |
| | | 14 kg / 31 lbs (CP600) |
| | | 115 - 230 V |

ORDERING INFORMATION

More information about product numbers and ordering, see page 87.

With the ever-growing requirement for higher sample throughput in many laboratories the Micronic Push Cap Decapper CP620 enables significant productivity gains to be made in the opening of multiple sample storage tubes. The internal cap grippers are developed in such a way that cross-contamination is minimized.



SUPERIOR FEATURES

- No adapters required
- Decaps 96 tubes in 8 seconds
- Simple and safe operation
- Friendly user interface for an easy selection of tube size
- CE-approved and conforms to UL
- Operates stand-alone or in a fully automated environment
- Minimum cross-contamination
- Remote control via RS232
- Compatible with the Micronic range of 96-format internal thread tubes (up to 1.40ml)
- Also compatible with other brands of tubes and caps



NEED ASSISTANCE?

We're committed to providing fast, customer focused support on all your technical enquiries. E-mail us at helpdesk@micronic.com or call +1 484 480 3372 (Micronic America) / +65 85 028 878 (Micronic Asia) / +31 (0)320 277 099 (Micronic Europe), we'll be right there with you.


SPECIFICATIONS

| | | |
|--|--|--|
| | | Compatible with tubes up to 1.40ml |
| | | Compatible with Micronic tubes and tubes from other brands |
| | | 10-15 seconds (whole rack) |
| | | 531 x 258 x 366 mm / 20.9 x 10.2 x 14.4 inch |
| | | 25 kg / 55 lbs |
| | | 115 - 230 V |

ORDERING INFORMATION

More information about product numbers and ordering, see page 88.

The Micronic Screw Cap Recapper CS100 line is developed to improve the manual (de)capping of tubes. The manual recappers are available for 96- (CS100 model), 48- (CS110 model) or 24- (CS120 model) well format tubes. The cost-effective tools are set to a precise torque value and stop turning when a perfect seal is established, ensuring an optimal sealing performance.

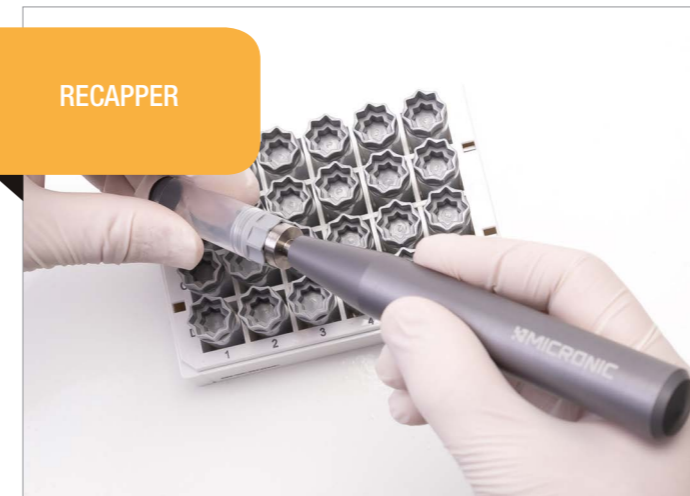

SUPERIOR FEATURES

- Quick and easy to use
- Each screw cap is reproducibly tightened securely and uniformly
- Minimizes contamination
- Ergonomic lightweight solution for left and right handed use
- Set to a precise torque value and stops turning when the perfect seal is established
- No need for battery recharging
- Made from stainless steel
- Models available for Micronic 96-, 48- and 24-well format Screw Caps, as well as caps from other brands



SPECIFICATIONS

| | | |
|--|--|--|
| | | Compatible with tubes from 0.30ml to 6.00ml |
| | | Compatible with Micronic tubes and tubes from other brands |
| | | 130 x 19 mm / 5.12 x 0.75 inch |
| | | 0.095 kg / 0.21 lbs |

ORDERING INFORMATION

More information about product numbers and ordering, see page 88.

The Micronic Screw Cap Recapper CS500 line is developed in conjunction with several leading laboratories. The handheld device offers users the ability to cap, decap or recap a row of 8 (CS500 model), 6 (CS510 model) or 4 (CS520 model) tubes in less than 5 seconds - 10 times faster than by hand! Thereby, minimizing sample evaporation and significantly improving productivity.



SUPERIOR FEATURES

- Highly durable solution to cap, decap or recap a row of 8,6 or 4 tubes at once
- Each screw cap is reproducibly tightened securely and uniformly
- Caps a row of 8 tubes in 5 seconds and a whole rack of 96 tubes in only 1 minute - 10 times faster than by hand!
- Minimizes contamination
- Easy positioning of screw caps above the tubes
- Ergonomic lightweight solution for left and right handed use
- Caps can be easily discarded using the large ejection handle
- Very small footprint for bench top use
- No need for battery recharging
- Models available for Micronic 96-, 48- and 24-well format Screw Caps, as well as caps from other brands

ADDITIONAL OPTIONS

- Docking station for CS500 line



SPECIFICATIONS

| | | |
|--|--|--|
| | | Compatible with tubes from 0.30ml to 6.00ml |
| | | Compatible with Micronic tubes and tubes from other brands |
| | | <5 seconds (per row) |
| | | 178 x 161 x 710 mm / 7.01 x 6.34 x 27.95 inch |
| | | 5.5 kg / 12.0 lbs |
| | | 115 - 230 V |

ORDERING INFORMATION

More information about product numbers and ordering, see page 88.

The Micronic Docking Station for CS500 semi-automated recapping systems is a compact, benchtop unit designed for efficient tube recapping in labs with medium throughput. The combination can be used to recap 96-, 48- and 24-well format sample storage tubes. The Docking Station offers the consistency of an automated recapping system, but at a much lower cost.



SUPERIOR FEATURES

- Consistent capping and decapping action is assured
- User friendly
- Improves ergonomics
- Small footprint for bench top use
- Caps, decaps or recaps a row of 8-6-4 tubes in 5 seconds
- Each screw cap is reproducibly tightened securely and uniformly
- Compatible with the complete line of Micronic CS500 Recappers
- Compatible with ANSI / SLAS standard format storage racks from Micronic and other brands



COMBI-DEAL

Benefit from an attractive combi-deal when purchasing the Docking Station and Recapper together. Inquire now! E-mail us at sales@micronic.com or call +1 484 480 3372 (Micronic America) / +65 85 028 878 (Micronic Asia) / +31 (0)320 277 099 (Micronic Europe), we'll be right there with you.

SPECIFICATIONS

| | | |
|--|--|---|
| | | Compatible with tubes from 0.30ml to 6.00ml |
| | | Compatible with ANSI / SLAS standard format storage racks |
| | | <5 seconds (per row) |
| | | 420.5 x 370 x 503.5 mm / 16.56 x 14.57 x 19.82 inch |
| | | 9 kg / 19.8 lbs |

ORDERING INFORMATION


More information about product numbers and ordering, see page 88.

The Micronic Screw Cap Recapper CS700 line offers users the ability to cap, decap or recap 96 (model CS700), 48 (model CS710) or 24 (model CS720) tubes with Screw Caps in a single action. Designed to improve efficiency, the Micronic Screw Cap Recapper can be used as a stand-alone instrument or be simply integrated into a fully automated laboratory environment.



SUPERIOR FEATURES

- Extremely durable design for reliable operation
- Caps, decaps or recaps 96 Screw Cap tubes at once in less than 30 seconds
- Compatible with the complete Micronic range of 96-well format tubes (up to 2.00ml)
- Precise torque value for optimal sealing quality
- No adapters required
- Easy-to-use graphical touch screen interface
- Operates stand-alone or in a fully automated environment
- Sensor cap detection that enhances process safety for automation (optional)
- Ethernet connection provides many possibilities for different communication protocols
- Models available for Micronic 96-, 48- and 24-well format Screw Caps, as well as caps from other brands

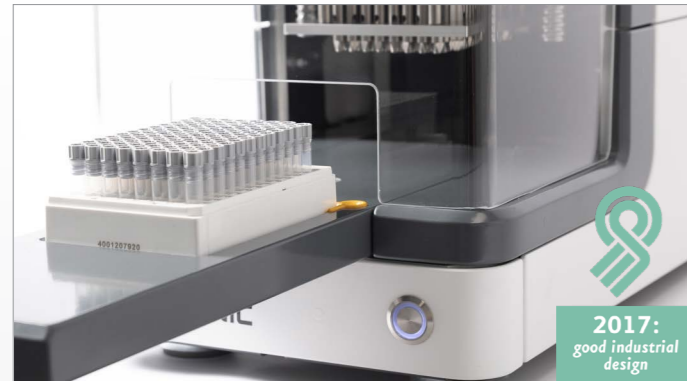

"It's a very smoothly running, easy to use and silent device. I also think it is a great benefit that the Screw Cap Recapper is contamination-free."

- Radboud UMC

SPECIFICATIONS

| | | |
|--|--|--|
| | | Compatible with tubes from 0.30ml to 6.00ml |
| | | Compatible with Micronic tubes and tubes from other brands |
| | | 30 seconds (whole rack) |
| | | 590 x 251 x 460 mm / 23.2 x 9.9 x 18.1 inch |
| | | 38 kg / 84 lbs |
| | | 115 - 230 V |

ORDERING INFORMATION

More information about product numbers and ordering, see page 89.

The Tube Handler HT500 is an affordable sample management system that sorts your valuable liquid and solid samples. The device can be used to pick specific specimens from a large group of samples. The HT500 has a total rack capacity of 20 racks, specify at least on target rack and one source rack per run.



SUPERIOR FEATURES

- Total rack capacity: 20
- Fast pick & place operation of approx. 400 tubes per hour
- 2D code reading and 1D code reading (source and target rack)
- Easy to navigate and intuitive interface
- Semi-automatic linking between rack ID and rack position
- Loading CSV-file pick list from USB thumb drive
- Output CSV-file of rack content
- Separate safety stop ensures sample/product protection
- Compatible with Micronic racks (ANSI/SLAS standard)
- Compatible with 0.30ml, 0.40ml, 0.50ml, 0.75ml, 0.80ml, 1.10ml and 1.40ml Micronic tubes



ADDITIONAL OPTIONS

- Contact us for the compatibility with other brands

SPECIFICATIONS

| | | |
|--|--|--|
| | | Compatible with tubes from 0.30ml to 1.40ml |
| | | Compatible with Micronic tubes in 96-well format |
| | | Sort tubes from a total capacity of 20 racks |
| | | 666 x 1153 x 604 mm / 24.88 x 23.78 x 23.74 inch |
| | | 60 kg / 132 lbs |
| | | 230 V |

ORDERING INFORMATION

More information about product numbers and ordering, see page 89.

The space efficient Tube Handler HT700 provides an easy, automated and reliable solution to sort sample storage tubes. The device can be used to pick specific specimens from a large group of samples. Automated sample identification and collection reduces human errors and reliefs researchers from repetitive manual labour.



SUPERIOR FEATURES








- Minimal footprint increase with doubling of capacity
- Source rack capacity: 24
- Target rack capacity: 1-3
- Fast pick & place operation of approx. 500 tubes per hour
- Picking and 2D code reading a high number of samples with a space efficient design
- 2D code reading and 1D code reading (source and target rack)
- Easy to navigate and intuitive interface
- Automatic linking between rack ID and rack position
- Loading CSV-file pick list from USB thumb drive
- Grouping field in pick list for sample clustering in target rack
- Output CSV-file of rack content
- Enclosed system for operator safety
- Separate safety stop ensures sample/product protection
- Compatible with Micronic racks (ANSI/SLAS standard)
- Compatible with 0.30ml, 0.40ml, 0.50ml, 0.75ml, 0.80ml, 1.10ml and 1.40ml Micronic tubes

ADDITIONAL OPTIONS

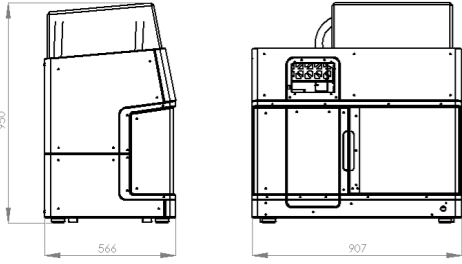- Contact us for the compatibility with other brands



SPECIFICATIONS

| | | |
|---|---|---|
|  |  | Compatible with tubes from 0.30ml to 1.40ml |
| |  | Compatible with Micronic tubes in 96-well format |
| |  | Sort tubes from 24 source racks into 1-3 target racks |
| |  | 566 x 907 x 950 mm / 223 x 358 x 374 inch |
| |  | 120 kg / 265 lbs |
| |  | 230 V |

ORDERING INFORMATION

More information about product numbers and ordering, see page 89.

ORDERING INFORMATION - SAMPLE HANDLING EQUIPMENT & SOFTWARE

TUBE READERS

from page 74-75

| Catalog number | Description | Case Size |
|----------------------------|----------------------------|-----------|
| Single Tube Readers | | |
| MP55127 | Micronic Tube Reader DT300 | 1 unit |
| MP55128 | Micronic Tube Reader DT510 | 1 unit |

RACK READER DR505

from page 76

| Catalog number | Description | Case Size |
|----------------------------|---|-----------|
| Single Rack Readers | | |
| MP35221 | Micronic Rack Reader DR505 | 1 unit |
| MP35240 | Micronic Side Barcode Reader for DR500 line | 1 unit |
| MP35261 | Micronic Rack Reader DR505 & Side Barcode Reader Pack | 1 unit |

RACK READER DR700 RANGE

from page 77

| Catalog number | Description | Case Size |
|----------------------------|---|-----------|
| Single Rack Readers | | |
| MP35230 | Micronic Rack Reader DR700 | 1 unit |
| MP35232 | Micronic Rack Reader DR710 | 1 unit |
| MP35250 | Micronic Side Barcode Reader for DR700 line | 1 unit |
| MP35262 | Micronic Rack Reader DR700 & Side Barcode Reader Pack | 1 unit |
| MP35263 | Micronic Rack Reader DR710 & Side Barcode Reader Pack | 1 unit |

PUSH CAP CAPPER CP400

from page 78

| Catalog number | Description | Case Size |
|-----------------------|---|-----------|
| Manual Cappers | | |
| MP35050 | Micronic Push Cap Capper CP400 (including 1 adapter of your choice) | 1 unit |
| MP35054 | Micronic 0.50ml Adapter for Push Cap Capper CP400 | 1 unit |
| MP35053 | Micronic 0.75ml Adapter for Push Cap Capper CP400 | 1 unit |
| MP35052 | Micronic 1.10ml Adapter for Push Cap Capper CP400 | 1 unit |
| MP35051 | Micronic 1.40ml Adapter for Push Cap Capper CP400 | 1 unit |

PUSH CAP CAPPER CP600/CP610

from page 79

| Catalog number | Description | Case Size |
|-------------------------|---|-----------|
| Push Cap Cappers | | |
| MP35020 | Micronic Push Cap Capper CP600 for tube sizes up to 1.40ml in 96-format | 1 unit |
| MP35021 | Micronic Push Cap Capper CP610 for tube sizes up to 2.00ml in 96-format | 1 unit |



SERVING SCIENTISTS
SINCE 1984

CERTIFIED CLASS 7 CLEAN ROOM
MANUFACTURING

SAMPLE STORAGE STARTER PACKS

Everything
you need

EVERYTHING YOU NEED TO GET STARTED WITH USING 2D CODED TUBES TO TRACK AND STORE YOUR VALUABLE SAMPLES. MICRONIC OFFERS A FULLY TRACEABLE STORAGE SOLUTION AT A LOW COST. BUYING THE STARTER PACK CAN SAVE YOU UP TO 20% COMPARED TO BUYING THE PRODUCTS INDIVIDUALLY. SINCE THE AMOUNT OF STORED SAMPLES DIFFERS LABORATORY BY LABORATORY, MICRONIC OFFERS STARTER PACKS IN A RANGE OF SIZES: BASIC, STANDARD, ADVANCED, PREMIUM AND PREMIUM+.

SAMPLE STORAGE WITH MICRONIC



COLLECT SAMPLES

Micronic products are essential during your sample storage process. When a liquid sample is obtained, it is often temporarily stored in a large tube, like vacuum containers. For long-term storage, the sample is transferred from the larger tube into several Micronic sample storage tubes.



CLOSE SAMPLE STORAGE TUBES

To securely close your sample storage tubes for long-term storage in -80°C, the TPE Cap-clusters from Micronic are an ideal and economical solution. Place a TPE Capcluster on the tubes in a 96-well format rack and the Micronic Push Cap Capper CP600 closes all 96 tubes in only 7 seconds! Remove the black foil to leave the tubes individually capped.



SCAN TUBE CODES

The Micronic Rack Reader reads the 2D Data-Matrix codes of sample storage tubes and the 1D barcodes of sample storage racks in just 5 seconds. The intuitive user interface of the Micronic Code Reader Software is designed for maximum functionality and enables full sample traceability.



ADD SAMPLES TO YOUR DATABASE

Micronic Rack Readers are the link between your Sample Management Software and every individual sample. To make sure that the Micronic sample storage solution works with Sample Management Software we have extensively tested our readers with a wide range of Software Systems.



PLACE SAMPLES IN ALLOCATED FREEZER

After the sample data has been saved, the rack cover can be placed on the rack. The samples are now ready to be stored at the allocated freezer position.

STARTER PACKS

SAMPLE STORAGE STARTER PACK OVERVIEW

| | SAMPLE STORAGE STARTER PACK BASIC | SAMPLE STORAGE STARTER PACK STANDARD | SAMPLE STORAGE STARTER PACK ADVANCED | SAMPLE STORAGE STARTER PACK PREMIUM | SAMPLE STORAGE STARTER PACK PREMIUM+ |
|--|--|---|---|--|---|
| Starter packs with Screw Caps | | | | | Not Available |
| 10 racks filled with 2D coded tubes (internal and external thread) | Yes (0.75ml / 1.40ml) | Yes (0.75ml / 1.40ml) | No | No | n/a |
| 40 racks filled with 2D coded tubes (internal and external thread) | No | No | Yes (1.00/3.00/3.50/4.00ml) | Yes (0.75ml / 1.40ml) | n/a |
| Screw Caps in Grey to cap 10 racks with tubes | Yes | Yes | No | No | n/a |
| Screw Caps in Grey to cap 40 racks with tubes | No | No | Yes | Yes | n/a |
| Micronic Tube Reader DT510 | Yes | No | No | No | n/a |
| Micronic Rack Reader DR505 | No | Yes | Yes | Yes | n/a |
| Micronic Side Barcode Reader for DR500 line | No | No | Yes | No | n/a |
| Micronic Tube Reader DT300 | Yes | No | No | No | n/a |
| Micronic Screw Cap Recapper CS500 | No | No | No | Yes | n/a |
| Manual decapping tools | Yes | Yes | No | Yes | n/a |
| Starter packs with Push Caps | | | Not Available | | |
| 10 racks filled with 2D coded tubes (internal thread) | Yes (0.75ml / 1.40ml) | Yes (0.75ml / 1.40ml) | n/a | No | No |
| 40 racks filled with 2D Data-Matrix (internal thread) | No | No | n/a | Yes (0.75ml / 1.40ml) | Yes (0.75ml / 1.40ml) |
| TPE Caps in Grey to cap 10 racks with tubes | Yes | Yes | n/a | No | No |
| TPE Caps in Grey to cap 40 racks with tubes | No | No | n/a | Yes | Yes |
| Micronic Tube Reader DT510 | Yes | No | n/a | No | No |
| Micronic Rack Reader DR700 | No | No | n/a | No | Yes |
| Micronic Rack Reader DR505 | No | Yes | n/a | Yes | No |
| Micronic Side Barcode Reader for DR700 line | No | No | n/a | No | Yes |
| Micronic Tube Reader DT300 | Yes | No | n/a | No | No |
| Micronic Push Cap Capper CP400 | No | No | n/a | Yes | No |
| Micronic Push Cap Capper CP600 | No | No | n/a | No | Yes |
| Manual decapper-1 | Yes | Yes | n/a | Yes | Yes |
| Manual decapper-8 | Yes | Yes | n/a | Yes | Yes |

Tissue storage pack

INNOVATION AT MICRONIC



Industry leaders since 1984, Micronic serves scientists with sample storage and automation solutions. At Micronic, advancing scientific research to enhance quality of life is our corporate goal and to achieve this we make innovation and quality a priority. We engage in various innovative processes within our organization, like product development and customer focused lean manufacturing systems, in order to create new and improved products, solutions and services that provide radically better customer experiences. Our premium labware and precision engineered equipment is developed and manufactured in-house. This core competence gives us the opportunity to offer top-of-the-line products, delivering exceptional quality, reliability and durability. Above all, we stand behind our solutions and are confident that you will too.

ACTIVITIES

RESEARCH & DEVELOPMENT

The primary goal of our Research & Development (R&D) department is to develop new products, and to optimize existing ones. Another fundamental task of the team is to discover and create new knowledge about scientific and technological topics for the purpose of uncovering and enabling development of valuable new products, processes and services. At Micronic, we utilize a technology driven system in which R&D is directed towards developing those products, indicated by means of market research, that will meet an unmet need. By developing and manufacturing our products in-house, we have all first-hand high-tech manufacturing and R&D knowledge necessary to quickly adapt to developments in the market.

CUSTOM PROJECTS

To help you find the best solutions for your sample storage needs, we are open to OEM developments and customer specific requests. If you have any suggestions about new or improved sample storage products, please contact us at sales@micronic.com. Micronic can offer you highly/academically trained personnel, knowledge in materials and injection mold development, technical and electrical knowledge, and a large network of suppliers. Our R&D department has also access to the following facilities: 3D CAD engineering, printing/plotting machines, a test facility with injection molding machines, a technical department for prototyping and modifications, and a laboratory testing facility.

QUALITY COMMITMENT

Manufactured following industry leading strict tolerances, the precision and tube-to-tube consistency of Micronic labware optimizes the compatibility with automated sample handling equipment. The materials and auxiliary agents used to manufacture our sample storage labware are FDA compliant and are selected on quality and the highest purity levels. In addition to our high manufacturing and material standards, we also engage in extensive product testing. When new or improved products are developed, they undergo several tests during the process to fine-tune the product and to create the most optimal solution for specific customer applications. A wide range of



tests are conducted at our laboratory testing facility, examples are: freeze-thaw cycles, protein binding tests, durability tests, evaporation/sealing tests, ultra-low temperature resistance tests, code readability tests and code scratch tests. Please be aware that Micronic takes great care to test its products in several usage conditions. However, as it is not possible to cover all applications, we highly recommend customers to test the products also within their own application before usage. If you need any assistance in performing these application tests, our technical support team is very happy to help.

DEDICATION TO THE ENVIRONMENT

Our corporate goal extends to our commitment to safeguarding the natural environment for future generations. We have implemented environmental activities into our daily operations, and continuously work toward goals to increase efficiencies and reduce waste. In addition, Micronic constantly strives to utilize and develop solutions that minimize the impact on the environment – whether it is an energy efficient injection molding machine, plastic recycling, waste sorting, eco-friendly packaging or high-density products that promote efficiency at our customers' sites. Micronic takes further responsibility with respect to the environment by its internal environmental management system which is certified by Lloyd's Register Quality Assurance (ISO 14001). We are committed to comply with and where possible, considerably exceed all applicable laws and regulations related to environmental protection.



The ACT label is the first environmental impact factor label for laboratory products, developed by the non-profit organization My Green Lab. Micronic participates in this great program so that purchasers can make informed decisions about a product's environmental impact and to promote sustainability in the laboratory. Recently we have recertified our existing labels for the 1.40ml internally threaded tube and 4.00ml externally threaded tube. Both tube types now have an Environmental Impact Factor (EIF) of 32.8, which is a 12.8-point reduction compared to the last labels. With some steps we have taken at our manufacturing facility in reducing our use of water, energy and waste and making our packaging more sustainable we have improved our EIF-score. We are incredibly proud for our point reduction and it motivates us to further improve the score for the future.

| Environmental Impact Scale | |
|-------------------------------------|-------------|
| Decreasing Environmental Impact | |
| Manufacturing | |
| Manufacturing Impact Reduction | 1.0 |
| Renewable Energy Use | No |
| Responsible Chemical Management | 1.0 |
| Shipping Impact | 1.0 |
| Product Content | 10.0 |
| Packaging Content | 1.0 |
| User Impact | |
| Energy Consumption (kWh/day) | N/A |
| Water Consumption (liters/day) | N/A |
| Product Lifetime | 6.0 |
| End of Life | |
| Packaging | 4.0 |
| Product | 8.8 |
| Environmental Impact Factor: | 32.8 |
| Label Valid Through: | June 2023 |

| Environmental Impact Scale | |
|-------------------------------------|-------------|
| Decreasing Environmental Impact | |
| Manufacturing | |
| Manufacturing Impact Reduction | 1.0 |
| Renewable Energy Use | No |
| Responsible Chemical Management | 1.0 |
| Shipping Impact | 1.0 |
| Product Content | 10.0 |
| Packaging Content | 1.0 |
| User Impact | |
| Energy Consumption (kWh/day) | N/A |
| Water Consumption (liters/day) | N/A |
| Product Lifetime | 6.0 |
| End of Life | |
| Packaging | 4.0 |
| Product | 8.8 |
| Environmental Impact Factor: | 32.8 |
| Label Valid Through: | June 2023 |

To the best of our knowledge, the information contained herein is accurate and reliable. However, Micronic does not make any express or implied representation or warranty as to the accuracy or completeness of the information. It is the user's responsibility to inspect and test the Micronic products in order to satisfy itself as to the suitability of the products for its particular purpose. The user is responsible for the appropriate, safe and legal use, processing and handling of the Micronic products.

WHERE TO BUY

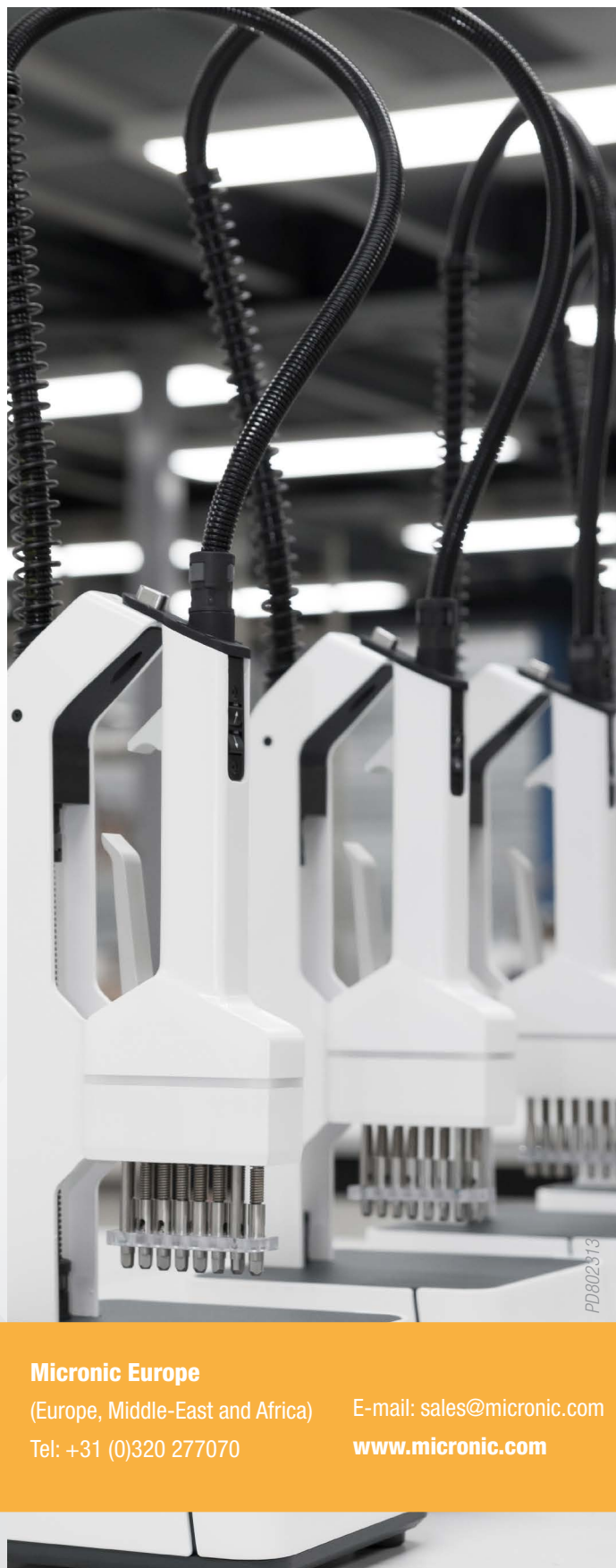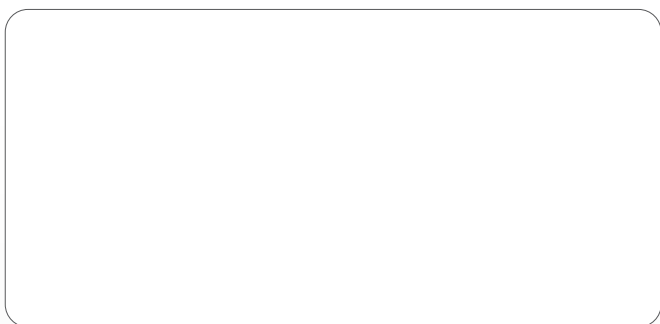

Micronic America

(US, Canada and Latin America)

Tel: +1 484 480 3372

Micronic Asia

(Asia and Pacific)

Tel: +65 85 028 878

Micronic Europe

(Europe, Middle-East and Africa)

Tel: +31 (0)320 277070

E-mail: sales@micronic.com

www.micronic.com